Millipore®

SigmaAldrich.com

User Guide

Millex®-FG Filter Unit (Non-Sterile)

REF

SLFG0250M

English

Français

Italiano

Deutsch

Español

Português

Nederlands

Dansk

Svenska

Latviski

Lietuviškai

Magyar

Česky

Polski

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CE

0459

R_xonly

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

PR05380

Rev 06/23

Date of issue: 2023-06

Symbol

Definition

Symbol

Definition

Non-sterile

Use-by date

Non-pyrogenic

Date of manufacture

Consult instructions for use

Manufacturer

Do not use if package is damaged and consult instructions for use

CE conformity marking

Do not re-use

Federal (USA) law restricts this device to sale by or on the order of a licensed healthcare professional.

REF

Catalogue number

CH

REP

Swiss authorised representative

LOT

Batch code

UDI

Unique device identifier

Temperature limit

Made in Ireland

Merck Millipore Ltd.
Tullagreen, Carrigtwohill,
Co. Cork, Ireland
T45 KD29
an affiliate of Merck KGaA,
Darmstadt, Germany

CH

REP

MDSS CH GmbH
Laurenzenvorstadt 61
5000 Aarau
Switzerland

The initial M, Millipore, Millex, Fluoropore, and Sigma Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2003-2024 Merck KGaA, Darmstadt, Germany and/or its affiliates. All rights reserved.

English

Millex®-FG Filter Unit (Non-Sterile)

SLFG0250M Qty: 1000 Pack

• 25 mm • Non-sterile • Single use only • Non-pyrogenic

Intended Purpose
The Millex®-FG Filter Unit (Non-Sterile) is a medical device that is supplied in a non-sterile state. Once sterilized, it is intended for use as a gas vent filter for ventilation of medical devices. Device is intended for use by medical and/or laboratory professionals only.

Introduction
This filter is intended for use in vacuum line protection, gas filtration, and tank or bioreactor venting where a sterile filter is NOT required. Clinical benefits include the reduction of microorganisms, particles, precipitates, and undissolved powders larger than the membrane's rated pore size from gaseous solutions to limit cross-contamination. Aqueous solutions are blocked from filtration through the hydrophobic membrane of the device. This non-sterile filter is compatible with ethylene oxide (EO) sterilization. Sterilization by any other method (e.g., gamma irradiation, electron beam, steam) is not recommended. For applications where a sterile filter is required, use the sterile equivalent, Cat. No. SLFG025LS.

Chemical Compatibility
This filter is compatible with most gaseous solutions. Compatibility of gaseous solutions should be tested before use. Visit [SigmaAldrich.com/MedMillex](https://www.sigmaaldrich.com/MedMillex) for more information.

WARNINGS

- Sterilize with EO only.
- Filter gaseous solutions with a temperature limit of 45 °C (113 °F).
- Do not use the same vent filter to filter gaseous solutions in both directions.
- Sudden loss of pressure could indicate failure of the filter.
- Single use only; do not re-use.
- This product is not suitable for filtration of aqueous solutions.
- Discard appropriately after single use. See "Disposal" below.

Directions for Use
To use the Millex®-FG Filter with a vacuum or air system, attach the appropriate Luer adapters to the vacuum or air system tubing.

If the Millex®-FG Filter becomes momentarily inoperable due to aqueous fluid completely filling the housing, clear the filter by disconnecting it from the upstream Luer connection and aseptically applying air pressure with a syringe.

If using the Millex®-FG Filter as an air vent, attach a needle to the Luer-slip connection of the filter, then insert the needle through the cap of the container you wish to vent.

Specifications and Performance Characteristics	
Materials	
Membrane	Hydrophobic Fluoropore™ polytetrafluoroethylene (PTFE), type FG
Pore size	0.2 µm
Housing	Polyvinyl chloride (PVC)
Dimensions	
Inlet to outlet	25 mm (0.98 in.)
Diameter	29 mm (1.14 in.)
Filtration area	4 cm² (0.62 in²)
Temperature limit	45 °C (113 °F)
Pressure limit at 21 °C	5.2 bar (75 psi) inlet and differential maximum
Connections	Female Luer-Lok™ inlet; male Luer-slip outlet

Notice
We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

Contact Information
For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

Disposal
Follow precautions for disposal of items contaminated with potentially infectious or hazardous biological material according to all applicable international, federal, state, and local regulations.

Technical Assistance
Visit the tech service page on our web site at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice). Any serious incident of this device should be reported to manufacturer and competent authority of country where user is established.

Standard Warranty
The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

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