

SAFC

3050 Spruce Street
Saint Louis, MO 63103 USA
Phone: 1-800-244-1173 • 1-314-534-4900

Fax: 1-888-289-8449 Global Email: safcglobal@sial.com

Product Information

COMPOZR® KNOCKOUT ZFN KIT - CHO SLC35A1

CATALOG NO. ZFNSLC35A1

Product Information

Zinc Finger Nucleases (ZFNs) are a class of engineered DNA-binding proteins which facilitate targeted genome editing by binding to a user-specified locus ,subsequently causing a double-strand break (DSB). Post cleavage of the DNA, the cell employs endogenous DNA repair processes, either non-homologous end joining (NHEJ) or homology-directed repair (HDR), to repair this targeted DSB. These repair processes can be channeled to generate precisely targeted genomic edits resulting in an cell line or an organism with specific gene disruptions (knockouts), integrations, or modifications. To provide the best potential of creating a functional gene knockout, CompoZr Knockout ZFNs are targeted to the 5' exon sequences which represent the first 2/3 of the coding region for the gene of interest.

Gene Target Background

The protein produced by the SLC35A1 gene resides on the membrane of the Golgi and is responsible for transporting the nucleotide sugar CMP-sialic acid into the Golgi. Knocking out function of this gene in a cell results in the production of asialylated glycan structures due to the loss of available sialic acid as a substrate to downstream sialic acid transferases.


Attachment to a cell surface receptor is a key first step in the infection cycle for many viruses (**Figure 3**). It has been shown that viral capsids derived from Minute Virus of Mice (MVM) bind to specific glycan structures with a terminal sialic acid. Knocking out the function of this gene in a cell results in asialylated glycan structures (**Figure 1**), thus eliminating the ability of MVM to bind to and enter the cell (**Figure 2**). Therefore, a SLC35A1 knockout host cell line can be engineered leveraging the ZFN technology in order to create a MVM resistant cell line.

SAFC has identified a ZFN pair targeting the CHO SLC35A1 gene (Gene ID 100689322). Upon genetic disruption at this loci, asialylated cell lines can be isolated. SAFC has validated this phenotypic affect, as well as resistance to MVM infection , by making SLC35A1 knockouts using the CHOZN GS-/- cell line (Product No. CHOGS).

Kit Components

CompoZr Knockout ZFNs are provided as aliquots (10 per kit) of mRNA encoding for the ZFNs along with plasmid DNA for both ZFN pairs. Also included in the kit are custom-designed PCR primers that allow for determination of the rate of mutation and for screening of individual clones that have the desired mutational event, and a positive control (genomic DNA that has been modified using the CompoZr Knockout ZFN).

Wildtype CHO

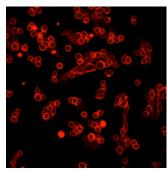

SLC35A1 knockout CHO

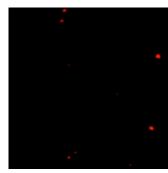


Figure 1. Wildtype and knockout cells were incubated with MAL II-biotin, washed and then incubated with Streptavidin-Alexa-647. Cells were washed and observed under a microscope. SLC35A1 knockout cells do not bind MallI lectin indicating the absence of surface sialic acid.

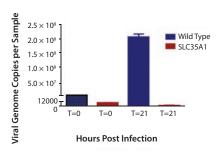


Figure 2. Wildtype and SLC35A1 knockout cells were incubated with MVM for 3 hours and then washed to remove non-specifically bound virus. Samples were then collected at time=0 hours and time=21 hours and analyzed by qPCR using a probe against the MVM NS2 gene. Time 0 indicates levels of viral attachment. Time 21 indicates levels of internalization and single cycle replication.

Features and Benefits

- First commercially available gene editing tool for viral resistance applications
 - Enables the generation of a MVM resistant host CHO cell line
 - Can be used to engineer a host cell line or a producing cell line
- Full kit provided with controls
 - Ten aliquots of transfection ready mRNA
 - Plasmid DNA provided for additional RNA generation or expansion
 - Primers for screening and sequencing
- Positive control DNA for CEL I assay validation
- Technical bulletin provided with detailed protocols for transfection and ZFN activity confirmation
- Technical support provided by the CHOZN R&D scientists
- Custom Cell Line Engineering services are available – inquire at CHOZN@sial.com



SAFC

3050 Spruce Street Saint Louis, MO 63103 USA Phone: 1-800-244-1173 • 1-314-534-4900

Fax: 1-888-289-8449 Global Email: safcglobal@sial.com

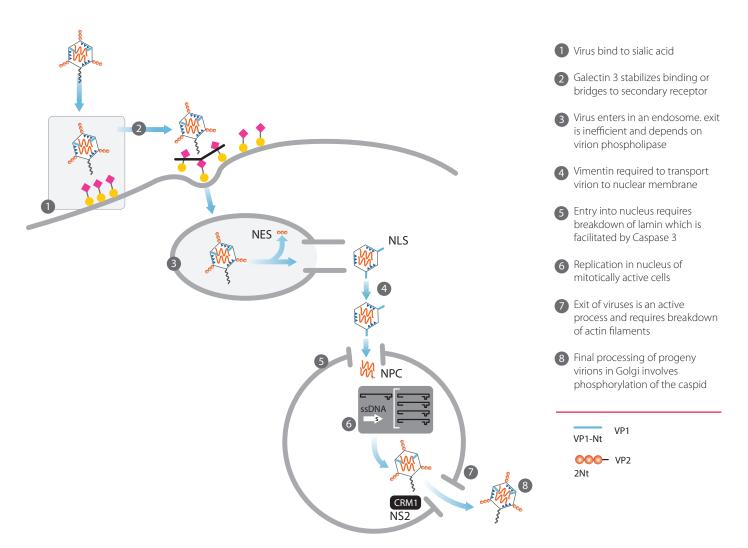


Figure 3. Infection and replication of Minute Virus of Mice in a host cell line

By eliminating the presence of surface sialic acid via ZFN mediated knockout of SLC35A1, MVM is unable to infect and replicate within a host cell line.

ZFN License Agreement

This Product and its use are the subject of one or more of the following patents controlled by Sangamo BioSciences, Inc.: U.S. Patent Nos. 6,534,261, 6,607,882, 6,746,838, 6,794,136, 6,824,978, 6,866,997, 6,933,113, 6,979,539, 7,013,219, 7,030,215, 7,220,719, 7,241,573, 7,241,574, 7,585,849, 7,595,376, 6,903,185, 6,479,626, US20030232410, US20090203140 and corresponding foreign patent applications and patents.

Before opening or using this product, please read the terms and conditions set forth in this license agreement. Your use of this product shall constitute acknowledgment and acceptance of these terms and conditions. If you do not agree to use this Product pursuant to the terms and conditions set out in this

License Agreement, please contact Sigma Technical Services within ten days of receipt to return the unused and unopened Product for a full refund; provided, however, that custom-made Products may not be returned for a refund.

The purchase of this Product conveys to you, the buyer, the non-transferable right to use the purchased Product for Licensed Research Use (see definition below) subject to the conditions set out in this License Agreement. If you wish to use this Product for any purpose other than Licensed Research Use, you must first obtain an appropriate license (see information set out below).

This Product may not be used for any purpose other than Licensed Research Use. Your right to use this Product for Licensed Research Use is subject to the following conditions and restrictions:



SAFC

3050 Spruce Street
Saint Louis, MO 63103 USA
Phone: 1-800-244-1173 • 1-314-534-4900

Fax: 1-888-289-8449 Global Email: safcglobal@sial.com

- 1. "Licensed Research Use" means any use for research purposes, other than:
 - (a) Licensing, selling, distributing, or otherwise providing Modified Animals to any third party other than Sigma and its affiliates as provided herein: provided however, that you may provide Modified Animals to researchers within your research organization located at the same research facility or campus. A "Modified Animal" means an animal having a genomic modification at the target site that results from Customer's use of the Product. Modified Animal includes but is not limited to (a) heterozygotes and mosaic animals, (b) the descendants of Modified Animals, (c) animals created from the breeding of Modified Animals with other animals, and (d) animals created by the Customer which contain and/or incorporate genetic information derived from Modified Animals.
 - (b) GMP production of therapeutic, diagnostic, prophylactic or other medicinal Products intended for use in humans or non-human animals, or any other industrial use solely to the extent involving commercial sale of a Product or service. If a molecule or any derivative of such molecule is used in or administered to humans, then the production of such molecule shall be deemed to be GMP production and therefore in violation of this License Agreement;
 - (c) use for gene targeting and/or gene regulation to modify the genome of a plant cell, plant, or plant cell culture (in each case, whether constituting or derived from a vascular or non-vascular plant), or alter the nucleic acid or protein expression in a plant cell, plant, or plant cell culture. "Non-vascular" plants shall include but not be limited to algae, moss, and fungi; and
 - (d) modification or reverse-engineering of the Product in any way or creating any derivatives or sequence variants thereof.
- 2. You may not transfer the Product, its components, or any materials made through the use of this Product, including Modified Animals, to any third party without prior written approval of Sigma and without the transferee entering into a material transfer agreement with Sigma. Notwithstanding the foregoing:

- (a) the Product or materials made through use of the Product may be transferred by you to your legal affiliates or bonafide third party contractors performing paid work on your behalf, with the exception of creation of Modified Animals, provided the use by such third party contractors is limited to performance of work for you; and
- (b) you may donate Mice that are Modified Animals as defined above ("Modified Mice") to The Jackson Laboratory, a licensed distributor of Modified Mice.
- You may not transfer the Product or materials made through use of the Product to third party contractors performing paid work on your behalf for the purposes of creation of Modified Animals.
- 4. Your right to use the Product will terminate immediately if you fail to comply with these terms and conditions. You shall, upon such termination of your rights, destroy all Product, Modified Animals, and components thereof in your control, and notify Sigma of such in writing.
- You may not use the Product to support the filing of a patent application in any country in the world that contains claims directed to the Product or its uses.

For information on purchasing a license to this Product for purposes other than Licensed Research Use, contact your local Sigma Sales representative, who will refer you to the proper licensing representative, or in the USA call 800-325-3010.

For information on donating Modified Mice to The Jackson Laboratory, visit their website at: http://www.jax.org/grc/index.html.

©2016 Sigma-Aldrich Co. LLC. All rights reserved. SAFC and SIGMA-ALDRICH are trademarks of Sigma-Aldrich Co. LLC, registered in the US and other countries. CompoZr is a registered trademark of Sigma-Aldrich Co. LLC. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see product information on the SAFC website at sigma-aldrich.com/safc and/or on the reverse side of the invoice or packing slip. Sigma-Aldrich Corp. is a subsidiary of Merck KGaA, Darmstadt, Germany.