

### **ProductInformation**

### SIGMA QUALITY CONTROL TEST

# Enzymatic Assay of PROTEIN KINASE CATALYTIC SUBUNIT Sigma Prod. Nos. P-2645 and P-8289 Phosphorylating Activity

### PRINCIPLE:

Casein +  $\gamma$ -<sup>32</sup>P-ATP  $\frac{Protein Kinase}{Catalytic Subunit}$  > [32P]-Phosphorylated Casein + ADP

Abbreviations used:  $\gamma^{-32}$ P-ATP = Adenosine 5'-Triphosphate  $\gamma^{-32}$ P label ADP = Adenosine 5'-Diphosphate

**CONDITIONS:** T = 30°C, pH = 6.5

**METHOD:** Radioactive

#### **REAGENTS:**

A. 1000 mM Potassium Phosphate Buffer, pH 6.5 at 30°C (Prepare 100 ml in deionized water using Potassium Phosphate, Monobasic, Anhydrous, Sigma Prod. No. P-5379. Adjust to pH 6.5 at 30°C with 2 M KOH.)

- B. 5.0% (w/v) Casein Solution (Casein)(Use Casein, from Bovine Milk, 5% (w/v) Solution, Sigma Prod. No. C-4765.)
- C. 500 mM Magnesium Acetate Solution (Mg(OAc)<sub>2</sub>)
   (Prepare 10 ml in deionized water using Magnesium Acetate, Tetrahydrate, Sigma Prod. No. M-9147.)
- D. 250 mM Aminophylline Solution (AP)
   (Prepare 10 ml in deionized water using Aminophylline, Hydrate, Sigma Prod. No. A-1755.)
- E. 330 mM Dithiothreitol Solution (DTT) (Prepare 10 ml in deionized water using DL-Dithiothreitol, Sigma Prod. No. D-0632. PREPARE FRESH.)
- F. 10.0 mM Adenosine 5'-Triphosphate Solution (ATP) (Prepare 1 ml in deionized water using Adenosine 5'-Triphosphate, Disodium Salt, Sigma Prod. No. A-5394. **PREPARE FRESH**.)

Revised: 12/27/94 Page 1 of 4

# Enzymatic Assay of PROTEIN KINASE CATALYTIC SUBUNIT Sigma Prod. Nos. P-2645 and P-8289 Phosphorylating Activity

### **REAGENTS:** (continued)

- G. γ-<sup>32</sup>P-Adenosine 5'-Triphosphate Solution (γ-<sup>32</sup>P-ATP)
   (Use product with minimum radioactive concentration of 30 curies/mmole and 2 millicuries/ml.)
- H. 13.5% (w/v) Trichloroacetic Acid Solution (TCA) (Prepare 20 ml in deionized water using Trichloroacetic Acid, 6.1 N Solution, approximately 100% (w/v), Sigma Stock No. 490-10.)
- Protein Kinase Catalytic Subunit Enzyme Solution
   (Immediately before use, prepare a solution containing 200 400 units/ml of Protein Kinase Catalytic Subunit in cold Reagent J.)
- J. 39 mM DL-Dithiothreitol Solution (Enzyme Diluent)
   (Prepare 10 ml in deionized water using DL-Dithiothreitol, Sigma Prod. No. D-0632.
   PREPARE FRESH.)
- K. 6.75% (w/v) Trichloroacetic Acid Solution (Wash Solution) (Prepare 20 ml in deionized water using Trichloroacetic Acid, 6.1 N Solution, approximately 100% (w/v), Sigma Stock No. 490-10.)
- Methylethyl Cellosolve
   (Prepare by adding equal volumes of Ethylene Glycol Monoethyl Ether, Sigma Prod.
   No. E-2632, to Ethylene Glycol Monomethyl Ether, Sigma Prod. No. E-5378.)
- M. Scintillation Cocktail (Use Sigma-Fluor Universal LSC Cocktail for Aqueous Samples, Sigma Prod. No. S-4273.)

### PROCEDURE:

Prepare a reaction cocktail by pipetting (in milliliters) the following reagents into a suitable container:

Deionized Water	2.00
Reagent A (Buffer)	0.50
Reagent C (Mg(OAc) <sub>2</sub> )	0.25
Reagent D (AP)	0.10
Reagent E (DTT)	0.10
Reagent F (ATP)	0.05

Revised: 12/27/94 Page 2 of 4

# Enzymatic Assay of PROTEIN KINASE CATALYTIC SUBUNIT Sigma Prod. Nos. P-2645 and P-8289 Phosphorylating Activity

**PROCEDURE:** (continued)

Mix by swirling. Transfer 1 ml to a suitable container and add enough Reagent G ( $\gamma$ -<sup>32</sup>P-ATP) to yield approximately 150,000-200,000 counts/minute (cpm) in 0.05 ml of the solution. Then add 0.50 ml of Reagent B (Casein). This is the Reaction Cocktail.

Pipette 0.05 ml aliquots of Reagent I (Enzyme Solution) into a multiwell disposable titerplate. Place in a 4°C ice bath.

Add 0.05 ml of the Reaction Cocktail to each well and mix by air injection. Immediately transfer the titerplate to a 30°C water bath. Incubate at 30°C for 10 minutes. Then add 0.10 ml of Reagent H (TCA) to each well.

Filter the material in the wells through 0.45  $\mu m$  Millipore HA Type filters. Wash 3 times with Reagent K (Wash Solution).

Transfer the filters to suitable 2 dram scintillation vials containing 2.00 ml of Reagent L (Methylethyl Cellosolve). To each scintillation vial, add 5 ml of Reagent M. Count the radioactivity in a suitable scintillation counter.

### **CALCULATIONS:**

The total number of picomoles (pMoles) of ATP in the reaction mixture is calculated as follows:

$$\frac{(0.05) (0.01) (10^9)}{(3.00) (1.5)} = 1.1 \times 10^5 \text{ pmole/ml Reaction Cocktail}$$

Find cpm/pmole by counting 0.05 ml (5555 pmoles) of the Reaction Cocktail.

0.05 = Volume (in milliliters) of ATP used in the Reaction Cocktail
 0.01 = Millimolar concentration of ATP (Reagent F)
 10<sup>9</sup> = Conversion of millimoles to picomoles
 3.00 = Intermediate volume (in milliliters) of Reaction Cocktail
 1.5 = Total volume (in milliliters) of Reaction Cocktail

CPM counted = Actual count - background on filters

Revised: 12/27/94 Page 3 of 4

# Enzymatic Assay of PROTEIN KINASE CATALYTIC SUBUNIT Sigma Prod. Nos. P-2645 and P-8289 Phosphorylating Activity

### **CALCULATIONS:** (continued)

10 = Time of Assay (in minutes) as per the Unit Definition 0.05 = Volume (in milliliters) of Protein Kinase Catalytic Subunit used

#### UNIT DEFINITION:

One unit will transfer 1.0 picomole ( $10^{-12}$  mole) of phosphate from  $\gamma$ -<sup>32</sup>P-ATP to hydrolyzed, partially dephosphorylated casein (C-4765) per minute at pH 6.5 at 30°C.

#### FINAL ASSAY CONCENTRATIONS:

In a 0.10 ml solution, the final assay concentrations are 55 mM potassium phosphate, 14 mM magnesium acetate, 3 mM aminophylline, 4 mM dithiothreitol, 0.055 mM adenosine 5'-triphosphate, 0.83% (w/v) casein, and 10 - 20 units protein kinase catalytic subunit.

#### **REFERENCES:**

Riemann, E.M., Walsh, D.A., and Krebs, E.G. (1971) Journal of Biological Chemistry 246, 1986-1995

Mayer, S.E., Stull, J.T., Wastila, W.B., and Thompson, B. (1974) *Methods in Enzymology*, XXXVIII, Part C, 66-73

### NOTE:

- 1. The concentration of dithiothreitol does not include that contributed by the protein kinase catalytic subunit diluent.
- 2. This assay is based on the cited references.
- 3. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that Sigma products conform to the information in Sigma publications. Purchaser must determine the suitability of the information and products for its particular use. Upon purchase of Sigma products, see reverse side of invoice or packing slip for additional terms and conditions of sale.

Revised: 12/27/94 Page 4 of 4