Safety Sheet

Thermocycler Module

Select language:

English.	2	Italiano (Italian)	30
العربية (Arabic)	4	日本語 (Japanese)	32
Български (Bulgarian)	6	한국어 (Korean)	34
中文 (Chinese, Simplified)	8	Latviski (Latvian)	36
Hrvatski (Croatian)	10	Lietuvių (Lithuanian)	38
Čeština (Czech)	12	Malti (Maltese)	40
Dansk (Danish)	14	Polski (Polish)	42
Nederlands (Dutch)	16	Português (Portuguese)	44
Eesti keel (Estonian)	18	Română (Romanian)	46
Suomi (Finnish)	20	Srpski (Serbian)	48
Français (French)	22	Slovenčina (Slovak)	50
Deutsch (German)	24	Slovenščina (Slovenian)	52
Ελληνικά (Greek)	26	Español (Spanish)	54
Magyar (Hungarian)	28	Svenska (Swedish)	56

Intended Use

The Opentrons Thermocycler Module GEN2 is a fully automated on-deck thermocycler, providing hands-free PCR in a 96-well plate format. Its heated lid and disposable seal fit tightly over the plate, ensuring efficient sample heating and minimal evaporation.

Specifications

Power Requirements

Power Adapter

Input voltage	100-240 VAC (+/- 10%)
Input frequency	50/60 Hz
Input current	8.5 A/115 VAC 5 A/230 VAC
Output voltage	36 VDC
Output current	17.5 A
Output wattage	630 W maximum
Power source Required	C13

Instrument

Outer dimensions	Length	244.95 mm (0.96 inches)
	Width	172 mm (6.77 inches)
	Height	310.1 mm (12.21 inches)
Weight		8.4 kg (18.52 lbs)

Environmental Conditions

For indoor use only. This module is not rated for wet locations or any other outdoor parameter. See the manual for recommended parameters for system operation.

	Acceptable for Operation	Storage/Transport
Ambient Temperature	2 °C to 40 °C 35.6 °F to 104 °F	-10 °C to 60 °C 14 °F to 140 °F
Relative Humidity	30-80% non-condensing (below 30 °C)	10-85% non-condensing (below 30 °C)
Altitude	Up to 2000 m above sea level	Up to 2000 m above sea level
IEC61010-1	Pollution Degree II	
Overvoltage	Category II	
Ingress Protection Rating	IP2X	

Precautions

Please read these safety instructions and manual carefully before using Flex and/or any module. Keep this Safety Sheet near the instrument for reference.

- This module generates enough heat to cause serious burns. Allow maximum clearance to avoid burns.
- The lid seal may be hot when opened.
- Safety goggles should be worn at all times during operation.
- Do not set the plate temperature below the ambient temperature for over two hours. The resulting condensation could damage the module.
- Do not operate near or with explosive substances. Do not use with substances that may generate explosive gases.
- Read this module's Quick Start Guide and Flex manual before operating.
- Unplug when not in use.

Electrical Warnings

Failure to follow any of these warnings may result in injury, fire, and electric shock.

- Plug into a grounded, Class 1 circuit.
- Only use the power cord provided.
- Do not use a power cord with inadequate current or voltage ratings. Do not replace the power cord unless at the direction of Opentrons Support.
- Keep the power cord free of obstructions so you can remove it if necessary.
- Do not connect (plug in), disconnect (unplug), or use AC power cables if:
 - The cable is frayed or damaged.
 - Other attached cables, cords, or receptacles are frayed or damaged.
- Using damaged power cords can cause an electric shock hazard resulting in serious injury or damage to the robot.
- Do not replace the AC power cable unless at the direction of Opentrons Support.
- Keep the power cord free of obstructions so you wcan remove it if necessary.
- Do not attempt to repair or remove the outer case of the module or its power supply unless directed by Opentrons Support.

Maintenance

- Power OFF and unplug the instrument before beginning cleaning, to prevent risk of electrical shock.
- No internal parts require cleaning. The modules should never be disassembled for cleaning.
- Do not use acetone. This may cause damage or dissolve parts of the instrument.
- See the manual, Cleaning section for detailed cleaning instruction.

Technical Service

Do not attempt to modify, revise, or repair the instrument. Only authorized Opentrons Support may service the instrument. Unauthorized repair or service may void the warranty and/or cause hazardous conditions.

Installation

- Read the manual for detailed installation directions before installing any module.
- Ensure the module is properly oriented in the caddy sized for this module.
- Firmly connect all cords to the module, as directed in the manual.
- The mains power cord and plug must be easily accessible in the case of a hazard in which the cord and plug need to be pulled to remove power from the product.

Equipment Operation

Read the manual before operating.

Power ON/OFF

Plug into the wall outlet. From the front of the power supply, press (-) on the left side of the power switch to turn on.

To turn off, press (o) the right side of the power switch. Unplug power cord from outlet when not in use.

Emergency Disconnect

Push down the Emergency Stop Pendant (E-stop) to quickly stop robot motion. The Flex requires a functional, disengaged E-stop to be attached at all times. The E-stop does not cut power.

In case of an emergency, unplug the power cord from the power source. Contact Opentrons Support before restarting the robot.

Symbol Definitions

	CAUTION, documentation must be consulted
	CAUTION, possibility of electrical shock.
	CAUTION, hot surface
	CAUTION, risk of pinching
	CAUTION, risk of explosion!
	Manufacturer/ Legal entity
	This product has been evaluated by an OSHA certified Nationally Recognized Testing laboratory (NRTL) in the United States. Conforms to UL Std 61010-1 Certified to CSA Std 61010-1
	European Council Directives conformity marking.
	United Kingdom conformity marking

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© OPENTRONS 2025
Opentrons Labworks, Inc.
Opentrons Flex™ (Opentrons Labworks, Inc.) Registered names, trademarks,