

Product Information

Lectin from *Galanthus nivalis* (snowdrop)

Catalog Number **L8275**

Storage Temperature 2–8 °C

Synonym: GNL

Product Description

Molecular mass: This 52 kDa lectin complex is made up to 4 identical 13 kDa subunits.

Extinction coefficient ($E_{1\%}$): 1.8 at 280 nm (in water)

Lectins are proteins or glycoproteins of non-immune origin that agglutinate cells and/or precipitate complex carbohydrates. Lectins are capable of binding glycoproteins even in presence of various detergents.¹ The agglutination activity of these highly specific carbohydrate-binding molecules is usually inhibited by a simple monosaccharide, but for some lectins, di, tri, and even polysaccharides are required.

Lectins are isolated from a wide variety of natural sources, including seeds, plant roots and bark, fungi, bacteria, seaweed and sponges, mollusks, fish eggs, body fluids of invertebrates and lower vertebrates, and from mammalian cell membranes. The precise physiological role of lectins in nature is still unknown, but they have proved to be very valuable in a wide variety of applications *in vitro*, including:

- blood grouping and erythrocyte polyagglutination studies
- mitogenic stimulation of lymphocytes
- lymphocyte subpopulation studies
- fractionation of cells and other particles
- histochemical studies of normal and pathological conditions

Sigma offers a range of lectins suitable for the above applications. Most Sigma lectins are highly purified by affinity chromatography, but some are offered as purified or partially purified lectins, suitable for specific applications.

Many of the lectins are available conjugated to (conjugation does not alter the specificity of the lectin):

- fluorochromes (for detection by fluorimetry)
- enzymes (for enzyme-linked assays)
- insoluble matrices (for use as affinity media)

Please refer to the table for general information on the most common lectins.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

This lectin is soluble in phosphate buffered saline, pH 6.8, (1 mg/ml).

Storage/Stability

Store the product at 2–8 °C.

In solution, aggregation is thought to occur in the presence of high concentrations of 2-mercaptoethanol.

Lectin	Molecular Mass (kDa)	Subunits	Blood Group	Specificity Sugar	Mitogenic Activity
<i>Abrus precatorius</i>			—	gal	+
Agglutinin	134	4		gal	
Abrin A (toxin)	60	2		gal	
Abrin B (toxin)	63.8	2(αβ)		gal	
<i>Agarius bisporus</i>	58.5	—	—	β-gal(1→3)galNAc	
<i>Anguilla anguilla</i>	40	2	H	α-L-Fuc	
<i>Arachis hypogaea</i>	120	4	T	β-gal(1→3)galNAc	
<i>Artocarpus integrifolia</i>	42	4	T	α-gal→OMe	+
<i>Bandeiraea simplicifolia</i>					
BS-I	114	4	A, B	α-gal, α-galNAc	
BS-I-A ₄	114	4	A	α-galNAc	
BS-I-B ₄	114	4	B	α-gal	
BS-II	113	4	acq, B, Tk, T	glcNAc	
<i>Bauhinia purpurea</i>	195	4	—	β-gal(1→3)galNAc	+
<i>Caragana arborescens</i>	60; 120 ^a	2/4	—	galNAc	
<i>Cicer arietinum</i>	44	2	—	fetuin	
<i>Codium fragile</i>	60	4	—	galNAc	
<i>Concanavalin A</i>	102	4	—	α-man, α-glc	+
<i>Succinyl-Concanavalin A</i>	51	2	—	α-man, α-glc	+ ^b
<i>Cytisus scoparius</i>	—	—	—	galNAc, gal	
<i>Datura stramonium</i>	86	2(αβ)	—	(glcNAc) ₂	
<i>Dolichos biflorus</i>	140	4	A ₁	α-galNAc	
<i>Erythrina corallodendron</i>	60	2	—	β-gal(1→4)glcNAc	+
<i>Erythrina cristagalli</i>	56.8	2(αβ)	—	β-gal(1→4)glcNAc	
<i>Euonymus europaeus</i>	166	4(αβ)	B, H	α-gal(1→3)gal	+
<i>Galanthus nivalis</i>	52	4	(h)	non-reduc. α-man	
<i>Glycine max</i>	110	4	—	galNAc	+ ^c
<i>Helix aspersa</i>	79	—	A	galNAc	
<i>Helix pomatia</i>	79	6	A	galNAc	
<i>Lathyrus odoratus</i>	40-43	4(αβ)	—	α-man	+
<i>Lens culinaris</i>	49	2	—	α-man	+
<i>Limulus polyphemus</i>	400	18	—	NeuNAc	
Bacterial agglutinin	—	—	—	galNAc, glcNAc	
<i>Lycopersicon esculentum</i>	71	—	—	(glcNAc) ₃	
<i>Maackia amurensis</i>	130	2(αβ)	O	sialic acid	+
<i>Maclura pomifera</i>	40-43	2(αβ)	—	α-gal, α-galNAc	
<i>Momordica charantia</i>	115-129	4(αβ)	—	gal, galNAc	
<i>Naja mocambique mocambique</i>	—	—	—	—	
<i>Naja naja kaouthia</i>	—	—	—	—	
<i>Narcissus pseudonarcissus</i>	26	2	(h)	α-D-man	
<i>Persea americana</i>	—	—	—	—	
<i>Phaseolus coccineus</i>	112	4	—	—	
<i>Phaseolus limensis</i>	247(II)	8	A	galNAc	+
	124(III)	4			
<i>Phaseolus vulgaris</i>					
PHA-E	128	4	—	oligosaccharide	+
PHA-L	128	4	—	oligosaccharide	+
PHA-P					
PHA-M					

----- Table continued on next page -----

Lectin	Molecular Mass (kDa)	Subunits	Blood Group	Specificity Sugar	Mitogenic Activity
<i>Phytolacca americana</i>	32	—	—	(glcNAc) ₃	+
<i>Pisum sativum</i>	49	4(αβ)	—	α-man	+
<i>Pseudomonas aeruginosa PA-I</i>	13-13.7	—	—	gal	+ ^c
<i>Psophocarpus tetragonolobus</i>	35	1	—	galNAc, gal	
<i>Ptilota plumosa</i>	65; 170	—	B	α-gal	
<i>Ricinus communis</i>					
Toxin, RCA ₆₀	60	2	—	galNAc, β-gal	
Toxin, RCA ₁₂₀	120	4	—	β-gal	
<i>Sambucus nigra</i>	140	4(αβ)	—	αNeuNAC(2→6)gal	+ ^c
				galNAc	
<i>Solanum tuberosum</i>	50; 100 ^a 1, 2		—	(glcNAc) ₃	
<i>Sophora japonica</i>	133	4	A, B	β-galNAc	
<i>Tetragonolobus purpureas</i>	120(A) 58(BA) 117(C)	4 2 4	H H H	α-L-fuc α-L-fuc α-L-fuc	
<i>Triticum vulgaris</i>	36	2	—	(glcNAc) ₂ , NeuNAc	+
<i>Ulex europaeus</i>					
UEA I	68	—	H	α-L-fuc	
UEA II	68	—	—	(glcNAc) ₂	
<i>Vicia faba</i>	50	4(αβ)	—	man, glc	+
<i>Vicia sativa</i>	40	4(αβ)	—	glc, man	+
<i>Vicia villosa</i>	139	4	A ₁ +T _n	galNAc	
A ₄	134	4	A ₁	galNAc	
B ₄	143	4	T _n	galNAc	
<i>Vigna radiata</i>	160	4	—	α-gal	
<i>Viscum album</i>	115	4(αβ)	—	β-gal	
<i>Wisteria floribunda</i>	68	2	—	galNAc	

^a Concentration-dependent molecular weight

^b Non-agglutinating and mitogenic

^c Mitogenic for neuraminidase-treated lymphocytes

References

1. Nagata, Y., and Burger, M.M., Wheat germ agglutinin. Molecular characteristics and specificity for sugar binding. *J. Biol. Chem.*, **249**, 3116-3122 (1974).

IRB,MWM,JRC,NSB,SAG,MAM 02/15-1