MILLIPORE

CHEMICON now part of Millipore

CHEMISCREEN[™] MEMBRANE PREPARATION RECOMBINANT HUMAN β₃ ADRENOCEPTOR

CATALOG NUMBER:	HTS159M	QUANTITY:	200 units
LOT NUMBER:		VOLUME/CONCENTRATION:	1 mL, 2 mg/mL
BACKGROUND:	The beta adrenergic rece epinephrine, by coupling predominantly in heart, ti adipose β_3 results in lipo is associated with weight lacking the β_3 -adrenocep (Revelli <i>et al.</i> , 1997). The treatments. Millipore's β preparations made from GPCR surface expression adrenoceptor interaction lodocyanopindolol (ICYF [¹²⁵ I]-(-)ICYP, a greater th	eptors mediate the effects of endog to G_s to stimulate cAMP. Whereas he β_3 receptor is found primarily in lysis and thermogenesis. A polymo t gain in obese patients (Clement e otor display increased total body fat ese observations indicate that β_3 is $_3$ adrenoceptor membrane prepara our proprietary stable recombinant on; thus, they are ideal HTS tools for s. The membrane preparations ex P). With 10 µg/well β_3 Adrenocepto han 8-fold signal-to-background rate	genous catecholamines, such as s β_1 and β_2 are found adipose tissue. Activation of orphism in the human gene for β_3 <i>et al.</i> , 1995). In addition, mice t, particularly on a high fat diet a possible target for obesity itions are crude membrane c cell lines to ensure high-level of or screening of antagonists of β_3 hibit a Kd of 0.69 nM for [¹²⁵ I]-(-) or Membrane Prep and 0.75 nM tio was obtained.

APPLICATIONS:

Radioligand binding assay





MILLIPORE

CHEMICON now part of Millipore

10ug/well

Chem1-WT @ 10ug/well



Figure 2. Competition binding for β_3 adrenoceptor. β_3 adrenoceptor Membrane Preparation (10µg/well) or Wild-Type Chem-1 membrane preparation (Millipore catalog # HTS000MC1) was incubated with 0.75 nM [¹²⁵]-(-)ICYP and increasing concentrations of unlabeled SR59230A, and more than 8-fold signal:background was obtained.

Table 1. Signal:background and specific binding values obtained in a competition binding assay with varying amounts of β_3 adrenoceptor membrane prep.

	10 μg/well
Signal:background	13.3
Specific binding (cpm)	11074.7

SPECIFICATIONS: 1 unit = 10 μ g membrane preparation Bmax: 1.3 pmol/mg K_d: ~0.69 nM

Species: Full-length human ADRB3 cDNA encoding the β_3 adrenoceptor (Accession number NM_000025)

HOST CELLS: Chem-1, a adherent mammalian cell line without any endogenous β_3 adrenoceptor expression.

RECOMMENDED ASSAY CONDITIONS: Membranes are mixed with radioactive ligand and unlabeled competitor (see Figures 1 and 2 for concentrations tested) in binding buffer in a nonbinding 96-well plate, and incubated for 1-2 h. Prior to filtration, an FC 96-well harvest plate (Millipore cat. # MAHF C1H) is coated with 0.33% polyethyleneimine for 30 min, then washed with 50mM HEPES, pH 7.4. Binding reaction is transferred to the filter plate, and washed 3 times (1 mL per well per wash) with Wash Buffer. The plate is dried and counted.

Binding buffer: 50 mM Hepes, pH 7.4, 5 mM MgCl_2, 1 mM CaCl_2, filtered and stored at 4 $^{\circ}\!\!\mathrm{C}$

Radioligand: [¹²⁵I]-(-) lodocyanopindolol (Perkin Elmer # NEX189)

Wash Buffer: 50 mM Hepes, pH 7.4, 500mM NaCl, filtered and stored at 4 $\!\!\!^{\circ}\!\!\!^{\circ}\!\!\!^{\circ}$.

MILLIPORE

CH	ΗEΛ	Λ	CON
now	part	of	Millipore

	One package contains enough membranes for at least 200 assays (units), where an unit is the amount of membrane that will yield greater than 8-fold signal:background with ¹²⁵ I-labeled (-)Iodocyanopindolol at 0.75 nM.
PRESENTATION:	Liquid in packaging buffer: 50 mM Tris pH 7.4, 10% glycerol and 1% BSA with no preservatives. Packaging method: Membranes protein were adjusted to the indicated concentration in packaging buffer, rapidly frozen, and stored at -80°C.
STORAGE/HANDLING:	Maintain frozen at -70° for up to 2 years. Do not freeze and thaw.
REFERENCES:	Clement K <i>et al.</i> (1995) Genetic variation in the beta3-adrenergic receptor and an increased capacity to gain weight in patients with morbid obesity. <i>N. Engl. J. Med.</i> 333: 352-354.
	Revelli JP <i>et al.</i> (1997) Targeted gene disruption reveals a leptin-independent role for the mouse beta3-adrenoceptor in the regulation of body composition. <i>J. Clin. Invest.</i> 100: 1098-1106.
	For research use only; not for use as a diagnostic.

Unless otherwise stated in our catalog or other company documentation accompanying the product(s), our products are intended for research use only and are not to be used for any other purpose, which includes but is not limited to, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses or any type of consumption or application to humans or animals.

©2007-2008: Millipore Corporation. All rights reserved. No part of these works may be reproduced in any form without permission in writing.