Safety Sheet

Auto2D[®] Two Dimensional Electrophoresis Device and Consumables

BM-100

Translations available on the following pages:

English	2	한국어 (Korean)	30
Български (Bulgarian)	4	Latviski (Latvian)	32
中文 (Chinese)	6	Lietuviškai (Lithuanian)	34
Hrvatski (Croatian)	8	Norsk (Norwegian)	36
Čeština (Czech)	10	Polski (Polish)	38
Dansk (Danish)	12	Português (Portuguese)	40
Nederlands (Dutch)	14	Română (Romanian)	42
Eesti keel (Estonian)	16	Srpski (Serbian)	44
Français (French)	18	Slovenský (Slovak)	46
Deutsch (German)	20	Slovenščina (Slovenian)	48
Ελληνικά (Greek)	22	Español (Spanish)	50
Magyar (Hungarian)	24	Svenska (Swedish)	52
Italiano (Italian)	26	Türkçe (Turkish)	54
日本語 (Japanese)	28		

Intended Use

The Auto2D® Electrophoresis Device (Auto2D® Device) separates proteins utilizing their specific isoelectric point differences (First Dimension) and their molecular weight differences (Second Dimension) and will be able to separate several hundreds of proteins at a time.

Specifications

Atmospheric Pressure	70–106 kPa (525–795 mmHg)
Rated Voltage	100–240 VAC
Rated Power	200 VA
Rated Frequency	50/60 Hz

Storage and Stability

Auto2D® Device

Operating Temperature	5 °C–35 °C
Operating Humidity	35%–85% (without condensing)
Operating Conditions	- Indoor use only - Maximum altitude 2,000 meters - Overvoltage Category II - Pollution Degrees 2
Storage Temperature	15 °C–30 °C

Electrode Chip and Electrode Chip Plus

Storage Temperature 15 °C–30 °C
If you heat to dry do not exceed 40 °C. Keep away from direct sunlight and high temperature. Use included case for storage.

PAGE Chip

Storage Temperature 2 °C–8 °C
Do not place near the cold air outlet of the refrigerator. There is a risk of freezing.

IEF Chip

Storage Temperature -20 °C

Solution Chip and Solution Chip Plus

Storage Temperature 15 °C–30 °C
Keep away from direct sunlight and high temperature. Use included case for storage.

Precautions

Electrical Warnings

- Do not immerse Auto2D® Device body in liquid or allow liquid to enter any part of the Auto2D® Device. In the case liquid is spilled into the tray, immediately unplug the power cord from the power outlet and clean inside of the tray.
- To avoid danger of electrical shock or injury, do not touch inside of the tray or high voltage probe at the depth of the tray while the Auto2D® Device is operation.
- To avoid potential shock and/or fire hazards please be sure to utilize the correct country plug and power cord for your region. Verify the that the main input voltage in which the product is connected to matches the ratings on the back of the unit before operating. See the rating label on the back of the Auto2D® Device for the voltage.
- Do not plug or unplug the power plug with wet hands. Do not damage, disassemble or alter power cord. Do not put heavy material on, pinch, or bend the power cord forcefully. Be sure to clean the dust off of the plug with a lint free dry cloth before use.
- Do not operate the Auto2D® Device with wet hands.
- If the Auto2D® Device becomes hot, smokes, smells unusual turn the power off and unplug the power cord immediately and contact technical service at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).
- Do not plug in or operate the Auto2D® Device if the power cord is damaged. This may result in injury, fire, and electric shock.
- To disconnect the power supply to the Auto2D® Device, either unplug the power cord from the power outlet or disconnect the power cord at the AC inlet of the Auto2D® Device. Do not pull on the power cord to unplug.
- Install Auto2D® Device so operating the disconnecting device (power switch) is easy and accessible.
- Failure to follow any of these may result in injury, fire, and electric shock.

Tray

- To avoid danger of electrical shock or injury, do not touch inside of the tray while the Auto2D® Device is operating. There is a high voltage probe at the depth of the tray.
- Do not open the tray cover during the operation of the Auto2D® Device.
- Do not allow inappropriate weight or force to be applied to the tray. Only use the Auto2D® accessories on the opened tray. Applying force or accidental impact could damage the tray.

IEF Chip

- Be sure to use rubber or plastic gloves when handling. If irritation occurs on your skin, immediately consult a physician.
- When handling the IEF Chip, follow general instructions for hazard prevention and safety instructions when handling chemicals.
- Only use Auto2D® IEF Chips with the Auto2D® Device. Failure to do so may result in damage to the Auto2D® Device.
- Do not place fingers near tray hinge when loading the IEF Chip into the Auto2D® Device.
- Let IEF chip adapt to room temperature before use.
- When opening IEF chip with scissors or a cutter, do not damage the gel attached to the tip. Also, be careful not to touch the gel directly during handling.

Solution Chip

Be sure to use rubber or plastic gloves when handling. If irritation occurs on your skin, immediately consult a physician.

Electrode Chip

- Be sure to use rubber or plastic gloves when handling. If irritation occurs on your skin, immediately consult a physician.
- Do not apply strong force or shock to the chip or the platinum part of the chip or there is a risk of damage.
- If the Electrode Chip platinum wire is cut, it will be a stabbing hazard and may cut you. Discontinue use of Electrode Chip with broken wire and contact technical service at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice) for additional help.
- Do not hold this chip too firmly as it may cut your hands.

PAGE Chip

- Fluid is enclosed with the gel. If fluid makes contact with eyes, immediately rinse with water and consult a doctor.
- Let the PAGE chip adapt to room temperature before use.

Maintenance and Service

Technical Service

Only authorized service personnel may service the Auto2D® Device. Unauthorized repair or service may void the warranty. Visit tech service at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Product Care And Handling

- Use only Auto2D® Device accessories. Do not attempt to make any technical alterations. Do not dismantle the instrument any further than is described in the user guide. If the instrument is broken or dropped, do not try to disassemble the instrument. Go to [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice) for assistance.
- Follow general instructions for hazard prevention and safety instructions when handling chemicals.
- Only operate the Auto2D® Device on a level and stable surface. Failure to do so may cause liquid to leak into the Auto2D® Device, resulting in injury or damage to the Auto2D® Device.

Cleaning

Be sure to unplug the power cord of Auto2D® Device before cleaning.

Auto2D® Device Exterior

- The Auto2D® Device is NOT autoclavable. Extreme heat will damage the display screen and other electronic components.
- The instrument body and control buttons can be sanitized by wiping with a soft cloth moistened with 70% ethanol.
CAUTION: When sanitizing, make sure that no liquid enters any part of the Auto2D® Device.

Electrophoresis Chamber

Please contact Technical Service at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice) before attempting to clean. Disassembling and cleaning the chamber may result in malfunction of the Auto2D® Device.

LCD Touch Panel

- Do not clean the display screen with sanitizing agents or other aggressive solutions. Wipe the screen with a soft, dry, nonabrasive cloth.
- To clean heavy dirt, please contact our Technical Service at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Fan Filter

Removed dirt on the filter with soft brush.

End Of Life Instructions
WEEE Directive

In accordance with European Union directive on the management of waste electrical and electronic equipment (WEEE), this product must not be disposed of in unsorted municipal waste at the end of its life. It must be taken to a collection and recycling center. For further information, go to [SigmaAldrich.com/weee](https://www.sigmaaldrich.com/weee).

This product utilizes consumable materials for its intended use. Any safety critical information regarding the consumable materials is in the Precaution section of this document.

User Guides

User guides and other documents are available online. Go to [SigmaAldrich.com](https://www.sigmaaldrich.com), enter the catalogue number in the search bar, select your product, and scroll to the documents section.

Symbol Definitions

	Danger. Warning. Caution Indicates the need for the user to consult the instructions for use for important cautionary information such as warnings and precautions that cannot, for a variety of reasons, be presented on the device itself.
	Affixed in accordance with the most recent European Council Directives which apply to this instrument. Please refer to the Declaration of Conformity for further details.
	Catalogue Number
	Manufacturer
	Serial Number

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The life science business of Merck operates
as MilliporeSigma in the U.S. and Canada.

The initial M, Millipore, Auto2D, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2021 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.