

AGGRECAN RECOMBINANT INTERGLOBULAR DOMAIN T₃₃₁ – G₄₅₈ (Aggrecan-IGD1)

CATALOG NUMBER:	CC1890	QUANTITY:	50 μg
LOT NUMBER:		CONCENTRATION:	XX mg/mL
BACKGROUND:	Aggrecan is a large aggregating discs and tendons [1,2]. The ag 130 glucosaminoglycan chains mass can reach $2.2 - 3.0 \times 10^6$ Within the aggrecan molecule 3 Domains G1 and G2 are connected domain (IGD), while the sequent regions for keratan sulfate and c G1 domain with hyaluronan and contain up to 50-100 aggrecant chain through 2 link proteins [1,7], which endowes cartilage with red Degradation of aggrecan appear molecules without the G3 doma aggrecan at 4 sites within the G1481, E1771 – A1772, E1871 – L1872) Cleavage at the latter site had breakdown products in rheumat To measure aggrecanase activitint region was first used b	g proteoglycan of articular carti- grecan core protein consists of are attached to the core pro- Daltons [4]. global domains G1, G2 and G cted by a rod-shaped polypept ice between domains G2 and G chondroitin sulfate chains. Agg link protein to form large aggre monomers noncovalently bour 2,4]. The aggregates form a hy esistibility to compression and G ars to initiate at the C-terminus ain increases with aging [5]. Is chrondroitin sulfate-rich regi) and 1 site within the interglol d been documented by analy oid and osteoarthritis [7]. ity, an artificial recombinant pu- nking FLAG-sequence and by Hughes et al. [8].	ilage. It is also found in aorta, of 2317 amino acids [2]. Up to obtain and the total molecular G3 can be distinguished. ide called interglobular G3 contains attachment precan interacts via the egates. Such aggregates can do to a single hyaluronan ydrated gel-like structurem, deformation. 5. The population of aggrecan solated aggrecanases cleave on (sites $E_{1667} - G_{1668}$, E_{1480} - bular domain ($E_{373} - A_{374}$) [6]. sis of cartilage proteoglycan rotein composed of aggrecan human immunoglobulin G1
DESCRIPTION:	Molecular form: The polypepti $(T_{331} - G_{458})$ is expressed in E contains cleavage sites for agg $-F_{342}$). It comprises the followinT A E D F V D I P E N F F G V GA R G S V I L T V K P I F E V SG E A T R P W G F P T P G L G(His-tag)Main cleavage sites are indicated15 493 Da.Purity: The recombinant aggreed21 kDa in SDS-PAGE. It represed	ide connecting human aggrec <i>E. coli</i> with a C-terminal His-t recanases ($E_{373} - A_{374}$) and mag amino acids: G G E E D I T V Q T V <u>T</u> W P I P S P L E P E E P F T F A P I S P A T A F T S E D L V V Q V ed by arrows. The calculated M can interglobular domain appe ents more than 90% of total pr	an globular domains 1 and 2 ag. The recombinant protein atrix metalloproteinases (N_{341} D M E L P L P R N I T E G E E I G A T A F A E V E N E T 'T A V P G Q P H L P G G M_r of the His-tagged protein is ars as a major band at about otein in the preparation.
APPLICATIONS:	Aggrecan interglobular domai metalloproteinases. For protein proteinase for various time inter Thereafter, aliquots of the incub Upon cleavage with aggrecan SDS-PAGE is reduced from 2	n is used as substrate for hase activity measurements, t vals. hation mixture are analyzed by ases the apparent M _r of agg 21 kDa to about 13 kDa. G	aggrecanases and matrix the protein is incubated with SDS-PAGE or by ELISA. recan interlobular domain in Quantitative measurement of



		aggrecanase cleavage requires a neoepitop antibody with specificity for the N-terminus A R G S V I L T appearing upon hydrolysis. The fragment with the newly formed N-terminus is fixed by the neoepitop antibody to a microplate and quantified with an anti-Histag antibody. In analogy, cleavage by matrix metalloproteinases can be measured with antibodies to neoepitopes appearing upon action of these enzymes.	
PRESENTATIO	N:	The calculated M_r of the His-tagged protein is 15 493 Da. The protein is solubilized in 50 mM Tris-HCI, pH 7.5, 150 mM NaCl, 5 mM CaCl ₂ .	
STORAGE/HAN	IDLING:	MT5-MMP is stable until the expiry date given on the label of stored at -70°C. The protein can be kept at -20 °C for several weeks and on ice for several days. Repeated freezing and thawing should be avoided.	
REFERENCES:		 Knudsen, C.B. and Knudsen, W. (2001) Seminars Cell & Developm. Biol. <u>12</u>, 69-78. Watanabe, H., Yamada, Y. And Kimata, K. (1998) J. Biochem. <u>124</u>, 687-693. Doege, K.J., Sasaki, M., Kimura, T. and Yamada, Y. (1991) J. Biol. Chem. <u>266</u>, 894-902. Hardingham, T.E. and Fosang, A.J. (1992) FASEB J. <u>6</u>, 861-870. Dudhia, J. Davidson, C.M., Wells, T.M., Vynios, D.H., Hardingham, T.E. and Bayliss, M.T. (1996) J. Biochem. <u>313</u>, 933-940. Tortorella, M.D. Pratta, M., Lin, RQ., Austin, J., Ross, O.H., Abbaszade, I., Burn, T. and Arner, E. (2000) J. Biol. Chem. <u>275</u>, 18566-18573. Lohmander, L.S., Neame, P.J. and Sandy, J.D. (1993) Arthritis Rheum. <u>36</u>, 1214-1222. Hughes, C.E., Buttner, F.H., Eidenmuller, B., Caterson, B. and Bartnik, E. (1997) J. Biol. Chem. <u>272</u>, 20269-20274. 	
Important Note:	During shipm products with centrifuging ta	lent, small volumes of product will occasionally become entrapped in the seal of the product vial. For I volumes of 200 μ L or less, we recommend gently tapping the vial on a hard surface or briefly he vial in a tabletop centrifuge to dislodge any liquid in the container's cap.	

FOR RESEARCH USE ONLY; NOT FOR USE IN DIAGNOSTIC PROCEDURES. NOT FOR HUMAN OR ANIMAL CONSUMPTION

Unless otherwise stated in our catalog or other company documentation accompanying the product(s), our products are intended for research use only and are not to be used for any other purpose, which includes but is not limited to, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses or any type of consumption or application to humans or animals.

©2002 - 2011: Millipore Corporation. All rights reserved. No part of these works may be reproduced in any form without permission in writing.