


## Sigma-Aldrich 80956 Timestrip® 12 Hours

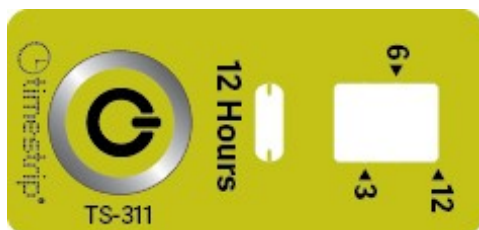


The Timestrip® 12 Hours indicator monitors time elapsed over a period of twelve hours. It is designed to be used at room temperature conditions. It is inert until activated and can be stored at room temperature until needed. It is self-adhesive, single use and irreversible. It has a pre-activation shelf life of two years from shipping.

### Instruction for use



**To Activate:** Fully squeeze the button on top of the indicator. A red line will appear to show that the indicator has been activated.

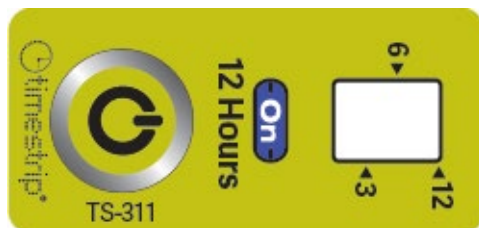

**Activated:** Peel off the liner and stick the indicator to the product or device to be monitored.



**Review:** The red dye will move through the viewing window and past the time markers. When the red dye reaches the end of the viewing window, one month has passed.

### Legal information

Timestrip is a registered trademark of Timestrip Limited

For full instructions, testing protocols, specifications and other information on this product, please visit [www.sigmaaldrich.com/timestrip](http://www.sigmaaldrich.com/timestrip)

The vibrant M, Supelco, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. Detailed information on trademarks is available via publicly accessible resources.  
© 2018 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the US and Canada.

