

3050 Spruce Street, St. Louis, MO 63103 USA
Tel: (800) 521-8956 (314) 771-5765 Fax: (800) 325-5052 (314) 771-5757
email: techservice@sial.com sigma-aldrich.com

Product Information

IgG from rat serum

reagent grade

Catalog Number I4131

Storage Temperature 2-8 °C

Product Description

Purified rat IgG is isolated from pooled normal rat serum by fractionation and ion-exchange chromatography.

Purified rat IgG may be used as a reference antigen, standard, blocking agent, or coating protein in a variety of immunoassays including ELISA, dot immunobinding, Western immunoblotting, immunodiffusion, and immunoelectrophoresis. Other applications include starting materials for the preparation of immunogens and solid phase immunoadsorbents.

Reagent

Supplied as an essentially salt-free, lyophilized powder.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

The product is essentially a salt-free preparation and should be reconstituted to a minimum of 1 mg/ml with 150 mM NaCl.

Protein concentration should be determined after reconstitution by measuring the absorbance at 280 nm ($E_{280}^{1\%}$ = 14).

Storage/Stability

Prior to reconstitution, store at 2–8 °C. After reconstitution, store in working aliquots at –20 °C. Repeated freezing and thawing is not recommended.

Product profile

Purity: ≥95% (SDS-PAGE)

ID by Immuno-Electrophoresis: Single major arc of precipitation in the gamma region versus anti-rat whole serum and anti-rat IgG. Minor gamma arcs may be present versus anti-rat whole serum

RC.PHC 06/12-1