

Product Information

Chorionic Gonadotropin Human

CG5

Storage Temperature: -20 °C

CAS RN: 9002-61-3

Synonym: hCG, Choriogonin

Product Description

pI: 2.95

Extinction Coefficient:³ $\epsilon^M = 1.41 \times 10^4$ (280 nm)

E1% = 3.6 (280 nm)

Human chorionic gonadotropin (hCG) is a glycoprotein hormone produced by the chorionic tissue of the placenta. The glycoprotein hormone family includes luteinizing hormone (LH), follicle-stimulating hormone (FSH), and thyroid-stimulating hormone (TSH), in addition to hCG. The function of hCG is to maintain the corpus luteum and stimulate steroid secretion from the ovary in the beginning stages of gestation. hCG appears in the blood and urine during the first trimester of early pregnancy, and levels decrease thereafter. It has been used for superovulation in animals.²

hCG consists of an α subunit of 92 amino acids and a β subunit of 145 amino acids. The α subunit is common among the family of glycoprotein hormones, whereas the hormone-specific β subunit, which exhibits different degrees of homology, may confer biologic specificity of the individual hormone. The amino acid sequences of the α subunit^{3,4} and the β subunit^{4,5} and the crystal structure of hCG⁷ have been reported. NMR studies of the α subunit have been reported.^{7,8}

The molecular mass of hCG is ~37.9 kDa, with ~31% carbohydrate by weight. The theoretical molecular mass of 37.9 kDa is based on the native form, which contains 2 subunits. The α subunit has a molecular mass of 14.9 kDa, with a polypeptide portion of ~10.2 kDa and a carbohydrate portion of ~4.7 kDa. The α subunit is N-glycosylated at Asn⁵² and Asn⁷⁸.⁷

The β subunit has a molecular mass of 23 kDa, with a polypeptide portion of ~16.0 kDa and a carbohydrate portion of ~7.0 kDa is for the carbohydrate.^{4,6,9,10} The β subunit is N-glycosylated at Asn¹³ and Asn³⁰.¹¹ The structure and location of the O-glycosidic carbohydrate units attached to the β subunit of hCG, at Ser¹²¹, Ser¹²⁷, Ser¹³², and Ser¹³⁸, have been investigated.¹² The β subunit has been characterized by LC-MS.¹³ A MALDI mass spectrometry report on the glycoforms of the hCG β -core fragment has been published.¹⁴

When hCG was used in combination with recombinant interferon- γ , there was a significant cooperative induction of nitric oxide synthesis (iNOS) in a dose-dependent manner in mouse peritoneal macrophages. This suggests that hCG may provide a second signal for synergistic induction of NO synthesis.¹⁵ hCG has been used as a biomarker for such conditions as epithelial cancer¹⁸ and germ cell tumors.^{17,18}

Chorionic gonadotropin human (CG5) is 0.2 μ m-filtered. A vial contains ~5,000 IU. The calculated value is based on the USP method for human Chorionic Gonadotrophin to meet 5,000 IU per vial, with a tolerance of not less than 80.0% (4,000 IU) and not more than 125.0% (6,250 IU) of the potency stated.

Preparation Instructions

When the contents of a vial of CG5 are reconstituted with water to an hCG concentration of 1,000 I.U./mL, the hCG solution will also contain 0.02 M sodium phosphate buffer, pH ~7.2, and mannitol (stabilizer) at a concentration of 20 mg/mL. hCG is also soluble in aqueous glycerol and glycols and is insoluble in ethanol.³ Solutions should be sterile-filtered and not autoclaved.

Precautions and Disclaimer

This product is for R&D use only, and not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

References

1. MORGANS, F., Canfield, R. E., Vaitukaitis, J. L., & Ross, G. T. (1974). Properties of the subunits of human chorionic gonadotropin. *Endocrinology*, 94(6), 1601-1606.
2. Hogan, B., et al., Manipulating the Mouse Embryo, A Laboratory Manual, 2nd ed., Cold Spring Harbor Laboratory Press (Cold Spring Harbor, NY), p. 130 (1994).
3. Bellisario, R., Carlsen, R. B., & Bahl, O. P. (1973). Human chorionic gonadotropin: linear amino acid sequence of the α subunit. *Journal of biological chemistry*, 248(19), 6796-6809.
4. Morgan, F. J., Birken, S., & Canfield, R. E. (1975). The amino acid sequence of human chorionic gonadotropin. The alpha subunit and beta subunit. *Journal of Biological Chemistry*, 250(13), 5247-5258.
5. Carlsen, R. B., Bahl, O. P., & Swaminathan, N. (1973). Human chorionic gonadotropin: linear amino acid sequence of the β subunit. *Journal of biological chemistry*, 248(19), 6810-6827.
6. Lapthorn, A., Harris, D. C., Littlejohn, A., Lustbader, J. W., Canfield, R. E., Machin, K. J., ... & Isaacs, N. W. (1994). Crystal structure of human chorionic gonadotropin. *Nature*, 369(6480), 455-461.
7. de Beer, T., van Zuylen, C. W., Hård, K., Boelens, R., Kaptein, R., Kamerling, J. P., & Vliegthart, J. F. (1994). Rapid and simple approach for the NMR resonance assignment of the carbohydrate chains of an intact glycoprotein Application of gradient-enhanced natural abundance 1H - 13C HSQC and HSQC-TOCSY to the α -subunit of human chorionic gonadotropin. *FEBS letters*, 348(1), 1-6.
8. De Beer, T., Van Zuylen, C. W., Leeftang, B. R., Hård, K., Boelens, R., Kaptein, R., ... & Vliegthart, J. F. (1996). NMR Studies of the Free α Subunit of Human Chorionic Gonadotropin: Structural Influences of N-Glycosylation and the β Subunit on the Conformation of the α Subunit. *European journal of biochemistry*, 241(1), 229-242.
9. Bahl, O. P. (1969). Human chorionic gonadotropin: I. Purification and physicochemical properties. *Journal of Biological Chemistry*, 244(3), 567-574.
10. Swaminathan, N., & Bahl, O. P. (1970). Dissociation and recombination of the subunits of human chorionic gonadotropin. *Biochemical and Biophysical Research Communications*, 40(2), 422-427.
11. Sisinni, L., & Landriscina, M. (2015). The role of human chorionic gonadotropin as tumor marker: biochemical and clinical aspects. *Advances in Cancer Biomarkers: From biochemistry to clinic for a critical revision*, 159-176.
12. Kessler, M. J., Mise, T., Ghai, R. D., & Bahl, O. P. (1979). Structure and location of the O-glycosidic carbohydrate units of human chorionic gonadotropin. *The Journal of biological chemistry*, 254(16), 7909-7914.
13. Black, R. S., & Bowers, L. D. (2000). Mass Spectrometric Characterization of the β -Subunit of Human Chorionic Gonadotropin. *Mass Spectrometry of Proteins and Peptides: Mass Spectrometry of Proteins and Peptides*, 337-354.
14. Jacoby, E. S., Kicman, A. T., Laidler, P., & Iles, R. K. (2000). Determination of the glycoforms of human chorionic gonadotropin β -core fragment by matrix-assisted laser desorption/ionization time-of-flight mass spectrometry. *Clinical Chemistry*, 46(11), 1796-1803.
15. Kim, H. M., & Moon, Y. H. (1996). Human chorionic gonadotropin induces nitric oxide synthase mRNA in mouse peritoneal macrophages. *Biochemical and biophysical research communications*, 229(2), 548-552.
16. Iles, R. K. (2007). Ectopic hCG β expression by epithelial cancer: malignant behaviour, metastasis and inhibition of tumor cell apoptosis. *Molecular and cellular endocrinology*, 260, 264-270.
17. Mitsios, J. V., McClellan, A., Brown, S., & Gronowski, A. M. (2014). Human chorionic gonadotropin and α -fetoprotein in cerebral spinal fluid: method validation and retrospective review. *Clinical Biochemistry*, 47(7-8), 632-635.
18. Barlow, L. J., Badalato, G. M., & McKiernan, J. M. (2010). Serum tumor markers in the evaluation of male germ cell tumors. *Nature Reviews Urology*, 7(11), 610-617.

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2016-2025 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

23252148 Rev 08/25

**Millipore
Sigma**