




VHP Compatible Milliflex[®] PLUS Pump Hub User Guide

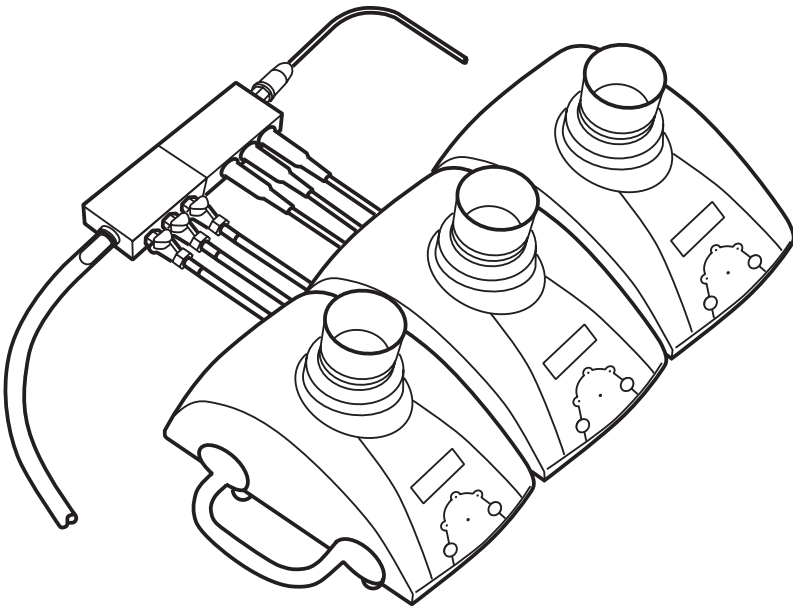


Introduction


The VHP Compatible Milliflex PLUS Pump Hub is used to connect up to three Milliflex PLUS pumps.

Fluid Line Connections


Push the Milliflex PLUS Pump fluid line connectors into the fluid hub until a click is heard. Install the other end of the tubing onto the pumps.

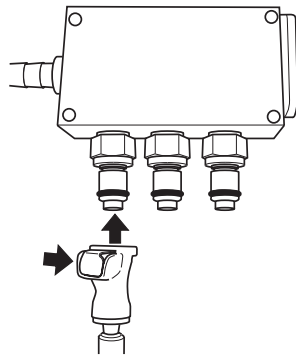

Insert the output tubing onto the stainless steel connector.

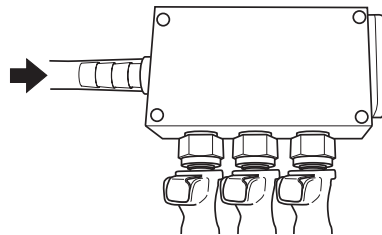


To disconnect the fluid lines, push on the button and pull the connector.

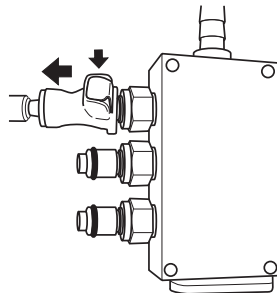


Electrical Connection


Push the Milliflex PLUS Pump electrical cables into the electrical hub until a click is heard. Connect the other end of the cables onto the pumps.

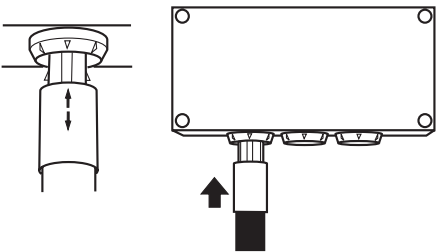

Connect the power supply to the electrical hub and tighten the connector.

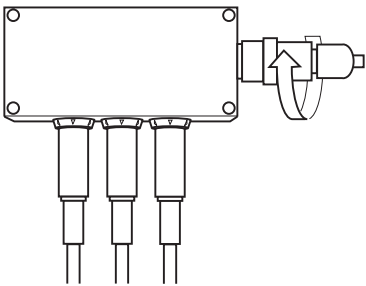


To disconnect the electrical cables from the hub, pull the cables out of the hub.

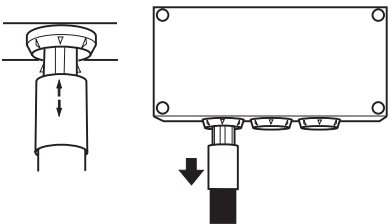

Connecting the Hubs


The fluid and electrical hubs may be connected or disconnected by sliding them together.

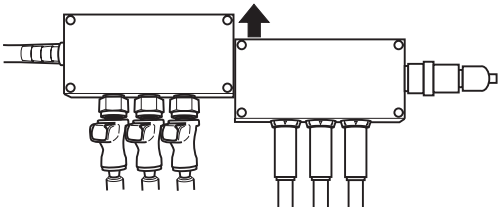

Decontamination

The hubs must be cleaned after each use. To decontaminate the fluid line hub and the tubing, use the sanitizing mode of the Milliflex PLUS Pump (refer to the Milliflex PLUS Pump User Guide). The external parts of both hubs may be sprayed with an alcohol solution or a standard laboratory surface decontamination solution.

Hang the hubs inside the isolator during the decontamination cycle as shown below.

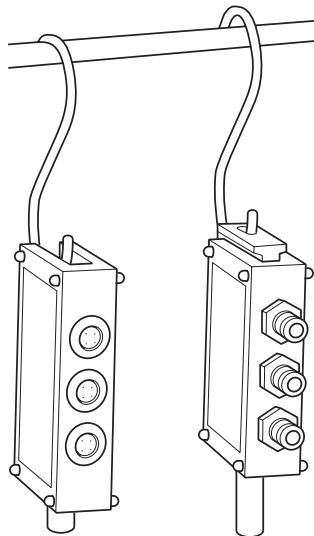

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

For more information:

Visit <http://www.millipore.com/techservice> or, in the U.S., call 1-800-MILLIPORE (1-800-645-5476).



Milliflex® is a registered trademark of Merck KGaA, Darmstadt, Germany.

The M mark is a trademark of Merck KGaA, Darmstadt, Germany.

© Copyright 2013 by Millipore SAS, Molsheim, France.
All rights reserved.