# MILLIPLEX® MAP Human Myokine Panel is an optimized, quantitative immunoassay that simultaneously measures 15 novel muscle-secreted factors

Skeletal muscle is actively involved in the synthesis and secretion of many proteins collectively termed "myokines." Myokines can act in autocrine and/or paracrine manners to regulate skeletal muscle metabolism and muscle proliferation and differentiation (muscle growth), as shown in Figure 1.

In addition, these myokines can act as endocrine hormones in altering cell metabolism, endothelial function and tumor growth/retardation. These myokines are also involved in inflammatory responses in a wide variety of tissues, including heart, adipose tissue, breast, colon, liver, intestine, endothelium, bone and skeletal muscle.

Emerging evidence supports the fact that the anti-inflammatory, disease-fighting effects of exercise may be mediated by these myokines, and that myokine signaling pathways intersect with innate immunity, neurological signaling and insulin response<sup>2,3</sup>.

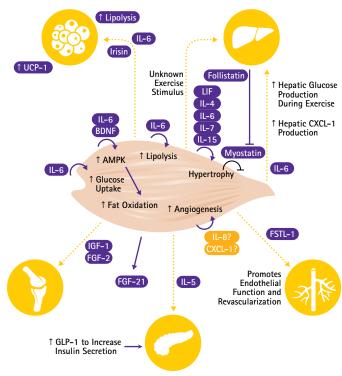


Figure 1.

Myokines, secreted by skeletal muscle, signal to diverse organs, including the liver, pancreas, bone and circulatory system. The precise mechanisms by which angiogenesis is upregulated are still being studied. Adapted from Pedersen BK, Febbraio MA. Muscles, exercise and obesity: skeletal muscle as a secretory organ. Nat Rev Endocrinol. 2012 Apr 3; 8(8): 457 – 65.



To meet the increasing need to quantify myokines in preclinical and translational research models, MilliporeSigma has developed the MILLIPLEX® MAP Human Myokine Panel. This 15-plex kit is designed for the simultaneous quantification of any or all of the following analytes in serum or plasma samples, which have the indicated biological functions:

- APLN: Apelin The function of apelin depends on the tissue in which it is being expressed. Apelin regulates blood pressure in vascular tissue, heart contractions in cardiac tissue, food and water intake in the brain, and glucose uptake and insulin inhibition in the digestive tract.
- BDNF: Brain-derived neurotrophic factor Associated with long-term memory, BDNF secretion increases in response to exercise.
- EPO: Erythropoeitin Because it stimulates red blood cell production, it may be administered to athletes to increase oxygen supply to the body, ostensibly improving performance.
- FABP3: Fatty acid-binding protein 3 This myokine is produced by heart muscle tissue and regulates cardiac uptake of long-chain fatty acids. It is a reliable and early biomarker of myocardial infarction and acute coronary events.
- **FGF21:** Fibroblast growth factor 21 FGF21 is activated by Akt and is a metabolism-regulating hormone and myokine.
- Irisin: The cleaved and secreted portion of fibronectin type III domain-containing protein 5 (FNDC5) The once-debated function of irisin has now been settled by a recent proteomics study using tandem mass spectrometry<sup>4</sup>. Irisin levels increase in response to exercise and may mediate the browning of white fat.
- FSTL1: Follistatin-related protein 1 Regulated by nitric oxide signaling, FSTL1 induces growth of new blood vessels in muscle and protects cardiac myocyte from ischemic-induced apoptosis.
- CX3CL1: Fractalkine CX3CL1 is a chemokine protein expressed on monocytes, natural killer cells, T cells, and smooth muscle cells. It has recently been associated with obesity, insulin resistance and Type 2 Diabetes.
- IL-6: Interleukin 6 IL-6 is, in a way, a "master myokine," in that it regulates multiple, divergent processes in response to exercise, ranging from glucose production, fat oxidation and lipolysis.
  IL-6 also regulates the immune response to exercise and triggers anti-inflammatory signaling.
- IL-15: Interleukin 15 IL-1 5 regulates T and NK cell activation and proliferation. As a myokine, it significantly reduces visceral fat.
- LIF: Leukemia inhibitory factor This protein inhibits differentiation and may activate cardiac and muscle stem cells in response to exercise.
- GDF8: Growth differentiation factor 8 (Myostatin) –
   In a feedback loop, myostatin inhibits the growth of
   muscle cells and blocks the differentiation of stem
   cells toward muscle lineages.
- OSM: Oncostatin M Oncostatin is in the IL-6 family of cytokines and likely shares functions of IL-6, particularly in the regulation of inflammation.

- **OSTN:** Osteocrin (Musclin) The main role of musclin is in the regulation of glucose metabolism.
- SPARC: Secreted protein acidic and rich in cysteine (Osteonectin) – SPARC is a myokine that blocks tumor progression in the colon by inducing apoptosis, in response to exercise.

# **Materials and Methods**

Multiplex assays using the standards included in the assay kit, as well as serum and plasma samples, were conducted using the MILLIPLEX® MAP Human Myokine Magnetic Bead Panel (Cat. No. HMYOMAG-56K), following the instructions in the included protocol. Capture antibody-coated beads were incubated with samples in order to immobilize the analyte. The biotinylated detection antibody was added; this detection antibody binds to the analyte to form a sandwich. Streptavidin-phycoerythrin (SA-PE) dye was added so it could bind to the sandwich and emit fluorescent light.

#### Results

Using standards for each analyte serially diluted in serum matrix, standard curves were prepared to determine the assay response (measured in Mean Fluorescence Intensity [MFI] with respect to analyte concentration). The standard curves established that the assay generated linear MFI response over 3 to 5 orders of magnitude for all analytes (Figure 2). No significant cross-reactivity was seen within the panel (Table 1).

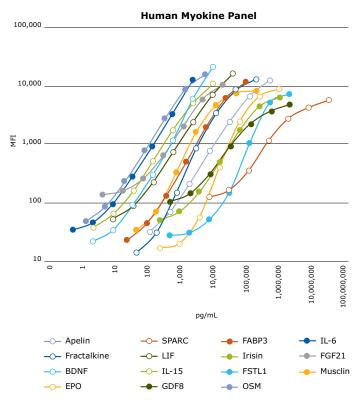


Figure 2.

Using standards for each analyte serially diluted in the validated serum matrix, standard curves were prepared to determine the assay response with respect to analyte concentration. The standard curves established linearity of the assay over three to four orders of magnitude for all analytes. The concentrations of analyte were measured in Mean Fluorescence Intensity (MFI).

# **Cross-Reactivity of Myokine Multiplexed Assays**

Fig.		MFI	Apelin	Fractalkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
Part	Frac		21	1108	17	15	109	43	27	91	15	45	25	32	29	122	27
Part		4	20	1115	16	16	113	43	27	100	15	44	26	35	30	133	27
Fig.			20	3985	17	16	113	43	30	95	14	44	27	33	31	135	26
Part		5	20	4039	17	15	116	43	30	97	15	44	26	34	30	135	27
Part			22	8821	18	17	115	46	29	91	17	47	24	36	30	130	27
BANK   Mart			20	8775	16	17	119	45	29	94	15	45	27	34	30	132	25
Part			21	13807	18	16	121	47	30	97	17	47	28	36	30	133	27
Fame			23		18	16	120	44	30	96	17	49	28	37	32	136	
Fig.   16	BDNF													-			
Second   S																	
Section   Sect																	
Fig.																	
Std																	
Pro																	
Second Part																	
Mathematical Part	FPO	std															
Std   19	2.0																
S		std										-				-	
Std   19																	
Secondary   Seco		std	19	12	23	7275	121	47	31	102	16	50	26	32	31	139	26
SPARC   Std   18		6	20	12	20	7080	120	48	28	99	17	48	26	33	30	136	28
SPARC   18			21	12	43	8686	117	43	27	97	15	44	26	33	29	127	27
Heat		7	19	12	37	8809	118	44	29	92	15	46	25	33	29	130	27
State   Stat	SPARC		18	12	17	16	1377	41	26	92	14	44	26	31	28	123	26
Start   Star		4	19	11	16	15	1359	42	29	94	16	47	22	31	28	133	25
Second   S			21	12	16	15	3061	41	29	91	14	44	24	31	29	127	26
Fig.																	
The color of the																	
Total Part																	-
Start   Star																-	
Part		-										-				-	
Std   20	LIF															-	
Fig.		std															
std     22     11     16     16     114     9555     27     90     14     42     24     32     28     117     25       std     22     12     15     14     116     9624     28     94     14     46     24     33     28     127     25       std     24     12     18     16     125     18040     28     89     15     44     25     36     28     128     25       7     26     12     19     17     116     18661     27     92     15     47     25     35     30     134     27       IL-15     44     19     10     14     15     102     38     590     90     14     43     25     30     29     124     26       44     19     11     16     16     112     41     572     101     15     44     24     31     31 <th< th=""><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th><th></th></th<>																	
6     22     12     15     14     116     9624     28     94     14     46     24     33     28     127     25       std     24     12     18     16     125     18040     28     89     15     44     25     36     28     128     25       7     26     12     19     17     116     18661     27     92     15     47     25     35     30     134     27       II-15     std     19     10     14     15     102     38     590     90     14     43     25     30     29     124     26       4     19     11     16     16     112     41     572     101     15     44     24     31     31     132     25       5td     18     12     15     16     112     41     2036     100     14     45     25     31     28		std															
The following color   The first color   The fi										94							
IL-15		std	24	12	18	16	125	18040	28	89	15	44	25	36	28	128	25
A     19     11     16     16     112     41     572     101     15     44     24     31     31     132     25       std 5     18     12     15     16     112     41     2036     100     14     45     25     31     28     133     25       std 5     19     11     15     16     108     40     1970     94     14     44     22     30     28     134     26       std 6     20     11     16     15     110     42     4888     95     15     43     25     31     29     134     25       std 7     19     12     16     15     113     51     10834     98     15     45     25     31     28     136     25       For 7     20     12     16     16     108     51     10730     97     14     45     24     34     30     135		7	26	12	19	17	116	18661	27	92	15	47	25	35	30	134	27
Std     18     12     15     16     112     41     2036     100     14     45     25     31     28     133     25       5     19     11     15     16     108     40     1970     94     14     44     22     30     28     134     26       std     20     11     16     15     110     42     4888     95     15     43     25     31     29     134     25       6     18     11     16     17     107     42     5039     94     13     43     24     31     29     133     25       std     19     12     16     15     113     51     10834     98     15     45     25     31     28     136     25       GDF8     std     19     11     14     15     100     39     26     425     14     43     24     30     27	IL-15		19	10	14	15	102	38	590	90	14	43	25	30	29	124	26
5     19     11     15     16     108     40     1970     94     14     44     22     30     28     134     26       std 6     20     11     16     15     110     42     4888     95     15     43     25     31     29     134     25       std 7     19     12     16     15     113     51     10834     98     15     45     25     31     29     133     25       GDF8     std 19     12     16     16     108     51     10730     97     14     45     24     34     30     135     25       GDF8     std 19     11     14     15     100     39     26     425     14     43     24     30     27     115     25       std 20     19     11     15     15     106     38     25     421     13     46     25     31		4 	19	11	16	16	112	41	572	101	15	44	24	31	31	132	25
std     20     11     16     108     40     1970     94     14     44     22     30     28     134     26       std     20     11     16     15     110     42     4888     95     15     43     25     31     29     134     25       std     19     12     16     15     113     51     10834     98     15     45     25     31     29     133     25       GDF8     std     19     12     16     16     108     51     10730     97     14     45     24     34     30     135     25       GDF8     std     19     11     14     15     100     39     26     425     14     43     24     30     27     115     25       std     19     11     15     15     106     38     25     421     13     46     25 <td< th=""><th></th><th></th><th>18</th><th>12</th><th>15</th><th>16</th><th>112</th><th>41</th><th>2036</th><th>100</th><th>14</th><th>45</th><th>25</th><th>31</th><th>28</th><th>133</th><th>25</th></td<>			18	12	15	16	112	41	2036	100	14	45	25	31	28	133	25
6     18     11     16     17     107     42     5039     94     13     43     24     31     29     133     25       std     19     12     16     15     113     51     10834     98     15     45     25     31     28     136     25       GDF8     std     19     11     14     15     100     39     26     425     14     43     24     30     27     115     25       std     19     11     14     15     106     38     25     421     13     46     25     31     30     125     25       std     19     11     15     15     106     38     25     421     13     46     25     31     30     125     25       std     19     11     15     15     106     39     25     1269     13     45     23     31     28																	
18     11     16     17     107     42     5039     94     13     43     24     31     29     133     25       std 7     19     12     16     15     113     51     10834     98     15     45     25     31     28     136     25       GDF8     std 19     11     14     15     100     39     26     425     14     43     24     30     27     115     25       std 4     19     11     14     15     106     38     25     421     13     46     25     31     30     125     25       std 19     11     15     15     106     39     25     1269     13     45     23     31     28     127     24       std 5     19     11     15     15     104     39     25     1199     13     48     25     31     27     120					-												
Formula     7     20     12     16     16     108     51     10730     97     14     45     24     34     30     135     25       GDF8     std 4     19     11     14     15     100     39     26     425     14     43     24     30     27     115     25       std 5     19     12     17     15     106     38     25     421     13     46     25     31     30     125     25       std 5     19     11     15     15     106     39     25     1269     13     45     23     31     28     127     24       std 5     18     11     15     15     104     39     25     1199     13     48     25     31     27     120     24       std 20     12     15     14     102     38     27     2605     14     60     23     31     30 </th <th></th>																	
GDF8					-												
4 19 12 17 15 106 38 25 421 13 46 25 31 30 125 25   std 19 11 15 15 106 39 25 1269 13 45 23 31 28 127 24   5 18 11 15 15 104 39 25 1199 13 48 25 31 27 120 24   std 20 12 15 14 102 38 27 2605 14 60 23 31 30 124 25   6 19 11 15 16 107 38 29 2693 14 63 23 31 29 126 25   std 20 11 17 16 114 41 25 3847 14 126 23 32 29 123 25	- GDE8	etd															
std     19     11     15     15     106     39     25     1269     13     45     23     31     28     127     24       5     18     11     15     15     104     39     25     1199     13     48     25     31     27     120     24       std     20     12     15     14     102     38     27     2605     14     60     23     31     30     124     25       6     19     11     15     16     107     38     29     2693     14     63     23     31     29     126     25       std     20     11     17     16     114     41     25     3847     14     126     23     32     29     123     25	3DF0															-	
5 18 11 15 15 104 39 25 1199 13 48 25 31 27 120 24   std 20 12 15 14 102 38 27 2605 14 60 23 31 30 124 25   6 19 11 15 16 107 38 29 2693 14 63 23 31 29 126 25   std 20 11 17 16 114 41 25 3847 14 126 23 32 29 123 25		std															
std     20     12     15     14     102     38     27     2605     14     60     23     31     30     124     25       6     19     11     15     16     107     38     29     2693     14     63     23     31     29     126     25       std     20     11     17     16     114     41     25     3847     14     126     23     32     29     123     25																	
std 20 11 17 16 114 41 25 3847 14 126 23 32 29 123 25		std															
7		6	19	11	15	16	107	38	29	2693	14	63	23	31	29	126	25
18 12 18 15 105 41 27 <mark>3897</mark> 16 125 23 30 28 125 26			20	11	17	16	114	41	25	3847	14	126	23	32	29	123	25
		/	18	12	18	15	105	41	27	3897	16	125	23	30	28	125	26

Table 1.

Minor cross-reactivity (<5%) was observed between the musclin standard and all other analytes' antibodies. In addition, the musclin sample values fall in the low end of standard curve. Considering the musclin sample values fall in the low end of the standard curve, this amount of cross-reactivity would not be expected to affect the assay accuracy.

# **Cross-Reactivity of Myokine Multiplexed Assays (continued)**

	MFI	Apelin	Fractalkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
FABP3	std	18	13	19	16	92	48	32	105	568	49	28	34	34	150	28
	4	22	17	20	20	97	51	31	107	591	46	28	32	31	147	28
	std	20	14	19	17	98	49	32	104	2204	48	29	32	31	144	28
	5	22	15	21	17	103	47	33	110	2337	50	29	32	33	150	26
	std	21	15	20	17	104	52	32	110	6903	50	31	36	34	150	32
	6	19	16	20	18	101	52	35	112	6886	53	31	34	35	153	29
	std 7	23	14	20	21	108	51	30	113	12255	51	31	36	34	156	28
		20	15	20	18	104	51	28	108	12404	52	29	33	31	146	28
Irisin	std 4	26	20	22	22	108	56	36	110	22	709	32	37	39	159	36
		30	19	24	22	104	52	39	111	20	733	39	40	46	163	37
	std 5	25	16	23	27	101	56	37	124	22	2360	38	41	43	161	35
		28	19	26	22	114	56	34	123	24	2339	33	40	40	165	33
	std 6	27	16	24	20	111	56	35	112	22	5265	35	40	38	150	34
		25	18	27	21	115	55	36	114	23	5304	35	39	39	166	35
	std 7		18 15	24	22 24	113	55 54	38	107	25 21	6802	35	40 37	40	161 159	35 36
FSTL1	std	27	19	24	22	115	54	34	111	19	6814 56	36 226	36	40	146	30
13111	4	26	16	27	20	111	59	37	117	19	58	238	37	37	169	33
	std	25	18	25	23	118	51	33	116	19	63	1384	37	34	163	35
	5	26	18	25	23	112	54	34	111	21	59	1377	38	39	157	36
	std	26	16	22	20	104	58	37	112	21	67	5904	38	42	155	35
	6	26	17	24	21	116	51	37	111	22	70	6156	38	44	166	34
	std	26	18	26	20	115	58	35	111	24	105	8090	39	37	174	33
	7	26	21	26	22	124	55	33	110	26	105	8025	40	38	177	36
OSM	std	23	20	22	21	111	56	36	106	22	54	31	1201	39	143	33
	4	20	18	24	22	107	53	35	108	23	55	32	1192	39	161	31
	std	24	17	24	21	106	54	40	103	21	54	31	3857	40	154	33
	5	25	19	24	22	110	53	37	112	24	54	32	3711	39	155	33
	std	25	18	21	23	106	51	34	103	23	56	31	10305	41	156	34
	6	27	19	23	21	111	55	35	113	26	58	30	10428	40	159	33
	std 7	26	19	23	24	122	53	39	104	23	54	31	17246	39	154	30
		26	19	22	25	105	55	36	120	22	57	35	17557	43	160	34
IL-6	std 4	31	17	20	23	101	52	36	121	21	56	35	40	403	169	33
		26	18	22	20	110	55	32	115	22	52	35	38	381	166	35
	std 5	25	16	25	22	108	52	38	128	23	57	32	40	1314	175	33
		26	18	24	23	112	55	38	130	22	57	31	44	1325	178	36
	std 6	26	16	20	21	107	50	38	108	24	52	31	40	4245	155	33
		22	21	25	23	117	53	33	118	21	56	35	41	4445	159	36
	std 7	26	17	24	24	113	51	41	105	24	52	34	43	15480	161	32
FGF21	std	24 25	17 17	26 22	24 23	114 107	53 52	38	113	22	55 52	32 32	44 39	15857 38	161 1298	33 34
10121	4	24	18	22	21	107	51	35	120	24	56	32	35	37	1251	33
	std	23	16	23	25	111	51	37	117	20	59	31	42	38	4078	35
	5	26	18	23	23	108	53	37	114	20	55	32	39	40	3879	35
	std	26	23	21	26	111	52	35	114	23	63	32	46	40	9184	37
	6	24	21	25	26	108	55	36	116	25	63	33	40	41	8977	40
	std	37	37	34	32	120	58	39	130	30	97	45	64	41	13589	52
	7	34	37	33	29	126	63	41	123	29	102	40	68	40	13740	56
Musclin	std	34	18	24	24	108	63	34	112	27	58	35	48	42	157	2097
	4	30	20	22	26	110	67	40	113	25	56	39	43	36	161	2056
	std	53	27	42	42	129	127	42	146	42	83	63	78	41	196	5795
	5	52	34	50	42	134	125	43	145	40	85	56	85	45	207	5751
	std	180	97	200	138	278	392	75	286	154	276	184	306	64	352	8683
	6	193	92	206	147	283	407	80	300	155	291	211	316	73	374	8251
	std	676	353	719	514	801	1332	212	810	577	923	712	1176	138	1001	9107
	7	754	393	824	552	880	1487	235	892	605	1036	791	1223	158	1082	9132
	%	1.4	0.7	3.0	1.6	0.0	2.8	0.5	1.1	1.5	2.6	4.4	1.9	0.5	2.2	

# Table 1 (continued).

Minor cross-reactivity (<5%) was observed between the musclin standard and all other analytes' antibodies. In addition, the musclin sample values fall in the low end of standard curve. Considering the musclin sample values fall in the low end of the standard curve, this amount of cross-reactivity would not be expected to affect the assay accuracy.

Most of the antibodies in this human myokine panel can recognize monkey samples, except Fractalkine, LIF and IL-6 (Table 2). Apelin cross-reacts with mouse, dog, rabbit and minipig. BDNF cross-reacts with mouse and rabbit. EPO cross-reacts with dog and rabbit. IL-15 cross-reacts with horse. Myostatin cross-reacts with dog, minipig and horse. FABP3 cross-reacts with rat,

rabbit, minipig and horse. FSTL1 cross-reacts with mouse, rabbit, minipig and horse. OSM cross-reacts with dog and minipig. FGF21 cross-reacts with dog, minipig and horse. Note that only limited amount of normal samples were tested in this study, and we could not rule out the possible cross-reactivity in other normal or disease samples.

# Myokine Concentration [pg/mL]

	Apelin	Fractalkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
Human Exercise Sample 1	198	396	10844	1847	307150	48	28		40040	10652	16352	126	6143	5996	90
Human Exercise Sample 2	45	35	8238		289560				12736	2913	11651	16	187	2657	76
Human Exercise Sample 3	64	192	95	6779	173870		11		1258	10292	6379		91		104
Human Exercise Sample 4	254	137	15964	1152	271440			18312	6835	1539	9705	9	10	434	157
Monkey 1	122		172	8475	262260		14	1136	10811		2100	7		404	246
Monkey 2	170		3180	2006	172540				9451	768				145	2678
Monkey 3	163		2335	28661	359810		25		6345		3308	14		588	1630
Monkey 4	290		1061	7003	210100		20		10343		6088			367	75
Rat 1			4800						2477						
Rat 2			269						1971					152	
Rat 3			627						2948						
Rat 4			4725						2732						
Mouse 1			11						8						
Mouse 2	3715		208			157			50		7102				
Mouse 3	1615		82			61			36		8665				
Mouse 4									7	6289					
Dog 1				862				11409	163			36		449	
Dog 2	1244		100		7110				36					239	
Dog 3				443				2890	6			31		171	
Dog 4								12413	8	768		33		394	
Rabbit 1	876		44		6110				317		17725				
Rabbit 2	672		36						232		12108				
Rabbit 3	906		33	416					229		14563				
Rabbit 4	918		37						206		15646				
Minpig 1								9578	8	1539	21471			95	
Minpig 2								9356	8	706	44759	21		104	
Minpig 3	118			735				4442	102			19			
Minpig 4	696	797		1290	22700	94	17	478343	529	1918	18559	345		2066	
Horse 1							19	10505	451		6814			115	
Horse 2							33	7474	218		7747		11		
Horse 3							22	11606	491		7960			134	
Horse 4							18	1043	165		3598				

**Table 2.**Samples, taken from various species, show not all myokines are expressed to the same degree in all subjects or tissues.

	Apelin	Frac- talkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
							pg/m	L							
std 1	122	49	2	244	7324	10	2	488	24	244	488	1	1	5	49
std 2	488	195	10	977	29297	39	10	1953	98	977	1953	6	2	20	195
std 3	1953	781	39	3906	117188	156	39	7813	391	3906	7813	23	9	78	781
std 4	7813	3125	156	15625	468750	625	156	31250	1563	15625	31250	94	38	313	3125
std 5	31250	12500	625	62500	1875000	2500	625	125000	6250	62500	125000	375	150	1250	12500
std 6	125000	50000	2500	250000	7500000	10000	2500	500000	25000	250000	500000	1500	600	5000	50000
std 7	500000	200000	10000	1000000	30000000	40000	10000	2000000	100000	1000000	2000000	6000	2400	20000	200000
							MFI								
std 1	33	14	22	17	124	54	38	104	23	51	28	48	35	137	34
std 2	69	31	35	21	163	89	65	143	44	71	31	88	46	161	71
std 3	212	150	94	58	354	227	157	310	133	155	53	243	95	258	333
std 4	750	849	307	411	1,154	736	524	906	495	493	155	806	282	629	1,625
std 5	2420	3328	1175	2431	2770	2401	1767	2169	1999	1717	1029	2839	930	2029	4693
std 6	6581	8392	5990	6602	4213	7084	4981	3603	6238	4441	5059	8730	3321	5997	7515
std 7	12355	12910	21111	8705	5622	16056	10651	4744	11610	6275	7170	15794	12158	10415	8098
Sensitivity	63.9	40.4	8.4	814.6	7481.1	4.1	4.9	258.6	9.7	281.2	736.7	5.7	0.9	14.1	59.2
Intra-Assay CV% QC (10 assays)	2.1	2.7	3.6	1.6	2.5	2.2	1.9	2.8	2.6	3.9	3.1	5.6	4.2	4.1	7.2
Inter-Assay CV% QC (10 assays)	10.7	7.9	14.2	5.9	15.3	6	9.4	7.7	9	8.8	7.6	9.5	13	10	11.9

**Table 3.**Assay precision, as determined by measuring signals from included standards.

Known amounts of each analyte (points 3, 4, 5 on the standard curves) were spiked into normal human serum/plasma samples. Percent recovery was calculated using the following equation: Calculated concentration of analyte / Expected concentration \*100 (Table 4).

# Spike and Recovery in serum [n=5] and plasma [n=4] samples

	Apelin	Fractalkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
							Average								
+std 3	100	96	105	71	90	95	109	95	97	105	119	84	100	82	106
+std 4	92	104	113	77	90	101	105	128	93	115	114	77	130	92	109
+std 5	85	94	135	72	75	101	99	150	90	99	97	72	155	90	156
Average	93	98	118	74	85	99	104	124	93	106	110	77	129	88	124

**Table 4.**Recovery of all spiked samples was between 70% and 130%.

Next, linearity tests were performed to determine the extent to which the dose-response of each analyte was linear in a particular diluent (Table 5). Calculations rely on dividing the Observed Values by the Expected Values. Unfortunately, some analytes were barely detectable in the human sepsis samples we tested. We were, therefore, compelled to spike known amounts

of protein standards into normal serum/plasma samples and test the dilution linearity. The linearity range was 70% to 130% (Table 4). The kit components were stable over a wide range of temperatures. Storing the kit at 37°C for 7 days did not significantly alter the assay performance (data not shown).

# Average Dilution Linearity in Human Serum/Plasma Samples [n=10 samples]

	Apelin	Fractalkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
1:4 dil			64	79	139		94	110	89		67	99	79	114	
1:8 dil			49	54	150		98	111	86		79	89	66	124	_
1:16 dil			93	79	120		118	89	95		109	84	98	93	
Average			69	71	136		103	103	90		85	90	81	110	

Table 5.

Average dilution linearity in human serum/plasma samples.

### Average Dilution Linearity in Spiked Serum/Plasma Samples [n=10 samples] (continued)

	Apelin	Fractalkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
1:4 dil	111	85	58	105	100	97	95	74	92	106	73	106	77	91	109
1:8 dil	123	84	48	108	98	91	94	63	87	105	70	110	66	84	111
1:16 dil	129	84	42	108	87	90	94	61	83	100	64	114	61	75	116
Average	121	84	49	107	95	93	94	66	87	104	69	110	68	83	112

#### Table 5 (continued).


Average dilution linearity in human serum/plasma samples.

To test the consistency and accuracy between MILLIPLEX® assays, we compared the quantitation of analytes existing in both this human myokine panel and other MILLIPLEX® panels. All the standard curves showed overlapping response and analyte concentrations calculated in biological samples correlated very well.

For example, FGF21 and FABP3 sample values were measured using both MILLIPLEX® Human Myokine Panel, Human Liver Protein Panel and Human CVD1 Panel. Standard curve created using standards of

purified FGF21, and FABP3 from two panels showed overlapping assay response and linear range of the assays. Assay correlation was excellent, with slope and R value approaching unity (Figure 3).

In addition to FGF21 and FABP3, we have compared the same analytes in this human myokine panel with other MILLIPLEX® panels. All the standard curves showed overlapping responses and analyte concentrations that correlated very well with biological samples (data not shown).

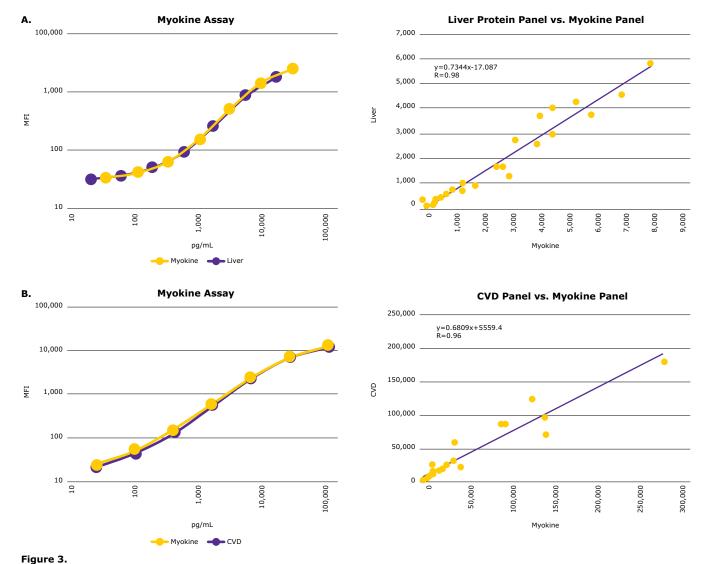

Comparison of the MILLIPLEX® MAP Human Myokine Panel with the Liver Panel and CVD Panel for the measurement of FGF21 (top) and FABP3 (bottom) respectively.

Finally, biological qualification of the assay panel was performed using serum/plasma samples from healthy subjects and from those with sepsis. As expected, 15 novel myokines were simultaneously quantified by the assay. FABP3, Irisin, FSTL1, OSM, IL-6, FGF21 and Musclin were significantly upregulated in sepsis samples (Table 6).

# Normal Sample Summary Data [pg/mL]

	Apelin	Fractalkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
N	8	9	10	11	12	13	14	15	16	17	18	19	20	21	22
Average	228	101	9548	1239	425343	10	3	8625	1943	495	6491	13	3	134	150
Min	0	0	350	0	169720	0	0	0	740	0	0	0	0	0	0
Max	420	260	20480	3460	739890	30	20	40110	3110	2370	17170	37	9	570	790
% detectable	88	88	100	63	100	50	13	38	100	38	88	75	88	63	50

# Sepsis Sample Summary Data [pg/mL]

	Apelin	Fractalkine	BDNF	EPO	SPARC	LIF	IL-15	Myostatin	FABP3	Irisin	FSTL1	OSM	IL-6	FGF21	Musclin
N	32	33	34	35	36	37	38	39	40	41	42	43	44	45	46
Average	122	78	9280	1298	544958	6	7	3777	38972	1905	14632	56	643	2016	174
Min	0	0	110	0	204800	0	0	0	240	0	0	0	0	0	0
Max	460	410	20180	12520	1258000	80	60	31600	249140	9060	74170	317	8679	12400	880
% detectable	69	69	100	41	100	19	25	28	100	60	84	97	81	75	59

#### Table 6.

Comparison of biomarkers expressed in normal and sepsis serum/plasma samples. FABP3, Irisin, FSTL1, OSM, IL-6, FGF21 and Musclin were shown to be upregulated in sepsis samples.

# Conclusion

The MILLIPLEX® MAP Human Myokine Panel is sensitive, accurate and reproducible. The sample values generated with this new panel are consistent with previous panels that include some of the same analytes. This panel provides an ideal immunoassay for diverse research areas, including metabolic, neuromuscular and idiopathic myopathy diseases.

# References

- Pedersen BK, Febbraio MA. Muscles, exercise and obesity: skeletal muscle as a secretory organ. Nat Rev Endocrinol. 2012 Apr 3; 8(8): 457 – 65.
- Egan B, Zierath JR. Exercise metabolism and the molecular regulation of skeletal muscle adaptation. Cell Metab. 2013 Feb 5; 17(2):162 – 84.
- 3. Pedersen BK, Hoffman-Goetz L. Exercise and the immune system: regulation, integration, and adaptation. Physiol Rev. 2000 Jul; 80(3): 1055 81.
- Jedrychowski MP, Wrann CD, Paulo JA, Gerber KK, Szpyt J, Robinson MM, Nair KS, Gygi SP, Spiegelman BM. Detection and Quantitation of Circulating Human Irisin by Tandem Mass Spectrometry. Cell Metab. 2015 Aug 12. pii: S1550-4131(15) 00392-7.

# To Place an Order or Receive Technical Assistance

In the U.S. and Canada, call toll-free 1(800)-645-5476

For other countries across Europe, call +44 (0) 115 943 0840

For other countries across Europe and the world, please visit **EMDMillipore.com/offices** 

For Technical Service, please visit **EMDMillipore.com/techservice** 

MilliporeSigma 400 Summit Drive Burlington, MA 01803

