



E.coli UDG

Product Number: UDGEC

Product Description

E.coli uracil-DNA glycosylase (UDG) is obtained from the expression and multi-step purification of the recombinant *E.coli* strain cloned with *E.coli* UDG gene. *E.coli* UDG catalyzes the release of uracil from uracil-containing DNA. UDG can efficiently hydrolyze uracil on single- or double-stranded DNA but cannot hydrolyze uracil from oligonucleotides of 6 bases or less.

Components

Components	UDGEC-500U	UDGEC-5000U
<i>E.coli</i> UDG (5 U/μl)	100 μl	1 mL

Storage Buffer

10 mM Tris-HCl, pH7.4@25°C
50mM KCl
0.1 mM EDTA
1 mM DTT
0.1mg/ml BSA
50% Glycerol (v/v)

Storage

Store at -30 ~ -15°C and transport at ≤ 0°C

Unit Definition

The amount of enzyme required to catalyze the release of 60 pmol of uracil from uracil-containing double-stranded DNA per minute is defined as 1 U. Activity was determined by measuring the amount of [³H]-uracil released from a 50 μl reaction system containing 0.2 μg DNA within 30 min at 37°C.

Notes

For research use only. Not for use in diagnostic procedures.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

To place an order or receive technical assistance

In Europe, please call Customer Service:

France: 0825 045 645

Germany: 069 86798021

Italy: 848 845 645

Spain: 901 516 645 Option 1

Switzerland: 0848 645 645

United Kingdom: 0870 900 4645

For other countries across Europe, please call: +44 (0) 115 943 0840

Merck KGaA
Frankfurter Strasse 250
64293 Darmstadt, Germany

For other countries across Europe and the world, please visit:
SigmaAldrich.com/offices

For Technical Service, please visit:
SigmaAldrich.com/techservice

SigmaAldrich.com

Experiment Process

Reaction System	
ddH ₂ O	To 50 μl
10 x Taq Buffer (with 20 mM MgCl ₂)	5 μl
dUTP ^a	0.6 mM
DATP/dCTP/dGTP	0.2 mM each
Template DNA	optional
Primer1 (10μM)	2 μl
Primer2 (10μM)	2 μl
Taq DNA Polymerase (5 U/μl)	0.5 μl
<i>E. coli</i> UDG (5 U/μl) ^b	0.2 μl

a) According to experimental needs, the final concentration of dUTP can be adjusted between 0.2 - 0.6mM.

b) According to experimental needs, the amount of 50 μl reaction system is generally 0.1 - 1 U.

▲ According to experimental needs, the final concentration of MgCl can be adjusted between 2 - 3 mM.

Reaction System		
37°C	10 min	Degradation of uracil-containing templates
95°C	2 min	UDG inactivation, template denaturation
PCR		
94°C	30 sec	} 30 - 35 cycles
55°C	30 sec	
72°C	60 sec/kb	
72°C	7 min	Fully extention

▲ According to the experimental needs, the PRC reaction program can be adjusted.