

68067 Candida Selective Supplement

An antibiotic supplement recommended for selective isolation of *Candida* species from mixed cultures.

Composition

(per vial, sufficient for 500 ml medium)

Gentamicin 50 mg

Directions

Rehydrate the contents of 1 vial with 5.0 ml sterile distilled water. Mix well and aseptically add to 500 ml sterile molten, cooled to 50°C *Candida* Ident Agar, modified (Cat. No. 94382). Mix well and pour into sterile petri plates.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

