

Data Sheet

## YUMMER1G.1F Mouse Cell Line

Cancer Cell Line

**SCC246****Pack Size  $\geq 1 \times 10^6$  viable cells/vial****Store at: Liquid nitrogen****FOR RESEARCH USE ONLY****Not for use in diagnostic procedures. Not for Human or Animal Consumption.**

### Background

The great promise of immune-based therapies in cancer and recent progress in successful application of these approaches has brought to the fore the necessity of immune-competent models to evaluate immune system responses to cancer cells. Melanomas exhibit relatively high somatic mutation burden, and these mutations may act as neoantigens that generate anti-tumor immune responses. The development of immunocompetent cell models is critical to the advancement of cancer immunotherapy and understanding of immunity, although few tractable model systems are available.

The Yale University Mouse Melanoma (YUMM) lines are a comprehensive system of melanoma cell lines harboring defined genetic drivers of melanoma that are syngeneic with the immunocompetent C57BL/6 mouse model. YUMM cell lines were derived from the Braf/Pten genetically engineered mouse melanoma model, a highly penetrant genetically-engineered mouse model of melanoma (GEMM) in which melanomas develop with short latency. YUMM Exposed to Radiation (YUMMER) cell lines are clonal isolates of YUMM lines exposed to UVB radiation to reflect the high somatic mutation burden in melanomas. YUMMER cells harbor a high frequency of stable somatic mutations compared to YUMM cells; these mutations have been shown to stimulate host adaptive immune response.<sup>3</sup> The YUMMER1G.1F mouse melanoma cell line is both immunocompetent and reflective of the somatic mutations common in melanomas. YUMMER1.G1F carries three driver mutations of melanoma: Braf V600E, Pten -/- and Cdkn2 -/-.<sup>1,2</sup> YUMMER1.G1F cells are diploid, allowing for enhanced knockout frequency in CRISPR-based screens, are of female genotype, and have pheomelanin-producing background.<sup>4</sup> The characteristics of the YUMMER1.G1F cell line make it a valuable model for investigating sex-related differences in immune checkpoint inhibition and mechanisms of anti-tumor responses.

### Source

The YUMMER1.G1F mouse melanoma cell line is a diploid clonal isolate from YUMMER.G cells exposed to UVB radiation.<sup>4</sup> The original YUMM1.1 cell line was derived from a 4-hydroxytamoxifen-induced melanoma tumor in a C57BL/6 mouse into which mutations from the Braf/Pten genetically-engineered mouse model had been introduced via backcrossing.<sup>1</sup>


## Quality Control Testing

- Each vial contains  $\geq 1 \times 10^6$  viable cells.
- Cells are tested negative for infectious diseases by a Mouse Essential CLEAR Panel by Charles River Animal Diagnostic Services.
- Cells are verified to be of mouse origin and negative for interspecies contamination from rat, human, Chinese hamster, Golden Syrian hamster, and non-human primate (NHP) as assessed by a Contamination Clear panel by Charles River Animal Diagnostic Services.
- Cells are negative for mycoplasma contamination.

## Storage and Handling

YUMMER1.G1F cells should be stored in liquid nitrogen. The cells can be cultured for at least 10 passages after initial thawing without significantly affecting functionality.

## Representative Data

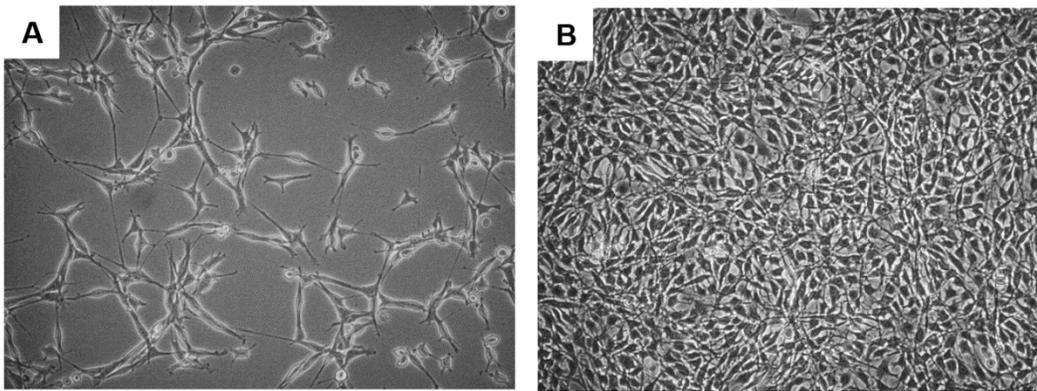

**Figure 1.** Brightfield images of YUMMER1.G1F one (A) and three (B) days after thaw in a T75 flask.

## Protocols

### Thawing Cells

1. Do not thaw the cells until the recommended medium is on hand. Cells can grow on normal tissue cultureware surfaces without any additional coating.  
Cells are thawed and expanded in YUMMER1.G1F Expansion Medium comprised of DMEM/F12 (Cat. No. DF-041-B) with 1X Non-Essential Amino Acids (Cat. No. TMS-001-C), 10% FBS (Cat. No. ES-009-B) and 1X penicillin/streptomycin (Cat. No. TMS-AB2-C, optional).
2. Remove the vial of frozen YUMMER1.G1F cells from liquid nitrogen and incubate in a 37 °C water bath. Closely monitor until the cells are completely thawed. Maximum cell viability is dependent on the rapid and complete thawing of frozen cells.  
**IMPORTANT:** Do not vortex the cells.
3. As soon as the cells are completely thawed, disinfect the outside of the vial with 70% ethanol. Proceed immediately to the next step.
4. In a laminar flow hood, use a 1- or 2-mL pipette to transfer the cells to a sterile 50 mL conical tube. Be careful not to introduce any bubbles during the transfer process.
5. Using a 10 mL pipette, slowly add dropwise 9 mL of YUMMER1.G1F Expansion Medium (Step 1 above) to the 15 mL conical tube.  
**IMPORTANT:** Do not add the entire volume of media all at once to the cells. This may result in decreased cell viability due to osmotic shock.

6. Gently mix the cell suspension by slowly pipetting up and down twice. Be careful not to introduce any bubbles.  
**IMPORTANT:** Do not vortex the cells.
7. Centrifuge the tube at  $300 \times g$  for 2-3 minutes to pellet the cells.
8. Decant as much of the supernatant as possible. Steps 5-8 are necessary to remove residual cryopreservative (DMSO).
9. Resuspend the cells in 15 mL of YUMMER1.G1F Expansion Medium.
10. Transfer the cell mixture to a T75 tissue culture flask.
11. Incubate the cells at 37 °C in a humidified incubator with 5% CO<sub>2</sub>.

### Subculturing Cells

1. YUMMER1.G1F cells should be passaged at ~80-85% confluency. Do not allow the cells to grow over 85% confluency.
2. Carefully remove the medium from the T75 tissue culture flask containing the 80% confluent layer of YUMMER1.G1F cells.
3. Rinse the flask with 10 mL 1X PBS. Aspirate after the rinse.
4. Apply 5-7 mL of Accutase™ and incubate in a 37 °C incubator for 3-5 minutes.
5. Inspect the flask and ensure the complete detachment of cells by gently tapping the side of the flask with the palm of your hand.
6. Add 10 mL of YUMMER1.G1F Expansion Medium to the plate.
7. Gently rotate the flask to mix the cell suspension. Transfer the dissociated cells to a 15 mL conical tube.
8. Centrifuge the tube at  $300 \times g$  for 3-5 minutes to pellet the cells.
9. Discard the supernatant, then loosen the cell pellet by tapping the tip of the tube with a finger.
10. Apply 2-5 mL of YUMMER1.G1F Expansion Medium to the conical tube and resuspend the cells thoroughly.  
**IMPORTANT:** Do not vortex the cells.
11. Count the number of cells using a hemocytometer.
12. Plate the cells to the desired density. Typical split ratio is 1:6.

### Cryopreservation of Cells

YUMMER1.G1F cells may be frozen in YUMMER1.G1F Expansion Medium and 10% DMSO using a Nalgene® slow freeze Mr. Frosty® container.

### References

1. Pigment Cell Melanoma Res 2016, 29(5): 590-597.
2. Nat Genet 2009, 41(5): 544-552.
3. Pigment Cell Melanoma Res 2017, 30(4): 428-435.
4. Ramseier JY, Charos A, Park K, Damsky W, Bosenberg MW (2019). Cancer Res 79(13 Suppl) Abstract #4622.

## Academic Use Agreement

Subject to local law

THIS PRODUCT MAY ONLY BE USED BY INDIVIDUALS EMPLOYED BY AN ACADEMIC INSTITUTION AND IS INTENDED SOLELY TO BE USED FOR ACADEMIC RESEARCH, WHICH IS FURTHER DEFINED BELOW. BY OPENING THIS PRODUCT, YOU ("PURCHASER") HEREBY REPRESENT THAT YOU HAVE THE RIGHT AND AUTHORITY TO LEGALLY BIND YOURSELF AND/OR YOUR EMPLOYER INSTITUTION, AS APPLICABLE, AND CONSENT TO BE LEGALLY BOUND BY THE TERMS OF THIS ACADEMIC USE AGREEMENT. IF YOU DO NOT AGREE TO COMPLY WITH THESE TERMS, YOU MAY NOT OPEN OR USE THE PRODUCT AND YOU MUST CALL MILLIPORESIGMA ("SELLER") CUSTOMER SERVICE (1-800-645-5476) TO ARRANGE TO RETURN THE PRODUCT FOR A REFUND.

"Product" means YUMMER1.G1F Mouse Cell Line (SCC246).

"Academic Research" means any internal in vitro research use by individuals employed by an academic institution. Academic Research specifically excludes the following uses of whatever kind or nature:

Re-engineering or copying the Product

Making derivatives, modifications, or functional equivalents of the Product

Obtaining patents or other intellectual property rights claiming use of the Product

Using the Product in the development, testing, or manufacture of a Commercial Product

Using the Product as a component of a Commercial Product

Reselling or licensing the Product

Using the Product in clinical or therapeutic applications including producing materials for clinical trials

Administering the Product to humans

Using the Product in collaboration with a commercial or non-academic entity

"Commercial Product" means any product intended for: (i) current or future sale; (ii) use in a fee-for-service; or (iii) any diagnostic, clinical, or therapeutic use.

Access to the Product is limited solely to those officers, employees, and students of PURCHASER's academic institution who need access to the Product to perform Academic Research. PURCHASER shall comply with all applicable laws in its use and handling of the Product and shall keep it under reasonably safe and secure conditions to prevent unauthorized use or access.

These use restrictions will remain in effect for as long as PURCHASER possesses the Product.

COMMERCIAL OR NON-ACADEMIC ENTITIES INTERESTED IN PURCHASING OR USING THE PRODUCT MUST CONTACT [licensing@emdmillipore.com](mailto:licensing@emdmillipore.com) AND AGREE TO SEPARATE TERMS OF USE PRIOR TO USE OR PURCHASE.

## Genetically Modified Organisms (GMO)

This product contains genetically modified organisms.

Este producto contiene organismos genéticamente modificados.

Questo prodotto contiene degli organismi geneticamente modificati.

Dieses Produkt enthält genetisch modifizierte Organismen.

Ce produit contient des organismes génétiquement modifiés.

Dit product bevat genetisch gewijzigde organismen.

Tämä tuote sisältää geneettisesti muutettuja organismeja.

Denna produkt innehåller genetiskt ändrade organismer.

---

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

## Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://SigmaAldrich.com/techservice).

## Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at [SigmaAldrich.com/terms](https://SigmaAldrich.com/terms).

## Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://SigmaAldrich.com/offices).

The life science business of Merck operates  
as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates.  
All other trademarks are the property of their respective owners. Detailed information on  
trademarks is available via publicly accessible resources.

© 2023-2024 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.  
Document Template 20306518 Ver 6.0  
20634623 Ver 3.0, Rev 10May2024, CJ

