



Rapid and Robust Protein Chromatography

Chromolith® WP 300 RP-18 2 mm I.D. HPLC Columns

The most hydrophobic of the Chromolith® WP 300 line, the RP-18 column is useful for the resolution of peptides and smaller proteins. One critical quality attribute (CQA) required by regulatory bodies is the peptide map of a biotherapeutic. Peptide maps generated by RP-HPLC provide valuable information about protein structure, stability, and purity. To be effective, the RP-HPLC column must be able to resolve a high percentage of the peptides in the sample. The more peptides, the better the information. The Chromolith® WP 300 RP-18 column gives unsurpassed RP-HPLC resolution of peptide maps from enzymatic digests. The improvements in silica and bonded-phase chemistry incorporated into the Chromolith® WP 300 line improve resolution by increasing efficiency and reducing peak tailing.

Key Benefits:


- Monolithic skeleton permits high flow rates to maximize throughput
- 300 Å mesopores permit large molecules to enter without fear of size-exclusion effects
- Matrix-loaded samples can be injected onto the column with little to no sample prep.

Protein Analysis on Chromolith® WP 300 RP-18

For large molecule separations, high efficiency separations are necessary in order to achieve resolution and good peak shape. Moving to sub-2 μm FPP-packed columns or 2.0 μm SPP-packed columns can deliver that desire; however, this comes at the cost of elevated backpressure. Chromolith® WP 300 RP-18, 2 mm I.D. columns provide UHPLC efficiencies, but at nearly 1/10th the backpressure.

Chromatographic conditions:

Columns:	Chromolith® WP 300 RP-18 100-2 mm (1.52370.0001) SPP, C18, 160 Å, 2.0 μm, 100-2.1 mm FPP, C18, 300 Å, 1.7 μm, 100-2.1 mm	
Mobile phase:	A: water (0.1% TFA) B: acetonitrile (0.08% TFA)	
Gradient:	4% B to 60% B in 10 minutes	
Flow rate:	0.38 mL/min	
Detection:	UV, 220 nm	
Column temperature:	30 °C	
Injection volume:	0.5 μL	
Sample:	HPLC Protein Mix 1 mg/mL, water 1) Ribonuclease 2) Cytochrome C 3) Holo-Transferrin 4) Apomyoglobin	

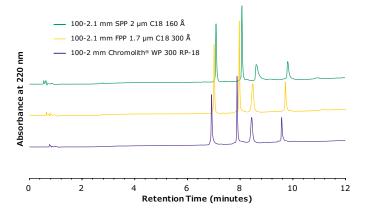



Excellent Lot-to-Lot Reproducibility

Chromolith® WP 300 RP-18 columns exhibit excellent batch-to-batch reproducibility, as demonstrated below with the same peptide map generated for Cytochrome C using three different Chromolith® WP 300 RP-18 columns across three different batches.

Chromatographic conditions:


Column	Chromolith® WP (1.52370.0001)	300 RP-18 100-	2 mm	
Mobile Phase	A: acetonitrile 0 B: water 0.1%	(, ,		
Gradient:	Time	%A	%B	
	0	5	95	
	25.0	30	70	
	30.0	30	70	
Flow rate:	0.190 mL/min			
Pressure:	18 bar			
Detection:	Vanquish DAD 20 Hz, UV, 214 nm			
Detector cell:	LightPipe 10 mm	1		
Temperature:	30 °C			
Injection volume:	0.2 μL			
Sample:	Rapid Trypsin Digestion with SOLu-Trypsin Rapid Digestion Kit 2.5 mg Cytochrome C was added in a PCR vial and dissolved in 320 μ L Rapid Trypsin Digestion Buffer. In the solution was 80 μ L SOLu-Trypsin added and incubated at 60 °C for 1 hour in a Thermomixer. The digestion was quenched by adding 12 μ L of hydrochloric acid 32 %.			


Antibody Fragment Analysis

Fragment analysis of a monoclonal antibody (mAb), also called middle-up analysis, is a useful technique in characterizing mAb domains without the inherent complexity of a peptide map. High efficiency is needed here to resolve subtle, structural variants of the mAb domains. The Chromolith® WP 300 RP-18 column is able to achieve the same separation efficiency and sensitivity as sub-2 μm FPP and 2.0 μm SPP-packed columns but at only 20% of the backpressure of those columns.

Chromatographic conditions:

Chromatographic conditions:					
Column	Chromolith® WP 300 RP-18, 2 mm I.D. (1.52370.0001) SPP, C18, 160 Å, 2.0 μ m, 100-2.1 mm FPP, C18, 300 Å, 1.7 μ m, 100-2.1 mm				
Mobile Phase	A: Water (0.1% (v/v) TFA)				
	B: Acetonitrile (0.08% (v/v) TFA)				
Gradient:	Time (Min)	%B			
	0	20			
	1	20			
	9	45			
Flow rate:	380 µL/min				
Detection:	UV, 220 nm				
Temperature:	80 °C				
Injection	1.0 µL				
volume:					
Sample:	SigmaMAb, 2 mg	/mL (SiLu™ Lite	Universal Antibody)		
DTT digest:	60 μL of 40 mM Dithiothreitol (DTT) solution was				
	added in a PCR vial, 40 µL mAb was added and				
	incubated at 37 °C for 30 minutes creating light chain				
	(LC) and heavy chain (HC) parts of the antibody.				

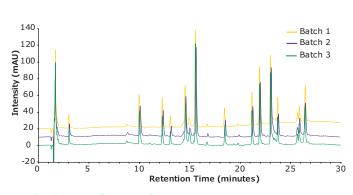

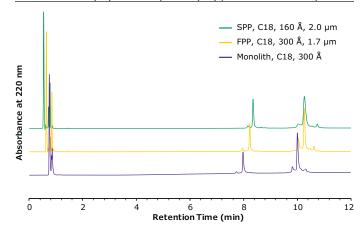

Ordering Information

Part Number	Description	Length (mm)	I.D. (mm)
1.52370.0001	Chromolith® WP 300 RP-18 Column	100	2
1.52371.0001	Chromolith® WP 300 RP-18 Column	50	2
1.52372.0001	Chromolith® WP 300 RP-18 Guard Columns (3 units)	5	2

To place an order or receive technical assistance

Order/Customer Service: SigmaAldrich.com/order Technical Service: SigmaAldrich.com/techservice

Safety-related Information: SigmaAldrich.com/safetycenter

MilliporeSigma 400 Summit Drive Burlington, MA 01803

SigmaAldrich.com

© 2021 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. MilliporeSigma, the vibrant M, Chromolith and Supelco are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

