

## 77981 CT Supplement (Tellurite Cefixime Supplement, Cefixime Tellurite Supplement)

A selective supplement recommended by ISO Committee for the isolation of *Escherichia coli* 0157.

### Composition

(one vial, sufficient for 500 ml medium)

Potassium tellurite 1.25mg

Cefixime 0.025mg

### Directions

Rehydrate the contents of 1 vial aseptically with 5 ml of sterile distilled water. Mix gently to dissolve the contents completely. Aseptically add the contents to 495 ml of sterile molten, cooled to 45-50°C MacConkey Sorbitol Agar or MacConkey Sorbitol *ChromoSelect* Agar (Cat. No. 88474). Mix well and pour into sterile petri plates.

### Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

