

3050 Spruce Street Saint Louis, Missouri 63103 USA Telephone 800-325-5832 • (314) 771-5765 Fax (314) 286-7828 email: techserv@sial.com sigma-aldrich.com

ProductInformation

Anti- phospho-Tau (pSer⁴⁰⁴)

Developed in Rabbit, Affinity Isolated Antibody

Product Number T 7444

Product Description

Anti-phospho-Tau (pSer⁴⁰⁴) is developed in rabbit using a synthetic phosphopeptide derived from the region of tau that contains serine 404 as immunogen. The serum is affinity purified using epitope-specific affinity chromatography. The antibody is preadsorbed to remove any reactivity towards a non-phosphorylated tau.

Anti-phospho-Tau (pSer⁴⁰⁴) recognizes human, mouse and rat tau (pSer⁴⁰⁴) (45-68 kDa). It has been used in immunoblotting applications.

Tau is a microtubule-associated phosphoprotein (MAP), localized in neuronal axons. It promotes tubulin polymerization and stabilizes microtubules.¹ The biological activity of tau is regulated by its degree of phosphorylation.^{1,2} Hyperphosphorylated tau is the major protein of the paired helical filaments (PHFs), which make up the pathological neurofibrillary tangles of Alzheimer's disease (AD). The PHFs are also found in the lesions of other central nervous system disorders.^{3,4}

Tau phosphorylation involves numerous kinases: glycogen synthase kinase 3 β (GSK-3 β), MARK kinase, MAP kinase, protein kinase A and C, cyclin-dependent kinase 5 (Cdk5), p38 kinase, c-Jun N-terminal kinase, and casein kinase II.^{1,2,5,6,7} Combined tau protein kinase II (TPKII), which consists of Cdk5 and GSK-3 β , is the most potent phosphorylation agent indirectly involved in the regulation of the phosphorylation state of tau in neuronal cells.^{6,8} In addition, tau is phosphorylated *in vitro* by osmotic cellular stress, which activates the stress-activated protein kinases (SAPKs).

To date, a total of 25 abnormal phosphorylation sites have been identified on hyperphosphorylated tau in an AD brain.¹⁰ Normal tau has approximately eight phosphorylation sites. The abnormal phosphorylation occurs usually on serine and threonine residues. Specifically, TPKII phosphorylates serines 202 and 404. GSK-3 β transfection phosphorylates serines 199, 202, 235, 396, 404 and 413, and threonines 205 and 231. These sites are among the major abnormal phosphorylation sites of tau.¹¹ Phosphorylation on these sites reduces the ability of a given tau species to promote microtubule self-assembly.^{11,12} Okadaic acid increases phosphorylation at threonine 231 and serines 235, 396 and 404. Phosphorylated serine 422 was found in the biopsies of brains from patients with Down syndrome, amyotropic lateral sclerosis, corticobasal degeneration, and Pick's disease. It was absent from control group of normal brains.¹³

The opposite process, tau dephosphorylation, is controlled by different protein phosphatases expressed in neurons. Protein phosphatases PP2A and PP2B efficiently dephosphorylate tau *in vitro* and restore biological activity in the assembly of microtubules.^{3,10,14}

Recently it was discovered that propyl isomerase (Pin 1) interacts with tau hyperphosphorylated on threonine 231 and restores the ability of tau to bind to microtubules.

Reagent

The antibody is supplied at approximately 0.5 mg/ml in Dulbecco's phosphate buffered saline (without Mg^{2+} and Ca^{2+}), pH 7.3, with 50% glycerol, 1.0 mg/ml BSA (IgG and protease free), and 0.05% sodium azide.

Precautions and Disclaimer

Due to the sodium azide content, a material safety data sheet (MSDS) for this product has been sent to the attention of the safety officer of your institution. Consult the MSDS for information regarding hazards and safe handling practices.

Storage/Stability

Store at -20 °C. For extended storage, upon initial thawing, freeze in working aliquots. Avoid repeated freezing and thawing to prevent denaturing the antibody. Working dilution samples should be discarded if not used within 12 hours.

Product Profile

One vial is sufficient for 10 immunoblots

A recommended working dilution of 1: 1000 is determined by immunoblotting using cell extracts from African green monkey kidney (CV-1). Data shows the specificity of Anti-phospho-Tau (pSer⁴⁰⁴) for the phosphorylated protein since no blocking is observed in the presence of the non-phosphorylated immunizing peptide.

Note: In order to obtain best results in different techniques and preparations we recommend determining optimal working concentration by titration test.

Results

1	kDa	2	3	4	5
	250				
	150				
	100	_	_	_	
	75				
	50				
	37				
				-	
None (-)		None	Tau [Gene	Tau [
÷		None (+)	pS40	Generic pS	Tau [pS ⁴⁰⁴])
			Tau [pS ⁴⁰⁴]NP	Š)4]₽

Figure 1 Peptide Competition

- Human recombinant tau added to background extracts and left untreated (lane 1) or treated with GSK (lanes 2-5) were resolved by SDS-PAGE on a 10% polyacrylamide gel and transferred to PVDF.
- 2. Membranes were blocked with a 5% BSA-TBST buffer for 1 hour at room temperature

- 3. After blocking, membranes were preincubated with different peptides as follow:
 - Lane 1, 2 no peptide
 - Lane 3 non phosphorylated peptide corresponding to the immunogen
 - Lane 4 a generic phosphoserine containing peptide
 - Lane 5 immunogen
- After preincubation membranes were incubated with 0.50 μg/mL Anti-tau [pSer⁴⁰⁴] antibody for two hours at room temperature in a 3% BSA-TBST buffer.
- 5. After washing, membranes were incubated with goat $F(ab')_2$ anti-rabbit IgG alkaline phosphatase and signals were detected.

The data in Figure 1 show that only the peptide corresponding to Tau [pSer⁴⁰⁴] blocks the antibody signal, thereby demonstrating the specificity of the antibody.

References

- 1. Haque, N., et al., Brain Res., 838, 69-77 (1999).
- 2. Jenkins, S.M., et al., Biochem. J., **345**, 263-270 (2000).
- 3. Gong, C.X., et al., J. Biol. Chem., **275**, 5535-5544 (2000).
- 4. Goedert, M., Nature, **399**, 739-740 (1999).
- 5. Maccioni, R.B., et al., Eur. J. Biochem., **268**, 1518-1527 (2001).
- 6. Arioka, M. et al., J. Neurochem., **60**, 461-468 (1993).
- 7. Anderton, B.H., et al., Biochem. Soc. Symp., **67**, 73-80 (2001).
- 8. Michel, G., et al., Biochem. Biophys. Acta, **1380**, 177-182 (1998).
- 9. Lundt, E.T., et al., J. Neurochem., **76**, 1221-1232 (2001).
- 10. Gong, C.X., et al., J. Neurochem., **62**, 803-806 (1994).
- 11. Takashima, A., et al., Neurosci. Res., **31**, 317-323 (1998).
- 12. Utton, M.A., et al., Biochem. J., **323**, 741-747 (1997).
- 13. Bussiere, T., et al., Acta Neuropathol. (Berlin), **97**, 221-230 (1999).
- 14. Wang, J.Z., et al., Mol. Brain Res., **38**, 200-208 (1996).

AH/PHC REV 11/04

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.

Sigma brand products are sold through Sigma-Aldrich, Inc.