

ProductInformation

SIGMA QUALITY CONTROL TEST PROCEDURE

Enzymatic Assay of TRYPSIN¹ (EC 3.4.21.4)

PRINCIPLE:

BAEE + $H_2O^{\frac{\text{Trypsin}}{\text{S}}}$ N α -Benzoyl-L-Arginine + Ethanol

Abbreviation used: BAEE = $N\alpha$ -Benzoyl-L-Arginine Ethyl Ester

CONDITIONS: $T = 25^{\circ}C$, pH = 7.6, A_{253nm} , Light path = 1 cm

METHOD: Continuous Spectrophotometric Rate Determination

REAGENTS:

- A. 67 mM Sodium Phosphate Buffer, pH 7.6 at 25°C (Prepare 100 ml in deionized water using Sodium Phosphate, Monobasic, Anhydrous, Sigma Prod. No. S-0751. Adjust to pH 7.6 at 25°C with 1 M NaOH.)
- B. 0.25 mM Nα-Benzoyl-L-Arginine Ethyl Ester Solution (BAEE) (Prepare 50 ml in Reagent A using Nα-Benzoyl-L-Arginine Ethyl Ester, Hydrochloride, Sigma Prod. No. B-4500.)
- C. 1 mM Hydrochloric Acid Solution (HCI) (Prepare 50 ml in deionized water using concentrated Hydrochloric Acid, Sigma Prod. No. H-7020.)
- D. Trypsin Enzyme Solution (Immediately before use, prepare a solution containing 500 BAEE units/ml of Trypsin in cold Reagent C.)

Enzymatic Assay of TRYPSIN¹ (EC 3.4.21.4)

PROCEDURE:

Pipette (in milliliters) the following reagents into suitable quartz cuvettes:

	Test	Blank
Reagent B (BAEE)	3.00	3.00
Equilibrate to 25° C. Monitor the A _{253nm} until constant, using a suitably thermostatted spectrophotometer. Then add:		
Reagent C (Hcl) Reagent D (Enzyme Solution)	0.20	0.20

Immediately mix by inversion and record the increase in A_{253nm} for approximately 5 minutes. Obtain the ΔA_{253nm} /minute using the maximum linear rate for both the Test and Blank.

CALCULATIONS:

 $(\Delta A_{253nm}/min \text{ Test} - \Delta A_{253nm}/min \text{ Blank})(df)$

BAEE Units/ml enzyme =

(0.001) (0.20)

df = Dilution factor

0.001 = The change in A_{253nm}/minute per unit of Trypsin at pH 7.6 at 25°C in a 3.2 ml reaction mix 0.20 = Volume (in milliliters) of enzyme used

units/ml enzyme

Units/mg solid = mg solid/ml enzyme

units/ml enzyme

Units/mg protein = -----

mg protein/ml enzyme

UNIT DEFINITION:

One BAEE unit will produce a ΔA_{253nm} of 0.001 per minute with BAEE as substrate at pH 7.6 at 25°C in a reaction volume of 3.2 ml.

FINAL ASSAY CONCENTRATION:

In a 3.2 ml reaction mix, the final concentrations are 63 mM sodium phosphate, 0.23 mM N α -benzoyl-L-arginine ethyl ester, 0.06 mM hydrochloric acid, and 100 units trypsin.

Enzymatic Assay of TRYPSIN¹ (EC 3.4.21.4)

REFERENCE:

Bergmeyer, H.U., Gawehn, K., and Grassl, M. (1974) in *Methods of Enzymatic Analysis* (Bergmeyer, H.U., ed) Volume I, 2nd ed., 515-516, Academic Press, Inc.,New York, NY

NOTES:

- 1. This assay procedure is not to be used to assay Sigma Prod. Nos. T-1763, T-4019, T-8386, T-8899, and T-9906.
- 2. This assay is based on the cited reference.
- 3. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that Sigma products conform to the information in Sigma publications. Purchaser must determine the suitability of the information and products for its particular use. Upon purchase of Sigma products, see reverse side of invoice or packing slip for additional terms and conditions of sale.