## User Guide

# Millex®-GP/HP Filter Unit (Sterile) with Millipore Express® PES Membrane

**REF** **SLGP033RS\*** **SLHP033RS**  
**SLGP033RB\*** **SLHP033RB**

### Available languages

|                                  |    |                                      |    |
|----------------------------------|----|--------------------------------------|----|
| English . . . . .                | 3  | Latviski (Latvian) . . . . .         | 30 |
| Français (French) . . . . .      | 6  | Lietuviškai (Lithuanian) . . . . .   | 33 |
| Italiano (Italian) . . . . .     | 9  | Magyar (Hungarian) . . . . .         | 36 |
| Deutsch (German) . . . . .       | 12 | Český (Czech) . . . . .              | 39 |
| Español (Spanish) . . . . .      | 15 | Polski (Polish) . . . . .            | 42 |
| Português (Portuguese) . . . . . | 18 | 日本語 (Japanese) . . . . .             | 45 |
| Nederlands (Dutch) . . . . .     | 21 | 繁體中文 (Traditional Chinese) . . . . . | 47 |
| Dansk (Danish) . . . . .         | 24 | Русский (Russian) . . . . .          | 49 |
| Svenska (Swedish) . . . . .      | 27 | Norsk (Norwegian) . . . . .          | 52 |

\* [Brazil Product Registration](#)  
[Number 81167709005.](#)  
[See page 2.](#)

PR05138w  
Rev 06/23  
Date of issue: 2023-06

[SigmaAldrich.com](https://SigmaAldrich.com)

**Millipore®**

## Definition of Symbols Used

| Symbol                                                                            | Definition                                                        | Symbol                                                                            | Definition                      |
|-----------------------------------------------------------------------------------|-------------------------------------------------------------------|-----------------------------------------------------------------------------------|---------------------------------|
|  | Sterilized using irradiation                                      |  | Temperature limit               |
|  | Non-pyrogenic                                                     |  | Use-by date                     |
|  | Consult instructions for use                                      |  | Date of manufacture             |
|  | Do not use if package is damaged and consult instructions for use |  | Manufacturer                    |
|  | Do not re-sterilize                                               |  | Caution                         |
|  | Do not re-use                                                     |  | CE conformity marking           |
|  | Catalogue number                                                  |  | Swiss authorized representative |
|  | Batch code                                                        |  | Unique device identifier        |

Made in Ireland  
 Merck Millipore Ltd.  
 Tullagreen, Carrigtwohill,  
 Co. Cork, Ireland  
 T45 KD29  
 an affiliate of Merck KGaA,  
 Darmstadt, Germany

MDSS CH GmbH  
 Laurenzenvorstadt 61  
 5000 Aarau  
 Switzerland

Brazilian Registration Holder (BRH)  
 Merck S/A  
 Rua Torre Eifel, nº 100  
 Lote 3ª, Gleba A  
 Parque Rincão  
 Cotia/SP

\* Brazil Product Registration Number  
 81167709005

## Millex®-GP/HP Filter Unit (Sterile) with Millipore Express® PES Membrane

|                   |                  |               |
|-------------------|------------------|---------------|
| Millex®-GP Filter | <b>SLGP033RS</b> | Qty: 50 pack  |
|                   | <b>SLGP033RB</b> | Qty: 250 pack |
| Millex®-HP Filter | <b>SLHP033RS</b> | Qty: 50 pack  |
|                   | <b>SLHP033RB</b> | Qty: 250 pack |

- 33 mm
- Sterile
- Single use only
- Non-pyrogenic

### Intended Purpose

The Millex®-GP/HP Filter Unit (Sterile) are medical devices intended for use as syringe filters to sterilize and/or clarify low volume solutions in direct patient care and pharmacy admixture applications. Devices are intended for medical and/or laboratory professional use only.

### Introduction

This document provides compatibility information, operating steps, and specifications for the Millipore Express® polyethersulfone (PES) family of sterile filters. These filters remove microorganisms, particles, precipitates, and undissolved powders larger than the membrane's rated pore size. These single-use filters are non-pyrogenic and non-toxic. They consist of a membrane filter sealed in a modified acrylic copolymer (MMA) housing.

### Applications

Typical research laboratory applications include the sterile filtration (GP) and/or clarification (GP/HP) of protein solutions, tissue culture media, additives, buffers, and water. Clinical benefits include the removal of microorganisms, particles, precipitates, and undissolved powders larger than the membrane's rated pore size from solutions for clinical use. In the hospital pharmacy, filters with Millipore Express® PES membrane can be used for the sterile filtration (GP) and/or clarification (GP/HP) of small volumes of protein pharmaceuticals, diagnostic imaging agents, chemotherapeutics, aqueous solutions, or water during admixture preparation. Direct patient care applications include sterilization (GP) and/or particulate removal (GP/HP) from epidural and other liquid anaesthetics as well as from irrigation solutions used in ophthalmic, otic, and other surgical procedures.

### Chemical Compatibility




These filters are compatible with most aqueous solutions. Compatibility of solution should be tested before use. Visit [SigmaAldrich.com/MedMillex](https://www.sigmaaldrich.com/MedMillex) for more information.

## WARNINGS

- Filter solutions with a temperature limit of 45 °C (113 °F).
- Syringe filters are for manual use only; do not use on automated systems.
- Not suitable for filtering high viscosity solutions or blood.
- Do not filter solutions containing 5 mg or less of active drug materials unless binding studies have been performed.
- To ensure sterility, do not use this product if the package is damaged.
- Use aseptic technique to maintain sterility removing product from primary package.
- Do not use with syringes smaller than 10 mL because pressures in excess of the maximum pressure rating may be reached, potentially causing damage to the filter and/or personal injury.
- Do not use for direct neuraxial applications which may come in contact with cerebrospinal fluid.
- Do not use the same syringe filter to filter solutions in both directions.
- Sudden loss of pressure could indicate failure of the filter.
- Single use only; do not re-use or resterilize.
- Make sure to wet the filter membrane thoroughly before injecting the solution; improperly wetted filters can become airlocked.
- Not suitable for use with medical pressure infusion equipment or gravity fed intravenous (IV) administration sets.
- Discard appropriately after single use. See "Disposal" below

## Directions for Use

1. Wearing gloves using aseptic technique: fill syringe with solution to be filtered.
2. Remove cover from package.
3. Attach suitable syringe to filter inlet and remove assembly from package.
4. Hold syringe with filter pointing up and top off by pushing a few drops through to wet filter.  
⚠ Excess solution may be hazardous and should be disposed with care.
5. Depress syringe plunger to filter solution into suitable container or compatible attachment on outlet.

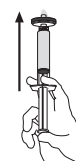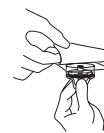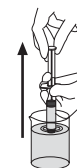## Specifications and Performance Characteristics

### Materials

|           |                                                          |
|-----------|----------------------------------------------------------|
| Membrane  | Hydrophilic Millipore Express® polyethersulfone (PES)    |
| Pore size | Millex®-GP filter: 0.22 µm<br>Millex®-HP filter: 0.45 µm |
| Housing   | Modified acrylic copolymer (MMA)                         |

### Dimensions

|                 |                     |
|-----------------|---------------------|
| Inlet to outlet | 27 mm (1.06 in.)    |
| Diameter        | 33 mm (1.30 in.)    |
| Filtration area | 4.52 cm² (0.70 in²) |

### Temperature limit

45 °C (113 °F)

### Pressure limit at 21 °C

10.3 bar (150 psi) inlet

### Filtration volume

10 mL to 200 mL

### Hold-up volume

≤0.1 mL after air purge

### Sterilization method

Gamma irradiation

### Connections

Female Luer-Lok™ inlet;  
male Luer-slip outlet

### Flow rate at 2.1 bar (30 psi), 21 °C

Millex®-GP filter: ≥ 150 mL/min  
Millex®-HP filter: ≥ 300 mL/min

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

## Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://SigmaAldrich.com/offices).

## Disposal

Follow precautions for disposal of items contaminated with potentially infectious or hazardous biological material according to all applicable international, federal, state, and local regulations.

## Technical Assistance

Visit the tech service page on our web site at [SigmaAldrich.com/techservice](https://SigmaAldrich.com/techservice).

Any serious incident of this device should be reported to manufacturer and competent authority of country where user is established.

## Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://SigmaAldrich.com/terms).

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---

Millipore, Millex, Durapore and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck operates  
as MilliporeSigma in the U.S. and Canada.