

## Product Information

## Caspase 3 Assay Kit, Colorimetric

## CASP3C

## Product Description


Caspases (Cysteine-requiring aspartate protease) are a family of proteases that mediate cell death and are important to the process of apoptosis. Caspase 3 (also referred to as CPP32, Yama, and apopain) is a member of the CED-3 subfamily of Caspases and is one of the critical enzymes of apoptosis.

Caspase 3, which is an effector Caspase, is the most studied of mammalian Caspases. Caspase 3 can process pro-Caspase 2, 6, 7, and 9 and specifically cleave most of the Caspase-related substrates known to date, including many key proteins such as:

- The nuclear enzyme poly(ADP-ribose) polymerase (PARP)<sup>1</sup>
- The inhibitor of Caspase-activated deoxyribonuclease (ICAD)<sup>2</sup>
- Gelsolin and fodrin, which are proteins involved in apoptosis regulation<sup>3</sup>

This cleavage is part of the mechanism leading to cell death. In addition, Caspase 3 plays a central role in mediating nuclear apoptosis, including chromatin condensation and DNA fragmentation, as well as cell blebbing.<sup>5</sup> Caspase 3 activity is tissue-, cell type-, or death stimulus-specific.<sup>5</sup>

The Caspase 3 colorimetric assay is based on the hydrolysis by Caspase 3 of the peptide substrate acetyl-Asp-Glu-Val-Asp *p*-nitroanilide (Ac-DEVD-pNA), to give the release of a *p*-nitroaniline (pNA) moiety:

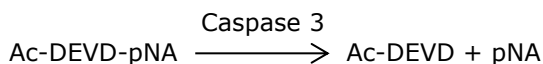

*p*-Nitroaniline has a high absorbance at 405 nm ( $E_{\text{mM}}^{\text{m}} = 10.5$ ). The concentration of the released pNA is calculated from the absorbance values at 405 nm or from a calibration curve prepared with defined pNA solutions.

This Caspase 3 Colorimetric Assay Kit provides the reagents needed for a quick and efficient detection of Caspase 3 activity in cell lysates and in purified preparations of Caspase 3.

The assay can be performed in either of these volumes:

- 1 mL volume and measured using a spectrophotometer
- 100  $\mu\text{L}$  volume in a 96-well plate using an ELISA reader.

This kit provides reagents sufficient for either:

- 100 standard 1 mL tests
- 1000 tests in 96-well plates.

## Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

## Reagents Provided

- 5 $\times$  Lysis Buffer (Component L2912): 5 mL (250 mM HEPES, pH 7.4, 25 mM CHAPS, 25 mM DTT)
- 10 $\times$  Assay Buffer (Component A0219): 20 mL (200 mM HEPES, pH 7.4, 1% CHAPS, 50 mM DTT, 20 mM EDTA)
- Caspase 3 (Component C5974, Lyophilized powder): 5  $\mu\text{g}$ 
  - Reconstitution with 50  $\mu\text{L}$  of 17 megohm water will give a solution of 100  $\mu\text{g}/\text{mL}$  Caspase 3 in 50 mM HEPES, pH 7.4, 0.1% CHAPS, 10 mM DTT, 100 mM NaCl, 1 mM EDTA and 10% sucrose.
  - Specific Activity: 1 unit per mg protein
  - Unit definition: One unit is the amount of enzyme that will cleave 1.0  $\mu\text{mol}$  of the substrate Ac-DEVD-pNA per minute at pH 7.4 at 25  $^{\circ}\text{C}$ .

- Ac-DEVD-pNA Substrate (Component A2559): 15 mg (Acetyl-Asp-Glu-Val-Asp *p*-nitroanilide)
- Ac-DEVD-CHO Inhibitor (Component A0835): 0.5 mg (Acetyl-Asp-Glu-Val-Asp-al)
- 4-Nitroaniline Standard (Component 185310): 1 mg
- Water (17 megohm) (Component W3888): 125 mL

## Reagents and Equipment Required

(Not provided)

Example Cat. Nos. are given where appropriate.

- Cells to undergo apoptosis. The example procedure here uses Jurkat E6-1 cells.
- Apoptosis inducer. Apoptosis may be either spontaneous or induced. The example procedure here uses staurosporine (Cat. No. S4400).
- Phosphate buffered saline (PBS), such as Cat. No. D8537
- DMSO, such as Cat. No. D8418
- Bovine serum albumin (BSA), such as Cat. No. A8022
- Spectrophotometer with quartz cuvettes
- ELISA reader
- Flat-bottom, 96-well plates suitable for ELISA
- Polypropylene test tubes and microcentrifuge tubes

## Storage/Stability

Store the kit at -20 °C.

## Preparation Instructions

**Note:** Use 17 megohm water only (Component No. W3888) in all the steps.

- 1× Assay Buffer: 20 mM HEPES, pH 7.4, 2 mM EDTA, 0.1% CHAPS, 5 mM DTT
  - Dilute 10× Assay Buffer 10-fold with 17 megohm water.
- Caspase 3 substrate (Ac-DEVD-pNA): 20 mM in DMSO.
  - Dissolve the vial contents (15 mg) in 1.2 mL of DMSO to prepare a stock solution.
  - Alternatively, dissolve 1 mg of substrate in 78.5 µL of DMSO. Store at -20 °C.
  - **For assays using 96-well plates**, dilute the 20 mM stock solution to **2 mM** with 1× Assay Buffer.
- Caspase 3 inhibitor (Ac-DEVD-CHO): 2 mM in DMSO.
  - Dissolve the vial contents (0.5 mg) in 500 µL of DMSO to prepare a stock solution. Store at -20 °C.
  - **For assays using 96-well plates**, dilute the 2 mM stock solution to **200 µM** with 1× Assay Buffer.
- Caspase 3 positive control:
  - Reconstitute the vial (5 µg) with 50 µL of 17 megohm water (100 µg/mL).
  - Store in aliquots at -70 °C.
  - Just before use, dilute an aliquot to 5 µg/mL 20-fold either in 1× Assay Buffer containing 1 mg/mL BSA, or 1× Lysis Buffer containing 1 mg/mL BSA.
- 1× Lysis Buffer: 50 mM HEPES, pH 7.4, 5 mM CHAPS, 5 mM DTT:
  - Dilute 5× Lysis Buffer 5-fold with 17 megohm water.
  - **Note:** In order to protect the cell lysate Caspases from non-specific proteolysis, protease inhibitor cocktails that omit particular inhibitors of cysteine proteases (such as E-64, or leupeptin) may be added.

- *p*-Nitroaniline Standard (for developing a calibration curve for assays in 96-well plate):
  - Dissolve the vial in 0.72 of mL DMSO. Store this stock solution at  $-20^{\circ}\text{C}$ .
  - To determine the actual concentration, dilute a sample of the stock solution 100-fold in  $1\times$  Assay Buffer.
  - Determine absorbance at 405 nm using a quartz cuvette.
  - Calculate the actual concentration of the stock solution, using the molar absorptivity of  $E^{\text{mM}} = 10.5$  at 405 nm.
- Staurosporine: 1 mg/mL in DMSO

## Procedure

The following procedure is an example using Jurkat cells induced to apoptosis using staurosporine and lysed prior to the determination of Caspase 3 activity.

Three controls are recommended for each Caspase 3 Colorimetric Assay:

- Inhibitor-treated cell lysate control (for measuring the nonspecific hydrolysis of the substrate)
- Caspase 3 positive control
- Reagent blank (negative control)

## Preparation of Cell Lysates from Apoptotic Cells

- 1.1.1. Induce apoptosis in a cell suspension of Jurkat cell (at least  $10^7$  cells) by addition of staurosporine to a final concentration of  $1\text{ }\mu\text{g/mL}$ . Reserve a sample of non-induced cells for a zero-time control.
- 1.1.2. Incubate for 2.5 to 3 hours at  $37^{\circ}\text{C}$  in a  $5\%$   $\text{CO}_2$  atmosphere.
- 1.1.3. Pellet the induced cells and the control cells by centrifugation at  $600 \times g$  for 5 minutes at  $4^{\circ}\text{C}$ .
- 1.1.4. Remove the supernatant by gentle aspiration.
- 1.1.5. Wash the cell pellets once with 1 mL of PBS. Centrifuge the cells and remove the supernatant completely by gentle aspiration.
- 1.1.6. Suspend the cell pellets in  $1\times$  lysis buffer at a concentration of  $100\text{ }\mu\text{L}$  per  $10^7$  cells.
- 1.1.7. Incubate the cells on ice for 15-20 minutes.

- 1.1.8. Centrifuge the lysed cells at  $16,000$  to  $20,000 \times g$  for 10-15 minutes at  $4^{\circ}\text{C}$ .
- 1.1.9. Transfer the supernatants to new tubes.
- 1.1.10. Analyze the lysates immediately, or freeze in liquid nitrogen and store in aliquots at  $-70^{\circ}\text{C}$ .

## 1 mL Volume Assay Method

- Equipment required:
    - Test tubes
    - Spectrophotometer
    - 1 mL quartz cuvettes
  - The positive control volume recommended in the reaction scheme is compatible with the expected activity found in  $0.5 \times 10^6 - 1.5 \times 10^6$  apoptotic Jurkat cells. The positive control and sample volumes can be increased if required.
  - Use quartz cuvettes **only**, since plastic cuvettes attenuate the absorption at 405 nm.
  - Yellowish color is visualized by the naked eye at approximately 0.2 Optical Density (OD) at 405 nm.
1. Place  $10\text{ }\mu\text{L}$  of cell lysate or Caspase 3 Positive Control in the appropriate tubes as indicated in Table 1.
  2. Add  $1\times$  Assay Buffer to each of the tubes as indicated in Table 1.
  3. Add  $10\text{ }\mu\text{L}$  of Caspase 3 inhibitor to the appropriate tubes.
  4. Start the reaction by adding  $10\text{ }\mu\text{L}$  of Caspase 3 substrate to each tube.
    - 4.1. Mix gently.
    - 4.2. Cover the tubes.
    - 4.3. Incubate at  $37^{\circ}\text{C}$  for 1.5 to 2 hours.
      - 4.3.1. If the signal is too low, continue the incubation overnight.
    - 4.4. Read the Absorbance at 405 nm.

- Calculate the Caspase 3 activity in  $\mu\text{mol}$  of pNA released per min per mL of cell lysate or positive control based on the formula:

- Activity, in  $\mu\text{mol pNA/min/mL} = (\text{OD} \times \text{D}) / E^{\text{mM}} \times \text{T} \times \text{V}$
- Where:
  - $E^{\text{mM}} = 10.5$
  - V = volume of sample in mL
  - D = dilution factor
  - T = reaction time in minutes

## 96-Well Plate Microassay Method

- Equipment required:
    - Flat bottom 96-well plate
    - ELISA reader
  - The positive control volume recommended in the reaction scheme is compatible with the expected activity found in  $0.25 \times 10^6$  to  $1 \times 10^6$  apoptotic Jurkat cells. The positive control and sample volumes can be increased if required.
  - Yellowish color is visualized by the naked eye at approximately 0.2 OD at 405 nm (OD405).
- Place 5  $\mu\text{L}$  of cell lysate or Caspase 3 Positive Control in the appropriate wells as indicated in Table 2.
  - Add 1 $\times$  Assay Buffer to each of the wells as indicated in Table 2.
  - Add the Caspase 3 Inhibitor to the appropriate wells.
  - Start the reaction by adding 10  $\mu\text{L}$  of Caspase 3 substrate to each well.
    - Mix gently by shaking.
    - Try to avoid forming bubbles in the wells.
    - Cover the plate.
    - Incubate at 37 °C for 70-90 minutes.
      - If signal is too low, continue the incubation overnight.
  - Read the Absorbance at 405 nm.
  - Calculate the results using a *p*-nitroaniline calibration curve. This method is recommended for accurate results to avoid miscalculations that stem from incompatibility of the ELISA reader and the plastic plates.

## *p*-Nitroaniline (pNA) Calibration Curve

- Prepare a series of *p*-nitroaniline solutions at the concentration range of 10  $\mu\text{M}$ –200  $\mu\text{M}$  by diluting the *p*-nitroaniline stock solution in 1 $\times$  Assay Buffer.
- Add 100  $\mu\text{L}$  of each dilution to a well.
  - Include 100  $\mu\text{L}$  of assay buffer as a blank.
- Read the Absorbance at 405 nm.
- Prepare a calibration curve of the absorbance values versus the concentrations of the *p*-nitroaniline solutions.
  - Alternatively, plot the OD405 values versus the amount of *p*-nitroaniline per well, in  $\mu\text{mol}$ , using the following values as a guideline.

## *p*-Nitroaniline Calibration Curve

$\mu\text{M}$ <i>p</i> -Nitroaniline	$\mu\text{mol}$ <i>p</i> -Nitroaniline per 100 $\mu\text{L}$
10	0.001
20	0.002
50	0.005
100	0.01
200	0.02

- Calculate the Caspase 3 activity in  $\mu\text{mol pNA}$  released per min per mL of cell lysate or positive control, based on the formula:

- Activity,  $\mu\text{mol pNA/min/mL} = (\mu\text{mol pNA} \times \text{D}) / (\text{T} \times \text{V})$
- Where:
  - V = volume of sample in mL
  - D = dilution factor
  - T = reaction time in minutes

## References

- Nicholson, D.W. *et al.*, *Nature*, **376(6535)**, 37-43 (1995).
- Sakahira, H. *et al.*, *Nature*, **391(9962)**, 96-99 (1998).
- Kamada, S. *et al.*, *Proc. Natl. Acad. Sci. USA*, **95(15)**, 8532-8537 (1998).
- Cohen, G.M., *Biochem. J.*, **326(Pt 1)**, 1-16 (1997).
- Porter, A.G., and Jänicke, R.U., *Cell Death Differ.*, **6(2)**, 99-104 (1999).

**Table 1.** Reaction scheme for 1 mL Volume Assay Method

	<b>Cell lysate</b>	<b>Caspase 3 (5 µg/mL)</b>	<b>Assay buffer</b>	<b>Caspase 3 inhibitor Ac-DEVD-CHO (2 mM)</b>	<b>Caspase 3 substrate Ac-DEVD-pNA (20 mM)</b>
Reagent blank	----	----	990 µL	----	10 µL
Non-induced cells	10 µL	----	980 µL	----	10 µL
Non-induced cells + inhibitor	10 µL	----	970 µL	10 µL	10 µL
Induced cells	10 µL	----	980 µL	----	10 µL
Induced cells + inhibitor	10 µL	----	970 µL	10 µL	10 µL
Caspase 3 positive control	----	10 µL	980 µL	----	10 µL
Caspase 3 positive control + inhibitor	----	10 µL	970 µL	10 µL	10 µL

**Table 2.** Reaction scheme for 96-Well Plate Microassay Method

	<b>Cell lysate</b>	<b>Caspase 3 (5 µg/mL)</b>	<b>1× Assay buffer</b>	<b>Caspase 3 inhibitor Ac-DEVD-CHO (200 µM)</b>	<b>Caspase 3 substrate Ac-DEVD-pNA (2 mM)</b>
Reagent blank	----	----	90 µL	----	10 µL
Non-induced cells	5 µL	----	85 µL	----	10 µL
Non-induced cells + inhibitor	5 µL	----	75 µL	10 µL	10 µL
Induced cells	5 µL	----	85 µL	----	10 µL
Induced cells + inhibitor	5 µL	----	75 µL	10 µL	10 µL
Caspase 3 positive control	----	5 µL	85 µL	----	10 µL
Caspase 3 positive control + inhibitor	----	5 µL	75 µL	10 µL	10 µL

---

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

## Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://SigmaAldrich.com/techservice).

## Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at [SigmaAldrich.com/terms](https://SigmaAldrich.com/terms).

## Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://SigmaAldrich.com/offices).

The life science business of Merck operates  
as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates.  
All other trademarks are the property of their respective owners. Detailed information on  
trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

CASP3Cpis Rev 07/22 LPG,SS,AC,HM,GCY,MAM

