PyroDetect Cryoblood

Artikel-Nr. 1.44155.0001

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PyroDetect Cryoblood

Article No. 1.44155.0001

Cryo-preserved human whole blood for detection of pyrogens with the Monocyte-Activation Test (MAT). To use with the PyroDetect Kit (Art. No. 1.44154.0001).

The PyroDetect Cryoblood is part of the PyroDetect System (Art. No. 1.44153.0001) for the detection of pyrogens in pharmaceuticals, medical devices and cosmetics using Interleukin-1 β response.

The cryo-preserved and pooled human whole blood is tested for sterility and for antigens of HIV, HAV, HCV and HBV.

Content and Storage

PyroDetect Cryoblood (Store at -80°C or lower) Art. No. 1.44155.0001	Content	Storage of opened reagents
Cryoblood	2 x 2 ml	Do not freeze after thawing

Expiration

For expiration of the PyroDetect Cryoblood please refer to the expiration date of the label. Don't use cryoblood past expiration date!

Warning and Precautions

For research use only, not for in vitro diagnostic use

Caution: The cryoblood should be handled by well-trained and authorized laboratory technicians in accordance with good laboratory practice.

Caution: We recommend the vaccination of personnel with direct contact to the cryoblood.

Caution: Do not use cryoblood past the expiration date printed on the label.

Caution: Do not heat cryoblood over 37°C .

Caution: After melting of the cryoblood mix with cell culture medium (RPMI) of the PyroDetect Kit for stabilization of monocytes in blood.

Preparation

The PyroDetect Cryoblood is to be used in combination with the PyroDetect Kit.

For the preparation of the cryoblood please refer to the user manual of the PyroDetect Kit.