

83055 BSM-Supplement

Supplement for the selective isolation and identification of Bifidobacteria like *Bifidobacterium longum* and *Bifidobacterium infantis*. The BSM-Supplement is a mixture of antibiotics and inhibits other bacteria.

Store in the dark. Some components are photosensitive.

Directions

Suspend 0.116g of the BSM Supplement in 1 to 5 ml sterile water, mix well. The supplement is faintly soluble but it can be added to the medium as a turbid and white suspension. Add the supplement to 1 litre sterile BSM agar (Cat. No. 88517) or BSM broth (Cat. No. 90273) at 55°C and mix well.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

