

OxyBlot[™] Protein Oxidation Detection Kit

Catalog No. S7150

FOR RESEARCH USE ONLY Not for use in diagnostic procedures

USA & Canada Phone: +1(800) 645-5476 In Europe, please contact Customer Service: France: 0825.045.645; Spain: 901.516.645 Option 1 Germany: 01805.045.645 Italy: 848.845.645 United Kingdom: 0870.900.46.45 For other locations across the world please visit <u>www.millipore.com/offices</u> This page left blank intentionally

TABLE OF CONTENTS

١.	INTRODUCTION	1
	Using this Manual	1
	Background	
	Principles of the Procedure	2
П.	KIT COMPONENTS	
	Materials Required But Not Supplied	
	Precautions	
Ш.	PROTOCOL	
	Fig. 1: Flow Diagram Of OxyBlot™ Protocol	
	Preparation of Protein Lysate	
	Derivatization of Protein Mixture	
	Reagents	
	Protocol	
	SDS-PAGE, Western Transfer and Immunodetection	7
	Reagents	7
	Protocol	8
IV.	DATA ANALYSIS	10
	Molecular Weight Protein Standards	
	Fig. 2: Molecular Weight Protein Standard	
V.	TROUBLESHOOTING	11
VI.	Appendix	13
	Preparation of Oxidized Proteins for Positive Controls	
	Reagent Preparation	
	Sensitivity of DNP Detection	
	Fig. 3: Sensitivity of DNP Detection Using the OxyBlot™ Kit	15
VII.	References	16
	References Cited in the Manual	
	Disclaimers	
	Warranty	19

For Research Use Only; Not for use in diagnostic procedures

I. INTRODUCTION

Using This Manual

OxyBlot[™] Protein Oxidation Detection Kit (S7150) provides the chemical and immunological reagents necessary to perform the immunoblot detection of carbonyl groups introduced into proteins by oxidative reactions with ozone or oxides of nitrogen or by metal catalyzed oxidation. Please read the entire manual before beginning the procedure. Should additional questions arise, contact Techical Support for assistance.

Background

Oxygen-derived free radicals have been implicated in important roles in aging, apoptosis, cancer, neurodegenerative diseases, chronic inflammatory diseases, pulmonary diseases, and cardiovascular diseases (for reviews, see ref. 1-4).

Oxygen free radicals are generated by environmental factors such as ionizing radiation and chemical substances. They are also produced during normal cellular metabolism by mitochondrial electron transport and the cellular redox system, and by immune responses (5-7). Although cells have developed various antioxidant defenses to protect against free radicals (7-9), free radicals that escape the defenses attack and modify subcellular components – nucleic acids, lipids and proteins (7, 10-12). In some cases, cells respond to the oxidative stimuli and allow the organism to adapt to the oxidative stress (13-15).

Proteins are one of the major targets of oxygen free radicals and other reactive species. Oxidative modification of proteins modifies the side chains of methionine, histidine, and tyrosine and forms cysteine disulfide bonds (16-19). Metal catalyzed oxidation of proteins introduces carbonyl groups (aldehydes and ketones) at lysine, arginine, proline or threonine residues in a site-specific manner (16, 20-22).

The oxidative modification of proteins can modulate biochemical characteristics of proteins such as enzymatic activity (21-23), DNA binding activities of transcription factors (26-28), and the susceptibility to proteolytic degradation (12, 27-30). While a relationship between protein oxidation and aging has been suggested (31-33), little is known about the importance of oxidative modification of individual proteins in the pathophysiology of free radical mediated processes.

OxyBlot[™] Protein Oxidation Detection Kit provides a sensitive and simple methodology for detection and quantification of proteins modified by oxygen free radicals and other reactive species.

Principles of the Procedure

Oxidative modification of proteins by oxygen free radicals and other reactive species such as hydroxynonenal occurs in physiologic and pathologic processes. As a consequence of the modification, carbonyl groups are introduced into protein side chains by a site-specific mechanism. The OxyBlot[™] Kit provides reagents for simple and sensitive immunodetection of these carbonyl groups, which is a hallmark of the oxidation status of proteins.

The carbonyl groups in the protein side chains are derivatized to 2,4-dinitrophenylhydrazone (DNPhydrazone) by reaction with 2,4-dinitrophenylhydrazine (DNPH). The DNP-derivatized protein samples are separated by polyacrylamide gel electrophoresis followed by Western blotting. Alternatively, the derivatized protein samples can be directly dot blotted onto a membrane filter. After either method of preparation, the filters are incubated with primary antibody, specific to the DNP moiety of the proteins. This step is followed by incubation with a horseradish peroxidase-antibody conjugate directed against the primary antibody (secondary antibody: goat anti-rabbit IgG). The filters are then treated with chemiluminescent reagents (luminol and enhancer). The luminol is converted to a light-emitting form at wavelength 428 nm by the antigen/primary antibody/secondary antibody/peroxidase complex (bound to the membrane) in an H₂O₂ catalyzed oxidation reaction. The light is detected by short exposure to blue-light sensitive films.

Under the conditions recommended in the kit, as little as 5 femtomoles of carbonyl residue can be detected. This sensitivity is at least 100 times greater than those obtained by other procedures such as radioisotope methodology utilizing ³H-labeled NaBH₄.

The detection of oxidatively modified proteins by immunoblotting has additional advantages. Individual oxidized proteins are separated and identified from a complex mixture by SDS-PAGE. The oxidative status of each protein can be analyzed quantitatively by comparison of the signal intensity of the same protein in different lanes on the same or different gels.

II. KIT COMPONENTS

Part #	Component	Volume	Storage Temp.
90448	1X 2,4-Dinitrophenylhydrazine (DNPH) Solution	4.0 mL	2° to 8°C
90449	Neutralization Solution	3.0 mL	2° to 8°C
90450	Mixture of Standard Proteins with attached DNP residues	75 µL	-25° to -15°C
90451	1° Antibody: Rabbit Anti-DNP Antibody	4.0 mL	2° to 8°C
90452	2° Antibody: Goat Anti-Rabbit IgG (HRP-conjugated)	2.0 mL	2° to 8°C
90463	1X Derivatization-Control Solution	4.0 mL	2° to 8°C

S7150 OxyBlot[™] Protein Oxidation Detection Kit

The reagents supplied in the kit are sufficient to process 200 samples (5 µL each) on 30 acrylamide gels (7 samples with derivatized and control reactions; 1.5 mm gel thickness), assuming 15 mL of diluted primary and secondary antibody solution are used for treating each membrane.

Materials Required But Not Supplied

(Refer to Sec. VI. *Appendix* for specific instructions regarding the preparation of some of these reagents).

- 1. Transfer Buffer (12 mM Tris, 96 mM Glycine, 20% Methanol)
- 2. 1X PBS-T (PBS, pH 7.2-7.5, containing 0.05% Tween 20)
- 3. Blocking/Dilution Buffer (1% BSA/PBS-T)
- 4. 1X Gel Loading Buffer
- 5. 12% SDS
- 6. 0.5-1.5 mL Eppendorf Tubes


- 7. Reagents necessary to perform SDS-PAGE
- 8. Transfer Apparatus
- 9. Chemiluminescent Reagent
- 10. Nitrocellulose or PVDF Membrane
- 11. Distilled water (dH₂O)

Precautions

- 1. Components 90448 and 90463 contain 2N HCL acid. Component 90448 also contains 2-4dinitrophenyl hydrazine (DNPH). Harmful if swallowed or inhaled; avoid contact with skin and eyes (wear gloves and eye protection); wash areas of contact immediately.
- 2. Store the protein standard (component 90450) at -25° to -15°C upon receipt. Store all other components at 2°C to 8°C.

III. PROTOCOL

Figure 1: Flow Chart of OxyBlot™ Procedure.

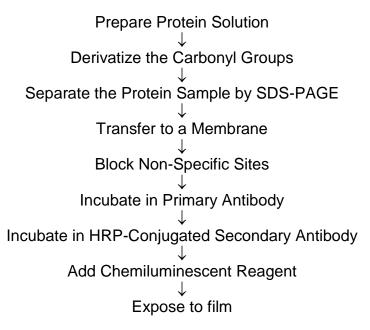

Preparation of Protein Lysate

The protein lysate may be prepared using a variety of protocols, such as those detailed in ref. 38 and 39. It is recommended that a reducing agent be added to the lysis buffer to prevent the oxidation of proteins that may occur after cell lysis. Lysis buffer containing either 1-2% 2-mercaptoethanol or 50 mM DTT sufficiently inhibits this oxidation, but has no adverse effect on the derivatization reaction in the OxyBlotTM protocol. Purified protein samples are also suitable for analysis with the OxyBlotTM Kit.

Derivatization of Protein Mixture

Reagents

- a. 1X DNPH Solution
- b. 1X Derivatization-Control Solution
- c. Neutralization Solution
- d. 12% SDS

Protocol

Two aliquots of each specimen to be analyzed will be treated. One aliquot will be subjected to the derivatization reaction. The other aliquot will serve as a negative control, by substituting the 1X Derivatization-Control Solution for the 1X DNPH Solution.

 Transfer 5 μL of a protein sample (crude or purified) into each of 2 eppendorf tubes (0.5-1.5 mL). The volume of protein samples can vary depending on the well size of the gel and the concentration of the sample. When a sample volume other than 5 μL is used, adjust the volume of other reagents accordingly.

Note: Using the protocol and reagents of this kit, 10 femtomoles of DNP residues in a protein will be readily detectable by a 30 second exposure (see Appendix). This is about equal to those in 2.5 ng of protein with M.W. 50,000 containing 0.2 mole carbonyl group/mole of protein.

Note: Generally, 15-20 µg of protein is recommended per derivatization reaction. Users may need to try several protein concentrations to obtain desirable results.

Note: Avoid high protein concentration (>10 μ g/ μ L) to assure solubility during the derivatization reaction.

- 2. Denature each 5 μ L aliquot of protein by adding 5 μ L of 12% SDS for a final concentration of 6% SDS.
- Derivatize the sample by adding 10 μL of 1X DNPH Solution to one of the tubes. To the aliquot designated as the negative control, add 10 μL of 1X Derivatization-Control Solution instead of the DNPH solution.
- 4. Incubate both tubes at room temperature for 15 minutes. Do not allow the reaction to proceed more than 30 minutes, as side reactions other than hydrazone linkage may occur.
- 5. Add 7.5 μ L of Neutralization Solution to both tubes. If reduction of the protein sample is desired and a reducing reagent was not present during lysis, add 2-mercaptoethanol to the sample mixture to achieve a final concentration of 0.74 M solution (1-1.5 μ L; 5%v/v).
- 6. Both the treated sample and the negative control are ready to load into a polyacrylamide gel.

Note: At this point, the samples can be stored at 4°C for up to one week. Thaw at room temperature.

SDS-PAGE, Western Transfer and Immunodetection

Reagents

- a. Transfer Buffer
- b. Blocking/Dilution Buffer: 1% BSA/PBS-T
- c. PBS-T: PBS (Phosphate buffered saline, pH 7.2-7.5) containing 0.05% Tween-20.
- d. DNP-lated molecular weight standard protein (90450).
- e. 1° Antibody: dilute 1° Antibody stock (90451) to 1:150 with Blocking/ Dilution Buffer
- f. 2° Antibody: dilute 2° Antibody stock (90452) to 1:300 with Blocking/ Dilution Buffer
- g. Chemiluminescent reagent: Refer to the manufacturer's instructions for preparation. Prepare the reagent just prior to use. About 0.125 mL of reagent per cm² of membrane is sufficient.

Protocol

 Upon first use, warm the standard protein mixture to room temperature. Transfer the amount necessary for the first experiment into one tube and divide the remaining mixture among several tubes. Store the aliquots at from excessive freeze-thaw and heating cycles. Combine 2.5 μL of the molecular weight standard with attached DNP residues and 20 μL of 1X Gel Loading Buffer prior to loading onto the gel. (See Sec. IV. *Data Analysis, Molecular Weight Protein Standards.*) Perform SDS-PAGE according to standard procedures (34).

Note: It is not necessary to add 1X Gel Loading Buffer to the samples prior to loading the gel, since the addition of the Neutralization Solution makes the sample dense enough to sink to the bottom of the well. However, the addition of Loading Buffer will not adversely affect the electrophoresis of the sample.

Note: Do not heat samples prior to loading into the gel.

Note: The DNP-BSA in 2.5 µL solution contains 100 femtomoles of DNP residues (determined spectrophotometrically).

Note: A portion or all of the derivatized and neutralized sample can be applied directly on SDS-PAGE or dot blotted onto a filter. The free DNPH in the derivatization mixture does not hamper the following detection reaction.

2. Electroblot proteins to a membrane (35).

Note: Refer to the transfer apparatus manual for specific transfer instructions and recommendations.

Note: If using PVDF membrane, pre-wet with 95% ethanol, then rinse with dH_2O .

Note: Optional: If non-reduced protein samples will be analyzed, soak the gel in 1% 2mercaptoethanol in 1X Gel Running Buffer prior to Western transfer. This pretreatment will enhance the detection sensitivity (36).

3. Incubate the membrane in Blocking/Dilution Buffer for 1 hour with gentle shaking. Use 0.3 mL/cm² of membrane.

- 4. Dilute the 1° Antibody stock 1:150 in Blocking/Dilution Buffer just before use. Use enough solution so that the membrane is completely immersed in the solution (15 mL for a 10 x 10 cm² membrane). An alternative method would be to place the membrane in a heat-sealable plastic bag (in this case, the amount of solution required will be less than that used for a membrane treated in a open container). Incubate the membrane in the 1° Antibody solution for ~1 hour at 18° to 25°C with gentle shaking.
- 5. Rinse the membrane 2 times with 1X PBS-T. Wash the membrane with 1X PBS-T once for 15 minutes, then twice for 5 minutes each at 18° to 25°C.
- Dilute the 2° Antibody stock 1:300 in Blocking/Dilution Buffer. Incubate the membrane in the 2° Antibody solution (15 mL for a 10 x 10 cm² membrane) for ~1 hour at 18° to 25°C with gentle shaking.
- 7. Wash the membrane as in step 5.
- 8. Drain the excess buffer from the membrane, place it on a plastic sheet with the protein side up, and cover the membrane. Prepare the chemiluminescent reagents according to manufacturer's directions just before use. Make just enough to cover the membrane with chemiluminescent reagent. The final reagent volume required is 0.125 mL/cm². Incubate for 1 minute at 18° to 25°C.
- 9. Drain off excess reagent by holding the membrane vertically and touching the edge against filter paper. As an alternative, the membrane may be dried by placing it between sheets of filter paper. Place the membrane inside a plastic page protector. Remove air pockets.
- 10. Place the membrane, protein side up, in a film cassette. Make sure that there is no free reagent in the cassette.
- 11. In a darkroom, carefully place an autoradiography film on top of the membrane.
- 12. Expose the film for 30 seconds, then develop. When all the procedures are correctly performed, the bands of molecular weight standard proteins should be readily visible after a 30 second exposure.
- 13. Place a second film on top of the membrane. Expose the film for the appropriate amount of time. Because each protein will differ in efficiency of transfer to a membrane and in the immunodetection of the attached DNP moieties, the extent of the carbonyl content in each protein is best investigated by comparison of the signal intensity of the same protein. The signal intensity of the molecular weight protein standards provides a convenient control for the experimental variations between separate experiments.

IV. DATA ANALYSIS

Proteins which have undergone oxidative modification will be identified by appearing as a band only in the lane containing the derivatized sample, but not in the lane containing the negative control. If proteins bands of the same molecular weight appear in both lanes, this is probably due to nonspecific binding of either of the antibodies and does not indicate a modified protein.

Molecular Weight Protein Standards

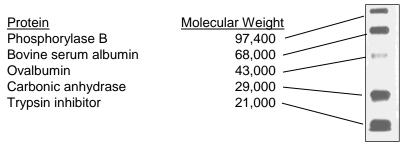


Figure 2: Molecular Weight Protein Standard.

2.5 µL of DNP-protein mixture was mixed with 20 µL of 1X SDS-Sample Buffer, separated on a 12% SDS polyacrylamide gel and electroblotted onto nitrocellulose membrane. The membrane was processed according to the procedure outlined in the OxyBlot™ Kit protocol, using reagents in the kit. The protein bands on the filter were detected using a chemiluminescent detection kit. The chemiluminescent signal was captured by a 15 second exposure on film.

The molecular weight protein standard mixture is composed of five proteins containing 1-3 DNP residues per molecule. The attached DNP residues permit the visualization of the proteins by the detection system of the kit. Since the proteins already have DNP residues, the protein standard mixture should NOT be subjected to the derivatization reaction. However, it is necessary to add 1X Gel Loading Buffer to the protein standards prior to gel loading (2.5 µL of protein standard plus 20 µL of 1X Gel Loading Buffer). In order to maintain consistency, gel loading buffer which contains a reducing agent must be added to the protein standard if the sample also contains a reducing agent. This standard serves as an internal control for the electrophoresis, western transfer and immunodetection steps of the OxyBlot™ procedure.

V. TROUBLESHOOTING



? No signal or poor signal is obtained.

1. Potential Problem: No transfer during Western Blot Procedure.

Recommendations:

- a. Check the power supply and electroblot apparatus.
- b. Check the orientation of the membrane and gel with respect to the electrodes of the transfer apparatus.
 - 2. Potential Problem: Poor transfer during Western Blot Procedure.

Recommendations:

- a. Re-evaluate the blotting procedure. (Check the transfer efficiency by staining procedure the blotting membrane (refer to ref. 37) or by transfer of pre-stained marker). Check the transfer time, current, transfer buffer, buffer temperature, gel concentration, thickness, etc.
- b. Check that the membrane was hydrated according to manufacturer's instructions.
 - 3. Potential Problem: Detection System.

Recommendations:

- a. Check that antibodies and other reagents are being stored correctly and properly used as recommended in the protocol.
- b. Make sure that chemiluminescent reagents are properly used and film developer and fixer are fresh.

? Diffused and/or distorted bands are observed.

1. Potential Problem: Electrophoresis.

Recommendations:

- a. Gel was overloaded with protein apply less electrophoresis sample to the gel.
- b. Make sure no air bubbles are trapped during electrophoresis.
- c. Check the electrophoresis conditions: make sure the buffer was made correctly.
 - 2. Potential Problem: Blotting.

Recommendations:

- a. Inspect the transfer cassette.
- b. Make sure that firm contact was made between the gel and the membrane.
- c. Make sure that any air bubbles that may have been trapped between the gel and the membrane are carefully removed prior to transfer.

High background is observed

1. Potential Problem: Inadequate Blocking of Membrane

Recommendations:

- a. Check that the Blocking/Dilution Buffer was properly prepared.
- b. Increase the incubation time of the blocking step.

VI. APPENDIX

Preparation of Oxidized Proteins for Positive Controls

1. Prepare a 5-10 mg/mL protein sample in the following buffer:

25 mM HEPES buffer, pH 7.2 25 mM ascorbate 100 μM FeCl₃

- 2. Incubate at 37°C for 5 hours.
- 3. Dialyze the sample vs. 50 mM HEPES, 1 mM EDTA.

Reagent Preparation

1. Electroblot Transfer Buffer (Half-strength Towbin Buffer)

To make 1 L:

2. 10X Phosphate Buffered Saline (PBS)

To make 1 L:

NaH₂PO₄ • H₂O (Monobasic monohydrate; FW 138.0) 26 g Na₂HPO₄ (Dibasic; FW 142.0) 11.5 g NaCl 85 g Adjust the pH to 7.2-7.5 using NaOH or HCl.

3.1X PBS-T

To make 1 L:

10X PBS	100 mL
Tween 20	0.5 mL
dH ₂ O	899.5 mL

4. Blocking/Dilution Buffer

To make 1 L:

1X PBS-T 1000 mL BSA 10 g Adjust the pH to 7.2-7.4.

5. 1X Gel Loading Buffer

To make 8 mL: 0.5 M Tris-HCl, pH 6.8 1.0 mL Glycerol 0.8 mL 14.3 M 2-Mercaptoethanol 0.1 mL (Reagent Grade) 10% SDS 1.6 mL 0.05% Bromophenol Blue0.32 mL dH₂O 4.18 mL

Divide into 1 mL aliquots and store at -25° to -15°C.

Sensitivity of DNP Detection

Figure 3: Sensitivity of DNP Detection Using the OxyBlot[™] Kit.

			1	2	3	4	5	6	7
D	NP in BSA Banc (femtomoles)	1		雨		1	-	-	
lane 1:	100					-	-	-	
lane 2:	30					-	-		
lane 3:	10					1			
lane 4:	3				-	-	-		
lane 5:	1								
lane 6:	0.3						-		
lane 7:	0.1								

Serial dilutions of DNP-protein mixture provided in the kit were analyzed as described in Sec. III Protocol. Exposure time was 1 minute.

References Cited in this Manual

- 1. Halliwell, B. and Gutteridge, J.M.C. (1990) Role of free radicals and catalytic metal ions in human diseases: an overview. *Methods in Enzymol.* **186**:1.
- 2. Halliwell, B. and Gutteridge, J.M.C. (1989) Free Radicals. *Biology and Medicine, 2nd Ed.* (Oxford, Clarendon Press).
- 3. Stadtman, E.R. (1992) Protein oxidation and aging. Science 257:1220.
- 4. Ching Kuang Chow, ed. (1988) Cellular Antioxidant Defense Mechanisms (CRC Press, FL).
- 5. Weizman, S.A. and Gordon, L.L. (1990) Inflammation and cancer: Role of phagocyte-generated oxidants in carcinogenesis. *Blood* **76**:655.
- 6. Babior, B.M. and Woodman, R.C. (1990) Chronic granulomatous disease. Semin. Hematol. 27:247.
- 7. Cance, B., et al. (1979) Hydroperoxide metabolism in mammalian organs. Physiol. Rev. 59:527.
- 8. Halliwell, B. (1994) Free radicals, antioxidants, and human disease: curiosity, cause, or consequence? *Lancet* **344**:721.
- 9. Miquel, J., et al., eds. (1989) Handbook of free radicals and antioxidant in biomedicine (CRC Press, FL).
- 10. Dizdaroglu, M. Chemistry of free radical damage to DNA and nucleoprotein in B. Halliwell and I. Aruomao, eds., (1993) *DNA and Free Radicals* (Chichester, Ellis Harwood), p19.
- 11. Ames, B.N., et al. (1991) Oxidative Damage and Repair: Chemical, Biological and Medical Aspects, K.J.A. Davies, ed. (Pergamon, Elmsford, NY), p181.
- 12.Davies, K.J.A. (1987) Protein damage and degradation by oxygen radicals. *J. Biol. Chem.* **262**:9895.
- 13.Schreck, R. *et al.* (1991) Reactive oxygen intermediates as apparently widely used messengers in the activation of the NF-kB transcription factor and HIV-1. *EMBO J.* **10**:2247.
- 14. Demple, B. and Halbrook, J.H. (1983) *Nature* **304**:466.
- 15. Christman, M. F., et al. (1985) Cell 41:753.
- 16. Stadtman, E. R. (1993) Oxidation of free amino acids and amino acid residues in proteins by radiolysis and by metal-catalyzed reactions. *Ann. Rev. Biochem.* **62**:797.
- 17.Davies, et al. (1987) Protein damage and degradation by oxygen radicals II. Modification of amino acids. J. Biol. Chem. **262**:9902.
- 18.Uchida, K. and Kawakishi, S. (1990) Site-specific oxidation of angiotensin I by copper(II) and Lascorbate: conversion of histidine residues to 2-imidazolones. *Arch. Biochem. Biophys.* **283**:20.
- 19.Heinecke, J.W., *et al.* (1993) Dityrosine, a specific marker of oxidation, is synthesized by the myeloperoxidase-hydrogen peroxidase system of human neutrophil and macrophages. *J. Biol. Chem.* **268**:4069.
- 20. Farber, J.M. and Levine, R.L. (1986) Sequence of a peptide susceptible to mixed function oxidation. Probable cation binding site in glutamine synthetase. *J. Biol. Chem.* **261**:4574.
- 21.Climent, I., *et al.* (1989) Derivatization of r-glutamyl semialdehyde residues in oxidized proteins by fluoresceninamine. *Anal. Biochem.* **182**:226.

- 22.Levine, R.L. (1983) Oxidative modification of glutamine synthetase I. Interaction is due to loss of one histidine residue. *J. Biol. Chem.* **258**:11823.
- 23.Oliver, C.N. (1987) Inactivation of enzymes and oxidative modification of proteins by stimulated neutrophils. *Arch. Biochem. Biophys.* **253**:62.
- 24.Abate, C., et al. (1990) Expression and purification of the leucine zipper and DNA-binding domaines of FOS and Jun: Both Fos and jun contact DNA directly. Proc. Natl. Acad. Sci. USA. 87:1032.
- 25.Hainaut, P. and Milner, J. (1993) Redox modulation of p53 conformation and sequence-specific DNA binding in vitro. *Cancer. Res.* **53**:4469.
- 26.Pognonec, P., *et al.* (1992) The helix-loop/leucine repeat transcription factor USF can be functionally regulated in a redox-dependent manner. *J. Biol. Chem.* **267**: 24563.
- 27.Rivett, A.J. (1986) Regulation of intracellular protein turnover: covalent modification as a mechanism of marking proteins for degradation. *Curr. Top. Cell Regul.* **28**:291.
- 28. Wolf, S.P. and Dean, R.T. (1986) Fragmentation of proteins by free radicals and its effect on their susceptibility to enzymic hydrolysis. *Biochem J.* **234**:399.
- 29. Davies, et al. (1987) Protein damage and degradation by oxygen radicals IV. Degradation of denatured protein. J. Biol. Chem. 262:9914.
- 30. Levine, R.L., *et al.* (1981) Turnover of bacterial glutamine synthetase: Oxidative inactivation proceeds proteolysis. *Proc. Natl. Acad. Sci. USA* **78**:2120.
- 31. Oliver, C.N., et al. (1987) Age related changes in oxidized proteins. J. Biol. Chem. 262:5488.
- 32. Smith, C.D., *et al.* (1991) Excess brain protein oxidation and enzyme dysfunction in normal aging and in Alzheimers disease. *Proc. Natl. Acad. Sci. USA.* **88**:10540.
- 33.Starke-Reed, P.E. and Oliver, C.N. (1989) Protein oxidation and proteolysis during aging and oxidative stress. *Arch. Biochem. Biophys.* **275**:559.
- 34.Laemmli, U.K. (1970) Cleavage of structural proteins during the assembly of the head of bacteriophge T4. *Nature* **227**:680.
- 35.Towbin, H., *et al.* (1979) Electrophoretic transfer of proteins from polyacrylamide gels to nitrocellulose sheets: procedure and some applications. *Proc. Natl. Acad. Sci. USA.* **76**:4350.
- 36.Lee, Y-J. and Shacter, E. (1995) Role of carbohydrates in oxidative modification of fibrinogen and other plasma proteins. *Arch. Biochem. Biophys.* **321(1)**:175-181.
- 37. Sasse, J. (1991) Detection of Proteins on Blot Transfer Membrane in *Current Protocols for Molecular Biology* (Ausubel, F.M., *et al*, eds.) p. 10.7.1 (John Wiley and Sons, New York).
- 38. Coligan, J. et al, eds. (1996) Current Protocols in Protein Science (John Wiley and Sons, New York).
- 39. Weissman, A.M. (1991) Solubilization of Cellular Proteins in *Current Protocols in Immunology* (Coligan, J.E. *et al*, eds.) p.8.1 (John Wiley and Sons, New York).

References Citing use of the OxyBlot kit

Wagner, K.R., *et al.* (2005) Plasma infusions into porcine cerebral white matter induce early edema, oxidative stress, pro-inflammatory cytokine gene expression and DNA fragmentation: Implications for white matter injury with increased blood-brain-barrier permeability. *Current Neurovascular Research* **2(2)**:149-155.

Agarwal, R. (2005) Smoking, oxidative stress and inflammation: Impact on resting energy expenditure in diabetic nephropathy. *BMC Nephrology* **6**:13.

Tezel, G., et al. (2005) Proteomic identification of oxidatively modified retinal proteins in a chronic pressure-induced rat model of glaucoma. *Investigative Ophthalmology & Visual Science* **46**:3177-3187.

Wu, A., et al. (2005) Dietary curcumin counteracts the outcome of traumatic brain injury on oxidative stress, synaptic plasticity, and cognition. *Experimental Neurology* in press.

Barreiro, E., et al. (2005) N-acetylcysteine increases manganese superoxide dismutase activity in septic rat diaphragms. *Eur Respir J* **26**:1032-1039.

Finn, P.F. and J.F. Dice (2005) Ketone bodies stimulate chaperone-mediated autophagy. *J. Biol. Chem.* **280**:25864-25870.

Bernhard, D., et al. (2005) Cigarette smoke metal-catalyzed protein oxidation leads to vascular endothelial cell contraction by depolymerization of microtubules. *The FASEB Journal* **19**:1096-1107.

Boyd-Kimball, D., et al. (2005) γ -Glutamylcysteine ethyl ester protection of proteins from A β (1-42)mediated oxidative stress in neuronal cell culture: A proteomics approach. *J. of Neuroscience Research* **79**:707-713.

Nagaoka, Y., et al. (2005) Specific inactivation of cysteine protease-type cathepsin by singlet oxygen generated from naphthalene endoperoxides. *BBRC* **331**:215-223.

Periquet, M. et al. (2005) Proteomic analysis of parkin knockout mice: alterations in eneary metabolism, protein handling and synaptic function. *J Neurochemistry* **95**:1259.

Elkon, H., et al. (2004) Oxidative stress, induced by 6-hydroxydopamine, reduces protease activities in PC12 cells: Implications for the pathogenesis of Parkinson's Disease. *J Molecular Neuroscience* 24:387-400.

Antony, J.M., et al. (2004) Human endogenous retrovirus glycoprotein-mediated induction of redox reactants causes oligodendrocyte death and demyelination. *Nature Neuroscience* 7:1088-1095.

Canton, M., et al. (2004) Evidence of myofibrillar protein oxidation induced by postischemic reperfusion in isolated rat hearts. *Am J Physiol Heart Circ Physiol* **286**:H870-H877.

Rowe, L.J., et al. (2004) Influence of early postmortem protein oxidation on beef quality. *J. Amin. Sci.* **82**:785-793.

Pocernich, C.B., et al. (2004) Effects of apolipoprotein E on the human immunodeficiency virus protein tat in neuronal cultures and synaptosomes. *J. of Neuroscience Research* **77**:532-539.

Cooksey, R.C., et al. (2004) Oxidative stress, beta cell apoptosis and decreased insulin secretory capacity in mouse models of hemochromatosis. *Endocrinology* **145**:5305-5312.

Kuzmenko, A.I. et al. (2004) Surfactant lipid peroxidation damages surfactant protein A and inhibits interactions with phospholipid vesicles. *J. Lipid Res.* 45:1061-1068.

Rezailki, L. et al. (2004). Respiration metabolism reduces oxidative and acid stress to improve long-term survival of Lactococcus lactis. *Molecular Microbiology* **53(5)**:1331-1342.

Palacino, J. et al. (2004). Mitochondrial Dysfunction and Oxidative Damage in parkin-deficient Mice. *J. of Biol. Chemistry* **279(18)**:18614-18922.

Ferri, A. et al. (2004). Activity of protein phosphatase calcineurin is decreased in sporadic and familial amyotrophic lateral sclerosispatients. *Journal Neurochemistry* **90 (5)**:1237.

Bakala, H., et al. (2003) Changes in rat liver mitochondria with aging. *Eur. J. Biochem.* 270:2295-2302.

Turchan, J., et al. (2003) Oxidative stress in HIV demented patients and protection ex vivo with novel antioxidants. *Neurology* **60**:307-314.

Graziewicz, M.A., et al. (2002) The mitochondrial DNA polymerase as a target of oxidative damage. *Nucleic Acids Research* **30**:2817-2824.

Lim, G.P., et al. (2001) The curry spice curcumin reduces oxidative damage and amyloid pathology in an Alzheimer transgenic mouse. *J Neuroscience* **21**:8370-8377.

Disclaimers

Tween[®] is a registered trademark of ICI Americas.

All trademarks, unless otherwise noted, are the property of Millipore.

Warranty

EMD Millipore Corporation ("EMD Millipore") warrants its products will meet their applicable published specifications when used in accordance with their applicable instructions for a period of one year from shipment of the products. **EMD MILLIPORE MAKES NO OTHER WARRANTY, EXPRESSED OR IMPLIED. THERE IS NO WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE.** The warranty provided herein and the data, specifications and descriptions of EMD Millipore products appearing in EMD Millipore's published catalogues and product literature may not be altered except by express written agreement signed by an officer of EMD Millipore. Representations, oral or written, which are inconsistent with this warranty or such publications are not authorized and if given, should not be relied upon.

In the event of a breach of the foregoing warranty, EMD Millipore Corporation's sole obligation shall be to repair or replace, at its option, the applicable product or part thereof, provided the customer notifies EMD Millipore Corporation promptly of any such breach. If after exercising reasonable efforts, EMD Millipore Corporation is unable to repair or replace the product or part, then EDM Millipore shall refund to the Company all monies paid for such applicable Product. **EMD MILLIPORE CORPORATION SHALL NOT BE LIABLE FOR CONSEQUENTIAL, INCIDENTAL, SPECIAL OR ANY OTHER DAMAGES RESULTING FROM ECONOMIC LOSS OR PROPERTY DAMAGE SUSTAINED BY ANY COMPANY CUSTOMER FROM THE USE OF ITS PRODUCTS.**

Unless otherwise stated in our catalog or other company documentation accompanying the product(s), our products are intended for research use only and are not to be used for any other purpose, which includes but is not limited to, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses or any type of consumption or application to humans or animals.

(c) 2009 – 2016: Merck KGaA, Darmstadt. All rights reserved. No part of these works may be reproduced in any form without permission in writing

This page left blank intentionally