





SIGMA QUALITY CONTROL TEST

## Product Information

### Enzymatic Assay of CHOLINE OXIDASE (EC 1.1.3.17)

#### PRINCIPLE:

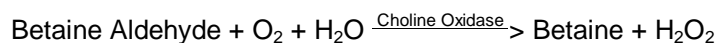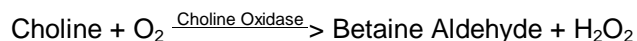

Abbreviation used:  
POD = Peroxidase

**CONDITIONS:** T = 37°C, pH = 8.0, A<sub>500nm</sub>, Light path = 1 cm

**METHOD:** Continuous Spectrophotometric Rate Determination

#### REAGENTS:

- A. 100 mM Tris HCl Buffer, pH 8.0 at 37°C  
(Prepare 100 ml in deionized water using Trizma Base, Sigma Prod. No. T-1503. Adjust to pH 8.0 at 37°C with 1 M HCl.)
- B. 2.1% (w/v) Choline Chloride Solution (Choline)  
(Prepare 100 ml in Reagent A using Choline, Chloride Salt, Sigma Prod. No. C-1879.)
- C. 1% (w/v) 4-Aminoantipyrine Solution (4-AAP)  
(Prepare 2 ml in deionized water using 4-Aminoantipyrine, Free Base, Sigma Prod. No. A-4382.)
- D. 1% (w/v) Phenol Solution (Phenol)  
(Prepare 5 ml in deionized water using Phenol, Sigma Prod. No. P-3653.)
- E. 10 mM Tris HCl with 2.0 mM Ethylenediaminetetraacetic Acid and 134 mM Potassium Chloride Solution (Enz Dil)  
(Prepare 100 ml in deionized water using Trizma Base, Sigma Prod. No. T-1503, Ethylenediaminetetraacetic Acid, Disodium Salt, Dihydrate, Sigma Stock No. ED2SS, and Potassium Chloride, Sigma Prod. No. P-4504. Adjust to pH 8.0 at 37°C with 1 M HCl.)

**Enzymatic Assay of CHOLINE OXIDASE  
(EC 1.1.3.17)**

**REAGENTS:** (continued)

- F. Peroxidase Enzyme (POD)  
(Use Peroxidase, Sigma Prod. No. P-8250.)
- G. Choline Oxidase Enzyme Solution (Choline Oxidase)  
(Immediately before use, prepare a solution containing 0.1 - 0.5 unit/ml of Choline Oxidase in cold Reagent E.)

**PROCEDURE:**

Prepare a reaction cocktail by pipetting (in milliliters) the following reagents into a suitable amber container:

Reagent B (Choline)	97.00
Reagent C (4-AAP)	1.00
Reagent D (Phenol)	2.00
Reagent F (POD, Purpurogallin units)	500

Mix by swirling.

Pipette (in milliliters) the following reagents into suitable cuvettes:

	<u>Test</u>	<u>Blank</u>
Reaction Cocktail	3.00	3.00

Equilibrate to 37°C. Monitor the  $A_{500\text{nm}}$  until constant, using a suitably thermostatted spectrophotometer. Then add:

Reagent E (Enz Dil)	-----	0.05
Reagent G (Choline Oxidase)	0.05	-----

Immediately mix by inversion and record the increase in  $A_{500\text{nm}}$  for approximately 5 minutes. Obtain the  $\Delta A_{500\text{nm}}/\text{minute}$  using the maximum linear rate for both the Test and Blank.

**CALCULATIONS:**

$$\text{Units/ml enzyme} = \frac{(\Delta A_{500\text{nm}}/\text{min Test} - \Delta A_{500\text{nm}}/\text{min Blank})(3.05)(\text{df})}{(12)(0.5)(0.05)}$$

3.05 = Volume (in milliliters) of assay

df = Dilution factor

12 = Millimolar extinction coefficient of Quinoneimine Dye at 500 nm under the conditions of the assay<sup>1</sup>

0.5 =  $\mu\text{mole}$  of Quinoneimine Dye formed per  $\mu\text{mole}$  of  $\text{H}_2\text{O}_2$

0.05 = Volume (in milliliter) of choline oxidase used in assay

**Enzymatic Assay of CHOLINE OXIDASE  
(EC 1.1.3.17)**

**CALCULATIONS:**

$$\text{Units/mg solid} = \frac{\text{units/ml enzyme}}{\text{mg solid/ml enzyme}}$$

$$\text{Units/mg protein} = \frac{\text{units/ml enzyme}}{\text{mg protein/ml enzyme}}$$

**UNIT DEFINITION:**

One unit will form 1.0  $\mu\text{mole}$  of  $\text{H}_2\text{O}_2$  from the oxidation of 1  $\mu\text{mole}$  of choline to betaine aldehyde per minute at pH 8.0 at 37°C.

**FINAL ASSAY CONCENTRATION:**

In a 3.05 ml reaction mix, the final concentrations are 96 mM Tris, 2.0% (w/v) choline, 0.01% (w/v) 4-aminoantipyrine, 0.02% (w/v) phenol, 15 units peroxidase, 0.03 mM ethylenediaminetetraacetic acid, 2 mM potassium chloride and 0.005 - 0.025 unit choline oxidase.

**REFERENCES:**

Okabe, H., Sagesaka, K., Nakajima, N., and Noma, A. (1977) *Clinica Chimica Acta* **80**, 87-94

Kessey, J. (1987) *Biochemica Information*, 1st ed., pp 19-20, Boehringer Mannheim Biochemicals, IN

**NOTES:**

1. The millimolar extinction coefficient is described in Kessey, J. (1982).
2. This assay is based on the cited references.
3. Peroxidase Unit Definition: One unit will form 1.0 mg purpurogallin from pyrogallol in 20 seconds at pH 6.0 at 20°C.
4. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that Sigma products conform to the information in Sigma publications. Purchaser must determine the suitability of the information and products for its particular use. Upon purchase of Sigma products, see reverse side of invoice or packing slip for additional terms and conditions of sale.