M Air T[™] Tester

A sieve-impaction sampler with ready-to-use agar cassettes



The M Air T is a lightweight, portable system for testing the microbial quality of air in clean rooms, barrier environments, and critical filling areas. The tester is small enough to be used in confined spaces, but powerful enough to sample up to 1000 liters in just 7 minutes. Programmable electronics and unique, ready-to-use agar cassettes make it the most convenient method for active air monitoring.

Fast

Sample up to 1000 liters (1 m^3) in less than 7 minutes.

Programmable

Reduce setup time by saving the air flow parameters in the tester's memory. You'll also reduce the risk of an operator selecting the wrong parameters. A delay timer allows the operator to leave the area before testing begins.

Convenient

Run up to 7 hours on a single charge. Depending on your usage, you may only have to re-charge the battery once a week.

No Overlapping Colonies

The sieve has nearly 1000 micro-perforations that evenly distribute air over the agar surface. Colony overlapping is minimized and results are easier to read.

Low Turbulence

The tester's small size and unique design will not disrupt the air currents inside a laminar hood.



For Use with Our Pre-filled Agar Cassettes

M Air T cassettes have a consistent filling level and flat surface. The distance between the sieve and the surface of the agar remains constant, which improves reproducibility. Other features of the cassettes, include:

- Double-sleeve packaging for barrier environments
- USP and EP compliance certificate
- Wings on the edge of the cassette for easy positioning on the tester head
- Integrated grid for easier colony counting
- The quantity of agar prevents desiccation of the medium and allows sampling up to 1000 liters (1 m³).
- Labels on each cassette for record-keeping and tracing
- Large choice of media
- Gamma-irradiated



Tester Specifications

Dimensions

255 mm (H); 105 mm (D); 125 mm (W) (10" x 4" x 5")

Weight (with batteries)

1.9 kg (4 lb)

Materials

Tester body: epoxy-coated aluminum Micro-perforated sampling sieve: 316 stainless steel (autoclavecompatible)

Sampling head: anodized aluminum

Sanitizing

Materials are compatible with most commonly-used sanitizing agents

Air flow rate

140 L/min for the first 500 liters, then 180 L/min (tester comes with a calibration certificate)

Regulatory compliance

Received CE mark

Ordering Information

Description	Qty/Pk	Catalogue No.
M Air T Tester		
M Air T Kit	1	ATAS05060
(Includes carrying case and the items below.		
Agar cassettes sold separately.)		
M Air T Tester	1	ATASPLRO 1
Battery charger, 50/60 Hz	1	ATAC05060
Micro-perforated sieve with cover	1	ATAHEADO 1
Allen-key	1	ATAKEYO01
Tripod	1	ATASUPO01

M Air T Agar Cassettes

Pre-filled, gamma-sterilized in double-sleeves (6 bags with 10 cassettes each)

Medium	Application	Qty/Pk	Catalogue No.
Tryptic Soy	Total count	60	ATSMTTD60
TSA with Pnase	Areas containing penicillin	60	ASTMPND60
Empty cassettes, EC) sterilized	114	ASTMC0114

To Place an Order or Receive Technical Assistance

For additional information call your nearest Millipore office:

In the U.S. and Canada, call toll-free 1-800-MILLIPORE

(1-800-645-5476) In the U.S., Canada and Puerto Rico, fax orders to 1-800-MILLIFX

(1-800-645-5439) On the Internet

http://www.millipore.com

To place an order call:

1-800-766-7000

E-mail: tech_service@millipore.com

Also Available From Fisher

Scientific in U.S. and Canada

Millipore Worldwide

AUSTRALIA

Tel. 1 800 222 111 or (02) 9888 8999 Fax (02) 9878 0788

AUSTRIA, CENTRAL EUROPE, C.I.S., AFRICA, MIDDLE EAST AND GULF Tel. (43) 1 877-8926

Tel. (43) 1 877-8926 Fax (43) 1 877-1654

BALTIC COUNTRIES

Tel. +358 9 804 5110 Fax +358 9 256 5660

BELGIUM AND LUXEMBOURG

Tel. (02) 726 88 40 Fax (02) 726 98 84

BRAZIL

Tel. (011) 548-7011 Fax (011) 548-7923

CANADA

Tel. 1-800-645-5476 Fax 1-800-645-5439

CHINA, PEOPLE'S

REPUBLIC OF Beijing:

Tel. (8610) 8519-1250 (8610) 8519-1058 Fax (8610) 8519-1255

Guangzhou: Tel. (8620) 8755-4021 Fax (8620) 8755-4350

Hong Kong: Tel. (852) 2803-9111 Fax (852) 2513-0313 Shanghai:

Tel. (8621) 5306-9100 Fax (8621) 5306-0838

CZECH REPUBLIC Tel. 02-2051 3841 02-2051 3842

02-2051 3842 Fax 02-2051 4298

DENMARK

Tel. 70 10 00 23 Fax 70 10 13 14

FINLANI

Tel. (09) 804 5110 Fax (09) 256 5660

FRANCE

Tel. (01) 30.12.70.00 Fax (01) 30.12.71.80

GERMANY

Tel. (06196) 494-0 Fax (06196) 43901

HUNGARY

Tel. 1-205 9784 Fax 1-205 9792

INDIA

Tel. (91) 80-839 46 57 Fax (91) 80-839 63 45

ITALY

Vimodrone (Milano): Tel. (02) 25.07.81 Fax (02) 26.50.324 Roma:

Tel. (06) 52.03.600 Fax (06) 52.95.735

JAPAN

Tel. (03) 5442-9711 Fax (03) 5442-9736 Analytical 9737 BioProcess 9734 Lab Water

KOREA

Tel. (822) 551-0990 Fax (822) 551-0228

MALAYSIA

Tel. 603-757 1322 Fax 603-757 1711

MEXICO

Tel. (525) 576-9688 Fax (525) 576-8706

THE NETHERLANDS Tel. 076-5022000

Fax 076-5022436

NORWAY

Tel. 22 67 82 53 Fax 22 66 04 60

POLAND

Tel. 22-669 12 25 22-663 70 31 Fax 22-663 70 33

PUERTO RICO

Tel. (787) 273-8495 Fax (787) 747-6553

SINGAPORE

Tel. (65) 842 1822 Fax (65) 842 4988

SPAIN AND PORTUGAL

Madrid: Tel. 917 283 960 Fax 917 292 909 Barcelona: Tel. 934 525 530

Fax 934 516 048

SWEDEN

Tel. 08-628 6960 Fax 08-628 6457

SWITZERLAND

Volketswil: Tel. (01) 908-30-60 Fax (01) 908-30-80 Lausanne:

Tel. (021) 641 2550 Fax (021) 641 2551

TAIWAN Taipei:

Tel. (886-2) 700-1742 Fax (886-2) 755-3267 Hsin Chu City: Tel. (886-3) 571-0178 Fax (886-3) 572-9520

U.K. AND IRELAND

Tel. (01923) 816375 Fax (01923) 818297

U.S.A.

Tel. (781) 533-6000 Fax (781) 533-3110

IN ALL OTHER COUNTRIES

Millipore Intertech (U.S.A.) Tel. +1 (781) 533-8622 Fax +1 (781) 533-8630

MILLIPORE

Millipore and M Air T are trademarks of Millipore Corporation or an affiliated company.

Lit. No. PF10176EN00 4/00 00-060

Printed in U.S.A.

Copyright © 2000 Millipore Corporation

or an affiliated company. All rights reserved.