

Data Sheet

BioTracker™ MB1 Tyrosinase Live Cell Dye

Live Cell Probe

SCT058**Pack Size 1 mg****Store at -20 °C****FOR RESEARCH USE ONLY****Not for use in diagnostic procedures. Not for human or animal consumption.**

Background

Melanoma is a form of skin cancer arising in melanocytes. Melanoma is an aggressive malignancy with rapid growth and early metastasis. Tyrosinase (TYR) is a potential biomarker for the detection of melanoma because its expression level is closely correlated with the level of malignancy.

BioTracker™ MB1 Tyrosinase live cell dye is a near-infrared dye that selectively reacts with TYR but not with various potential interfering agents, including inorganic salts (Na^+ , K^+ , Ca^{2+} , and Mg^{2+}), amino acids (Cys, Glu, and Arg), and other biological molecules including reduced GSH, BSA, glucose, hydrogen peroxide, vitamin C, lysozyme, glucose oxidase (GOD), and TYR.

Source

The BioTracker™ MB1 Tyrosinase Live Cell Dye (SCT058) does not contain genetically modified organisms.

Spectral Properties

Excitation max: 670 nm

Emission max: 690 nm

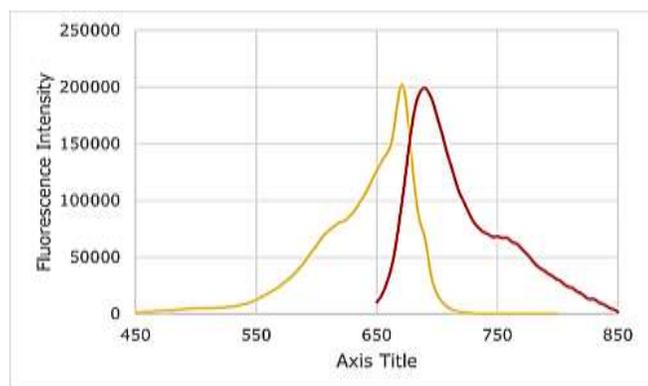


Figure 1: Probe excitation and emission data. 1 μL of probe at stock concentration (10 mM) was diluted in 1 mL of PBS buffer (20 mM, pH 6.5). Mushroom Tyrosinase (T3824) was added (10 U/mL) and incubated at 37 °C for 30 minutes to induce probe activation before undergoing excitation and emission scans. Spectral scans were conducted using a PerkinElmer FL8500 Fluorescence Spectrophotometer.

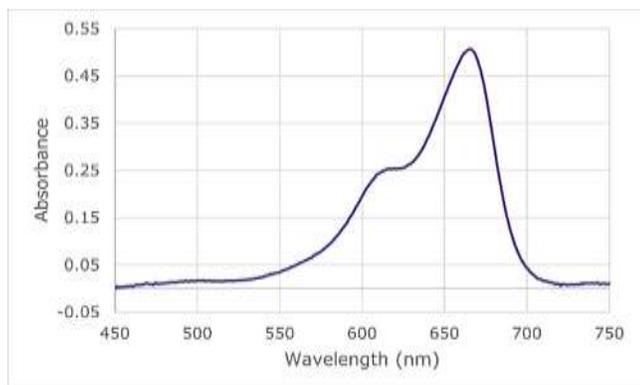


Figure 2: Probe absorbance data. 1 μL of probe at stock concentration (10 mM) was diluted in 1 mL of PBS buffer (20 mM, pH 6.5). Mushroom Tyrosinase (T3824) was added (10 U/mL) and incubated at 37 $^{\circ}\text{C}$ for 30 minutes to induce probe activation before undergoing absorbance scan. Spectral scans were conducted using a PerkinElmer FL8500 Fluorescence Spectrophotometer.

Quality Control Testing

Purity: $\geq 98\%$ confirmed by HPLC. Structure confirmed by HNMR, LC-MS and elemental analysis.

Storage and Handling

Store BioTracker™ MB1 Tyrosinase Live Cell Dye at $-20\text{ }^{\circ}\text{C}$, desiccated and protected from light.

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
Note: Centrifuge vial briefly to collect contents at bottom of vial before opening.

Presentation


Lyophilized

Representative Data

A. Control



B. +MB1

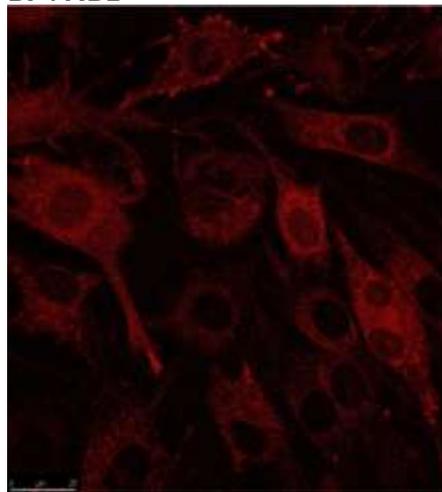

C. +MB1/+Kojic

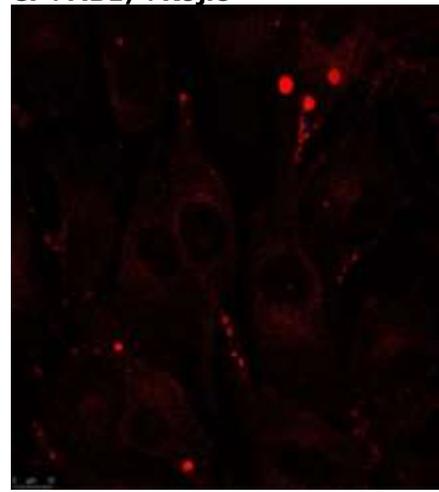


Figure 3: (B.) Imaging of MB1 Tyrosinase live cell probe in SK-MEL-28 cells. Cells were treated with 20 μM MB1 for 4 hours at 37 $^{\circ}\text{C}$. (C.) SK-MEL-28 cells were pre-treated with 1 mM Kojic acid (TYR inhibitor, K3125) for 1 hour before being treated with 20 μM MB1. (A.) Control cells were left unstained.

Protocols

Reagent Preparation

1. Before opening the vial, spin down the solid to the bottom by a microcentrifuge or by a desktop centrifuge.
2. Warm the vial to room temperature and prepare the MB1 probe (MW 448.58 g/mol) stock solution by dissolving the contents of one vial (1 mg) in 223 μ L DMSO to create a 10 mM solution (freeze aliquots at -20 °C).
3. Dilute in cell culture media to a final concentration of 10-20 μ M and add to cells in culture. Incubate at 37 °C for 4 hours.
4. Wash cells with PBS buffer before imaging.

Note: Optimal concentration must be determined by end user.

References

1. Li Z, Wang Y, Zeng C-C, Hu L, Liang X-J. 2018. Ultrasensitive Tyrosinase-Activated Turn-On Near-Infrared Fluorescent Probe with a Rationally Designed Urea Bond for Selective Imaging and Photodamage to Melanoma Cells. 90(6):3666–3669.

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates
as MilliporeSigma in the U.S. and Canada.

Merck BioTracker and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2024 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

Document Template 20306518 Ver 6.0

23062841 Ver 1.0, Rev 13MAY2024, AV

The Merck logo is displayed in a bold, red, sans-serif font.