## Safety Sheet

# DirectLoad™ Midi or Mini

## Horizontal Electrophoresis System

### DMIDI

### DMINI

|                                    |    |                                   |    |
|------------------------------------|----|-----------------------------------|----|
| English . . . . .                  | 2  | 한국어 (Korean) . . . . .            | 30 |
| Български (Bulgarian) . . . . .    | 4  | Latviski (Latvian) . . . . .      | 32 |
| 中文 (Chinese, Simplified) . . . . . | 6  | Lietuvių (Lithuanian) . . . . .   | 34 |
| Hrvatski (Croatian) . . . . .      | 8  | Norsk (Norwegian) . . . . .       | 36 |
| Čeština (Czech) . . . . .          | 10 | Polski (Polish) . . . . .         | 38 |
| Dansk (Danish) . . . . .           | 12 | Português (Portuguese) . . . . .  | 40 |
| Nederlands (Dutch) . . . . .       | 14 | Română (Romanian) . . . . .       | 42 |
| Eesti keel (Estonian) . . . . .    | 16 | Srpski (Serbian) . . . . .        | 44 |
| Français (French) . . . . .        | 18 | Slovenčina (Slovak) . . . . .     | 46 |
| Deutsch (German) . . . . .         | 20 | Slovenščina (Slovenian) . . . . . | 48 |
| Ελληνικά (Greek) . . . . .         | 22 | Español (Spanish) . . . . .       | 50 |
| Magyar (Hungarian) . . . . .       | 24 | Svenska (Swedish) . . . . .       | 52 |
| Italiano (Italian) . . . . .       | 26 | Türkçe (Turkish) . . . . .        | 54 |
| 日本語 (Japanese) . . . . .           | 28 |                                   |    |

## Intended Use

The DirectLoad™ Horizontal Electrophoresis System is used to perform high-resolution nucleic acid separation on agarose gels.

## Specifications

### Device

|               |                        |
|---------------|------------------------|
| Rated Voltage | Maximum 250 Vdc (±10%) |
| Rated Power   | Maximum 250 mA         |

### DMIDI Midi System

|            |        |         |              |
|------------|--------|---------|--------------|
| Dimensions | Height | 180 mm  | 7.0 inches   |
|            | Width  | 265 mm  | 10.25 inches |
|            | Depth  | 75 mm   | 3.0 inches   |
| Weight     |        | 1.65 kg | 3.64 lbs     |

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|            |        |         |            |
|------------|--------|---------|------------|
| Dimensions | Height | 100 mm  | 4.0 inches |
|            | Width  | 225 mm  | 9.0 inches |
|            | Depth  | 75 mm   | 3.0 inches |
| Weight     |        | 0.85 kg | 1.25 lbs   |

## Environmental Conditions

For indoor use only. This equipment is not rated for wet locations or any other outdoor parameter.

### Storage and Operating

|             |                           |
|-------------|---------------------------|
| Temperature | 10-65 °C                  |
| Humidity    | 20 to 80%, non-condensing |

Atmospheric Pressure Below 2000 M

This product is rated POLLUTION DEGREE 2 in accordance with IEC 664.

## User Guide

User guides and other documents are available online. Go to [sigmaaldrich.com](https://sigmaaldrich.com), enter the catalogue number in the search bar, select your product, and scroll to the documents section.

Accessories and related products (if applicable) can be ordered online at [sigmaaldrich.com](https://sigmaaldrich.com).

## Precautions

Please read these safety instructions carefully before using the device and/or any accessory supplied with this product. Keep a printed copy of the User Guide near the system for reference.

### General Warnings

- This equipment has been made as safe as possible for normal use. However, we do not guarantee safety if the equipment is used in a manner not specified.
- Any personnel from your organization who will use the system must be properly trained to ensure safe and effective use.
- Do not attempt to make modifications to the equipment. Contact us at [sigmaaldrich.com/techservice](https://sigmaaldrich.com/techservice) prior to repairs.
- Only use original manufacturer replacement parts. See the User Guide for catalogue numbers.
- Unauthorized changes to the system can endanger persons and damage the system and property of the company operating it.
- Check that there are no alarms or fault conditions before use.
- This equipment should only be used as specified as stated in the "Intended Use" section and described in the User Guide.
- The cover to this product only has a 1 joule of energy impact rating which can crack if dropped or exposed to falling objects. Please use caution handling this part to avoid a hazardous situation during operation and use.
- Do not attempt to use the tank without the safety lid properly in place.

### Electrical Warnings

- Do not touch plug or power supply with wet hands. This may result in injury, fire and electric shock.

### Usage Guidance and Restrictions

When used incorrectly, this product can deliver dangerous levels of electricity and are to be operated only by qualified personnel following the instructions laid out in the user guide. Anyone intending to use this equipment should read the complete manual thoroughly. The product must never be used without the safety lid correctly in position. The product should not be used if there is any sign of damage to the external tank or lid.

## Technical Service

For assistance, go to [sigmaaldrich.com/techservice](https://sigmaaldrich.com/techservice). Unauthorized repair or service to the device may void the warranty and/or cause hazardous conditions.

## Maintenance

- Unplug the instrument before beginning cleaning, to prevent risk of electrical shock.
- Use warm water and a mild detergent to clean. Water at temperatures above 60 °C can cause damage to the product and components.
- Compatible detergents include dish washing liquid, Hexane and Aliphatic hydrocarbons
- The tank should be thoroughly rinsed with warm water or distilled water to prevent build-up of salts, but care should be taken not to damage the enclosed electrode and vigorous cleaning is not necessary or advised.
- Do not submerge or soak in detergents for more than 30 minutes.
- Allow to air dry before use.
- The following cleaning agents will cause irreversible and accumulative damage, and should never be used:

|                      |                   |
|----------------------|-------------------|
| Acetone              | Methanol          |
| Phenol               | Ethanol           |
| Chloroform           | Isopropyl alcohol |
| Carbon tetrachloride | Alkalis           |
- RNAse Decontamination can be performed using the following protocol:
  1. Clean the units with a mild detergent as described above.
  2. Wash with 3% hydrogen peroxide (H<sub>2</sub>O<sub>2</sub>) for 10 minutes.
  3. Rinsed with 0.1% DEPC-(diethyl pyro carbonate) treated distilled water.

**Caution:** DEPC is a suspected carcinogen. Always take precautions when using. RNaseZAP™ cleaning agent can also be used. Please consult the instructions for use with acrylic gel tanks.

## Equipment Operation

Only operate on a level and stable surface.

### POWER ON

- Electrode cables should be inspected for defects prior to use.
- Connect the provided positive (red) and negative (black) cable from the tank lid to a certified power supply via the power supply's input terminals respectfully.

### POWER OFF

- Shut down any output being generated from the separately certified power supply. Once output power has ceased, remove the electrode cable connector from the power supply.
- Ensure the power is off prior to removing the lid.

## Symbol Definitions

|                                                                                   |                                                                |
|-----------------------------------------------------------------------------------|----------------------------------------------------------------|
|  | Consult this Safety Sheet for important cautionary information |
|  | Voltage or Shock Hazard                                        |
|  | Catalogue Number                                               |
|  | Serial Number                                                  |
|  | Manufacturer                                                   |

Declarations of Conformity are available on the product page at [sigmaaldrich.com](https://sigmaaldrich.com).

European Council Directives conformity marking.

United Kingdom conformity marking

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

## Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at [sigmaaldrich.com/terms](https://sigmaaldrich.com/terms).

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Made in the United Kingdom

EMD Millipore Corporation,  
Burlington, MA, U.S.A.,  
an affiliate of Merck KGaA,  
Darmstadt, Germany

UK Responsible Person:  
Millipore (UK) Ltd.  
Kirkton Campus  
Fleming Road,  
Livingston, UK  
EH 7BN

The life science business of Merck KGaA, Darmstadt, Germany  
operates as MilliporeSigma in the U.S. and Canada.

The initial M, Millipore, DirectLoad, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.  
© 2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.