

Product Information

DERMATO SUPPLEMENT

Product Number **D5185**

Product Description

An antibiotic supplement recommended for the selective isolation of pathogenic dermatophytes.

Components

Item

Cycloheximide	250.00 mg
Chlortetracycline	50.00 mg
Gentamicin	50.00 mg

Precautions and Disclaimer

For laboratory use only. Not for drug, household or other uses.

Preparation Instructions

Rehydrate the contents of one vial aseptically with 5 mls of 50% (v/v) acetone. Mix well and aseptically add to 500 mls of sterile, molten (D8059)D.T.M. Agar Base. Mix well and pour into sterile petri plates.

Storage

Store the supplement at 2-8°C.

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.