

Product Information

REDEExtract-N-Amp™ Seed PCR Kit

XNASS, XNAS**Product Description**

The REDEExtract-N-Amp™ Seed PCR Kit for direct PCR contains the reagents needed to rapidly extract and amplify genomic DNA from seeds (soybean, corn, wheat, etc.). Briefly, DNA is extracted from ground seed material by incubation in a mixture of Extraction Solution and Seed Preparation Solution at 55 °C for 10 minutes. There is no need for organic extraction, column purification, or precipitation of the DNA. After the extraction is stopped by incubation at 95 °C for 3 minutes, an equal volume of Neutralization Solution B is added, and the extract is ready for PCR.

An aliquot of the neutralized extract is then combined with the REDEExtract-N-Amp™ PCR Reaction Mix and user-provided PCR primers to amplify target DNA. The REDEExtract-N-Amp™ PCR Reaction Mix is a 2X ready mix containing buffer, salts, dNTPs, and *Taq* DNA polymerase. It is optimized specifically for use with the extraction reagents. This formulation also contains the JumpStart™ antibody for specific hot start amplification, and RED*Taq*® dye to allow direct loading of the PCR product onto an agarose gel.

Reagents Provided		XNASS	XNAS
	Cat. No.	10 Preps	100 Preps
Extraction Solution	E 7526	1.2 mL	6 mL
Seed Preparation Solution	S 1193	0.15 mL	0.9 mL
Neutralization Solution B	N 3910	1.2 mL	6 mL
REDEExtract-N-Amp™ PCR Reaction Mix, This is a 2X PCR mix containing buffer, salts, dNTPs, <i>Taq</i> DNA polymerase, RED <i>Taq</i> ® dye, and JumpStart™ antibody.	R 4775	0.15 mL	1.2 mL

Reagents and Equipment Required (Not Provided)

Items common to all procedures:

- Tubes or plates for PCR
- PCR primers, Cat. No. OLIGO
- Thermal cycler
- Water, PCR grade, Cat. No. W1754

For individual 1.5 mL tubes:

- 1.5 mL microcentrifuge tubes
- Heat block or thermal cycler
- Disposable plastic pestles

For individual 1.5 mL tubes with liquid nitrogen:

- 1.5 mL microcentrifuge tubes
- Liquid nitrogen
- Mortar and pestle
- Heat block or thermal cycler

For 96-well plates:

- Bead Mill
- 4 mm stainless steel grinding balls
- 2 mL square well block
- 96-well sealing mat
- 96-well PCR plate
- Thermal cycler
- Optional: Heat block with 96-well block

Precautions and Disclaimer

This product is for R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Storage

All components of the REDExtract-N-Amp™ PCR Kit can be stored at 2-8 °C for up to 3 weeks. For long-term storage, greater than 3 weeks, -20 °C is recommended. Do not store in a "frost-free" freezer.

Procedure

All steps are carried out at room temperature unless otherwise noted.

Grinding Seeds

Following are three different methods for grinding seeds.

Grind using a bead mill

1. Place 1 seed into each well of a 2 mL square well block.
Note: With *Arabidopsis* or similar sized seeds, approximately 50 seeds should be placed in a single well.
2. This grinding procedure is not recommended for corn seeds, because results with such large, tough seeds are inconsistent.

3. Pipette PCR grade water into the well according to the following volumes:

- 800 µL for soybean or similar sized seeds
- 600 µL for cotton or similar sized seeds
- 200 µL for canola, sorghum, wheat, or similar sized seeds
- 100 µL for *Arabidopsis* or similar sized seeds

4. Place a 4 mm stainless steel grinding ball in each well of the 2 mL 96-square well block and cover with sealing mat. Place block in the bead mill and shake at 1,500 rpm for 10 minutes. Continue to [Extraction of Seeds](#).

Grind individually using a plastic pestle

1. Place 1 seed into a 1.5 mL microcentrifuge tube.
Note: With *Arabidopsis* or similar sized seeds approximately 50 seeds should be placed in a single tube.
2. Pipette PCR grade water into the well according to the following volumes:
 - 800 µL for soybean or similar sized seeds
 - 600 µL for corn or similar sized seeds
 - 400 µL for cotton or similar sized seeds
 - 100 µL for *Arabidopsis*, canola, sorghum, wheat, or similar sized seeds
3. Incubate the seed with water for 1 hour at 55 °C.
4. Grind hydrated seeds in tube using a plastic pestle. Continue to [Extraction of Seeds](#).

Grind individually using liquid nitrogen

1. Grind seed into a fine powder in liquid nitrogen using a mortar and pestle.
Note: With small seeds, such as *Arabidopsis* and canola, more than one seed must be ground to collect enough ground seed material.
2. Transfer between 5 and 100 mg of ground seed material into a pre-weighed 1.5 mL microcentrifuge tube. Record mass of transferred seed material.
3. Pipette 4 µL of water for every mg of transferred ground seed material into the sample tube and vortex to mix. Continue to [Extraction of Seeds](#).

Extraction of Seeds

1. Pipette 45 µL of Extraction Solution into a 1.5 mL microcentrifuge tube or multi-well PCR plate. Add 5 µL of Seed Preparation Solution to the tube and pipette up and down to mix.

Note: If several extractions will be performed, sufficient volumes of Extraction and Tissue Preparation Solutions may be pre-mixed in a ratio of 9:1 up to 2 hours before use. The mixture should then be dispensed in 50 μ L volumes into tubes or multi-well plates.

- Pipette 5 μ L of the ground seed suspension from [Grinding Seeds](#) section into the Extraction Solution and Seed Preparation Solution mixture and vortex or pipette up and down to mix.
- Incubate the mixture at 55 °C for 10 minutes to extract DNA. Note that the ground seed will not appear to be digested at the end of this incubation; however, sufficient DNA will be released for PCR.
- Incubate the mixture at 95 °C for 3 minutes to stop the extraction.

Note: Steps 3 and 4 can be performed in a thermal cycler if using a 96-well PCR plate.

- Add 50 μ L of Neutralization Solution B to the mixture and vortex or pipette up and down to mix.
- Store the neutralized seed extract at 2-8 °C or continue to [PCR amplification](#) section.

PCR amplification

The REDExtract-N-Amp™ PCR Reaction Mix contains JumpStart™ antibody for specific hot start amplification. Therefore, PCR mixtures can be assembled at room temperature without premature *Taq* DNA polymerase activity.

Typical final primer concentrations are \sim 0.4 μ M each. The optimal primer concentration and cycling parameters will depend on the system being used.

- Add the following reagents to a thin-walled PCR microcentrifuge tube or plate:

Reagent	Volume
Water, PCR grade	Variable
REDExtract-N-Amp™ PCR reaction mix	10 μ L
Forward primer	Variable
Reverse primer	Variable
Seed extract	4 μ L*
Total volume	20 μ L

***Note:** The REDExtract-N-Amp™ PCR Reaction Mix is formulated to compensate for components in the Extraction, Seed Preparation, and Neutralization B Solutions. If less than 4 μ L of seed extract is added to the PCR reaction volume, use a 50:50 mixture of Extraction and Neutralization B Solutions to bring the volume of seed extract up to 4 μ L.

- Mix gently and briefly centrifuge to collect all the components to the bottom of the tube.
- For thermal cyclers without heated lids, add 20 μ L of mineral oil to the top of each tube to prevent evaporation.
- The amplification parameters should be optimized for individual primers, template, and thermal cycler.

Common cycling parameters:

Step	Temperature	Time	Cycles
Initial Denaturation	94 °C	3 minutes	1
Denaturation	94 °C	30 seconds	
Annealing	45 to 68 °C	30 seconds	30-35
Extension	72 °C	1-2 minutes (1 min/kb)	
Final Extension	72 °C	10 minutes	1
Hold	4 °C	Indefinitely	

- The amplified DNA can be loaded directly onto an agarose gel after the PCR is completed. It is not necessary to add a separate loading buffer/tracking dye.

Note: PCR products can be purified, if desired, for downstream applications such as sequencing with the GenElute™ PCR Clean-Up Kit, Cat. No. NA1020.

Troubleshooting Guide

Problem	Cause	Solution
Little or no PCR product is detected.	Seeds were not ground sufficiently.	For 96-well grind - Increase grinding time in bead mill. For seeds with a tough seed coat, it is helpful to break the seed before putting it into the plate to grind. For individual 1.5 mL tube - For seeds with a tough seed coat, it is helpful to break the seed before incubating it at 55 °C.
	PCR may be inhibited due to contaminants in the seed extract.	Dilute the extract with a 50:50 mixture of Extraction and Neutralization B Solutions. To test for inhibition, include a DNA control and/or add a known amount of template (100-500 copies) into the PCR along with the seed extract.
	A PCR component may be missing or degraded.	Run a positive control to ensure that components are functioning. A checklist is also recommended when assembling reactions.
	There may be too few cycles performed.	Increase the number of cycles (5-10 additional cycles at a time).
	The annealing temperature may be too high.	Decrease the annealing temperature in 2-4 °C increments.
	The primers may not be designed optimally.	Confirm the accuracy of the sequence information. If the primers are less than 22 nucleotides long, try to lengthen the primer to 25-30 nucleotides. If the primer has a GC content of less than 45%, try to redesign the primer with a GC content of 45-60%.
	The extension time may be too short.	Increase the extension time in 1-minute increments, especially for long templates.
	Target template is difficult.	In most cases, inherently difficult targets are due to unusually high GC content and/or secondary structure. Betaine (Cat. No. B0300) has been reported to help amplification of high GC content templates at a concentration of 1.0-1.7 M.
	JumpStart™ antibody is not working correctly.	Do not use DMSO or formamide with REDExtract-N-Amp™ PCR Reaction Mix. It can interfere with the enzyme-antibody complex. Other cosolvents, solutes (e.g., salts), extremes in pH, or other reaction conditions may reduce the affinity of the JumpStart™ antibody for Taq polymerase and thereby compromise its effectiveness.
	"Touchdown" PCR may be needed.	"Touchdown" PCR significantly improves the specificity of many PCR reactions in various applications. Touchdown PCR involves using an annealing/extension temperature that is higher than the Tm of the primers during the initial PCR cycles. The annealing/extension temperature is then reduced to the primer Tm for the remaining PCR cycles. The change can be performed in a single step or in increments over several cycles.
Negative control shows a PCR product or "false positive" results are obtained.	Reagents are contaminated.	Include a reagent blank without DNA template be included as a control in every PCR run to determine if the reagents used in extraction or PCR are contaminated with a template from a previous reaction.

References

1. Dieffenbach, C.W. and Dveksler, G.S. (Eds.) PCR Primer: A Laboratory Manual, 2nd ed., (Cold Spring Harbor Laboratory Press, New York, 2003).
2. Don, R.H., et al., 'Touchdown' PCR to circumvent spurious priming during gene amplification. *Nucleic Acids Res.*, 19, 4008 (1991).
3. Erlich, H.A. (Ed.) PCR Technology: Principles and Applications for DNA Amplification (Stockton Press, New York, 1989).
4. Griffin, H.G. and Griffin, A.M. (Eds.) PCR Technology: Current Innovations (CRC Press, Boca Raton, FL, 1994).
5. Innis, M.A., et al. (Eds.) PCR Strategies, Academic Press, New York, 1995.
6. Innis, M., et al. (Eds.) PCR Protocols: A Guide to Methods and Applications (Academic Press, San Diego, CA, 1990).
7. McPherson, M.J. et al. (Eds.) PCR 2: A Practical Approach (IRL Press, New York, 1995).
8. Newton, C.R. (Ed.) PCR: Essential Data (John Wiley & Sons, NY, 1995).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, REDExtract-N-Amp, REDTaq, Jumpstart and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

XNASpis Rev 09/22

Millipore
Sigma