# Sigma-Aldrich

#### Product Information

## 1-Fluoro-2,4-dinitrobenzene

≥ 98.5% (GC)

#### D1529

## **Product Description**

Molecular Formula: C<sub>6</sub>H<sub>3</sub>FN<sub>2</sub>O<sub>4</sub>

Molecular Weight: 186.1

CAS Registry Number: 70-34-8

Melting point: 26 °C¹ Density: 1.48 g/mL

Refractive Index: 1.5690 (20 °C)<sup>2</sup>

Molarity: 7.95 M

Synonyms: DNFB, 2,4-dinitrofluorobenzene, Sanger's Reagent, 2,4-dinitrobenzene fluoride, 2,4-DNFB, 1,2,4-Fluorodinitrobenzene, Fluoro-1,3-dinitrobenzene

1-Fluoro-2,4-dinitrobenzene (DNFB) is a halogenated and nitrated aromatic compound that modifies the *N*-terminal amino acids of proteins and peptides. DNFB is also commonly known as Sanger's Reagent, named after Frederick Sanger, who reported, in a classic 1945 paper, the use of DNFB for derivatization of proteins and peptides specifically at their *N*-termini.<sup>3</sup> In the Sanger procedure, DNFB is used to form dinitrophenyl derivatives,<sup>4</sup> for end group determination of proteins,<sup>3</sup> and to derivatize primary amines.<sup>5</sup>

The effect of DFNB on the enzymatic properties of mitochondrial b-c1 complex isolated from beef heart mitochondria indicated that the chemical modification by DNFB strongly inhibits the reductase activity of the complex.<sup>6</sup> DNFB has been reported to inactivate fructose diphosphatase by reacting with a cysteine residue.<sup>7</sup> DNFB can also cause reversible dinitrophenylation of thiols in muscle glyceraldehyde 3-phosphate dehydrogenase.<sup>8</sup> DNFB has been used as a sensitization reagent in contact hyperactivity studies.<sup>9</sup>

Several dissertations<sup>11-15</sup> have cited use of product D1529 in their protocols.

## Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

## Storage/Stability

The half-life of DNFB at 15 °C in a Tris buffer is:10

- 56.7 hours at pH 8.03
- 6.5 hours at pH 9.53

## **Preparation Instructions**

DNFB is soluble in chloroform (100 mg/mL). DNFB is also soluble in benzene, ether, and propylene glycol.

A saturated solution at 15 °C in water is 0.0086 M and in 8 M urea is 0.062 M.<sup>10</sup> The concentration of DNFB at saturation was determined spectrophotometrically as dinitrophenol after basic hydrolysis.

## References

1

- 1. The Merck Index, 12th ed., Entry# 4210 (1996).
- 2. Dean, J.A, (ed.), *Lange's Handbook of Chemistry*, 13th edition. McGraw-Hill (New York), p. 7-362 (1985).
- 3. Sanger, F., Biochem. J., **39(5)**, 507-515 (1945).
- 4. Greenstein, J.P., and Winitz, M., *Chemistry of the Amino Acids*, Vol. 2. Wiley (New York, NY), pp. 1552-1566 (1961).
- Fariss, M.W., and Reed, D.J., Methods Enzymol., 143, 101-109 (1987).
- Lorusso, M. et al., FEBS Lett., 195(1-2), 298-302 (1986).
- Pontremoli, S. et al., J. Biol. Chem., 240(9), 3464-3468 (1965).



- Shaltiel, S., and Soria, M., Biochemistry, 8(11), 4411-4415 (1969).
- Gimblet, C. et al., Cell Host Microbe, 22(1), 13-24.e4 (2017).
- 10. Murdock, A.L. *et al.*, *Arch. Biochem. Biophys.*, **114(2)**, 375-390 (1966).
- 11. Smookler, David, "The Role of TIMP3 in Models of Inflammation and Immunity". University of Toronto, Ph.D. dissertation, p. 57 (2010).
- 12. Cornely, Rhea, "The role of Annexin A6 in T cell activation". University of New South Wales, Ph.D. dissertation, p. 36 (2013).
- 13. Patra, VijayKumar, "Insight into the role of skin microbiome in UV-induced immune modulation and its potential link to pathogenesis of polymorphic light eruption". Medical University of Graz, Ph.D. dissertation, p. 33 (2018).
- 14. Gendrisch, Fabian, "The Role of Endoplasmic Reticulum Stress Responses in Contact Dermatitis". Albert-Ludwigs-Universität Freiburg im Breisgau, Ph.D. dissertation p. 36 (2019).
- Palmer, Brian C., "The Dermal Toxicity and Immunomodulatory Effects of Topically Applied Engineered Nanomaterials in a Model of Allergic Contact Dermatitis". University of Rochester, Ph.D. dissertation, p. 87 (2019).

### Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

#### Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

#### Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at <u>SigmaAldrich.com/terms</u>.

#### Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

