

Product Information

BCIP®/NBT Liquid Substrate System

Ready-to-use solution

B1911

Product Description

Storage temperature: 2-8 °C

The 5-bromo-4-chloro-3-indolyl phosphate (BCIP®)/nitro blue tetrazolium (NBT) Liquid Substrate System is a substrate system that is widely used with alkaline phosphatase conjugates.¹⁻³ This system produces a blue-purple product. The intense color can be observed visually, is very stable, and will not fade upon exposure to light.

The BCIP®/NBT Liquid Substrate System offers a convenient, ready-to-use substrate solution for visualizing alkaline phosphatase activity in immunoblotting and immunohistology. Immunohistochemical staining with BCIP®/NBT requires organic mounting media and can be counterstained with Nuclear Fast Red or Light Green. BCIP®/NBT is generally a more sensitive method than Fast Red. This product is not recommended for ELISA (multiwell) procedures.

Several publications,⁴ theses^{5,6} and dissertations⁷⁻²¹ have cited use of product B1911 in their protocols.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

The BCIP®/NBT Liquid Substrate System is a single solution, ready-to-use reagent.

Storage/Stability

Store the product at 2-8 °C for extended, long-term storage. Protect from light or UV sources.

Do not freeze. The product may be stored at room temperature for up to one year.

Usage

- Rinse specimens that have been incubated with an alkaline phosphatase conjugate in a non-phosphate wash buffer before treatment with the BCIP®/NBT Liquid Substrate System.
- Cover the entire specimen with the reagent during color development.
- Incubate the specimen at room temperature with the BCIP®/NBT reagent for ~10 minutes. Specimens and procedure may affect the length of time needed for color development.
- Monitor color development to avoid over-development.
- Stop color development by rinsing the specimen with water.

References

1. Blake, M.S. *et al.*, *Anal. Biochem.*, **136(1)**, 175-179 (1984).
2. Horowitz, J.P. *et al.*, *J. Med. Chem.*, **9(3)**, 447 (1966).
3. Smejkal, G.B., and Kaul, C.A., *J. Histochem. Cytochem.*, **49(9)**, 1189-1190 (2001).
4. Cabaleiro, C. *et al.*, *S. Afr. J. Enol. Vitic.*, **42(2)**, 165-174 (2021).
5. Unda, Faride, "Direct organogenesis and *Agrobacterium*-mediated transformation of *Exacum* styer group". University of British Columbia, M.Sc. thesis, p. 60 (2006).
6. Jang, Jungho, "Bioelectrochemical conversion of CO₂ to value added product formate using engineered *Methylobacterium extorquens*". Ulsan National Institute of Science and Technology, M.Sc. thesis, p. 27 (2018).

7. Tare, Rahul Shrikant, "Osteoblast-Stimulating Factor-1 (OSF-1): a peptide with diverse roles in bone development". University of Southampton, Ph.D. dissertation, p. 127 (2002).
8. Haupt, Borris, "Herpes simplex virus-1 (HSV-1) as a gene delivery vector for neural precursor cells". University College London, Ph.D. dissertation, p. 255 (2004).
9. van Boxtel, Evelien L., "Protein quaternary structure and aggregation in relation to allergenicity". Wageningen University, Ph.D. dissertation, pp. 91, 110 (2007).
10. Gunawardena, Harsha, "A clinical and serological study of adult and juvenile idiopathic inflammatory myopathy". University of Bath, Ph.D. dissertation, p. 61 (2010).
11. Kulkarni, Onkar P., "Role of therapeutic blockade of CCL2 in a mouse model of SLE and lupus nephritis". Ludwig-Maximilians-Universität München, Ph.D. dissertation, p. 50 (2010).
12. Senthilan, Pingkalai, "Identification and characterization of deafness genes in *Drosophila melanogaster*". Georg-August-Universität Göttingen, Dr. rer. nat. dissertation, p. 41 (2010).
13. Vance, Steven J., "The relationship between structure and function in natural surfactant proteins". University of Glasgow, Ph.D. dissertation, p. 246 (2012).
14. Mani, Jiju, "CD8⁺ T cell epitopes derived from the polyomavirus capsid protein VP1". University of Rostock, Dr. rer. nat. dissertation, p. 26 (2014).
15. Lercher, Lukas, "Chemical tools for the study of epigenetic mechanisms". University of Oxford, Ph.D. dissertation, p. 101 (2014).
16. Miri, Kamella, "The Imprinted Polycomb Group Gene *Sfmbt2* is Required for Trophoblast Maintenance and Placenta Development". University of Toronto, Ph.D. dissertation, p. 49 (2014).
17. Zhao, Qianru, "Regulation of starch biosynthesis in *Arabidopsis thaliana*". University of Guelph, Ph.D. dissertation, p. 69 (2015).
18. Eisel, David, "Reprogramming of M2-like macrophages by CD4⁺ T cells, transcription factor knockdown and miRNA transfection". Ruperto-Carola University Heidelberg, Dr. rer. nat. dissertation, p. 58 (2018).
19. Riera, Natalia Galindo, "Characterisation of a mouse model of Protocadherin 19". Cardiff University, Ph.D. dissertation, p. 58 (2019).
20. Sridhar, Krishna Chander, "Characterization of signaling events following single-cell death in myocardial cell clusters *in vitro*". Rheinischen Friedrich-Wilhelms-Universität Bonn, Dr. rer. nat. dissertation, p. 27 (2021).
21. Teixeira, Alyne Grasiela, "Development of microscale immunoassays using aqueous two-phase systems". Dalhousie University, Ph.D. dissertation, p. 158 (2021).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, BCIP, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2009-2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

B1911dat Rev 08/22 CS, RBG, GCY, MAM

**MILLIPORE
SIGMA**