English
Non-sterile Millex® 25 mm Filter Unit
Millex®-OR, Cat. No. SFGLO2500 (0,22 µm, 1,000/box)
Millex®-GS, Cat. No. SLGS02510 (0,22 µm, 1,000/box)
Millex®-HA, Cat. No. SLHA02510 (0,45 µm, 1,000/box)
● Single use only ● Non-sterile ● Non-pyrogenic

Intended Use
SLGS02510 and SFGLO2500 are intended for clarifying small volumes of aqueous solutions, and once sterilized, can be used as a sterilizing filter for aqueous solutions. SLHA02510 is intended for clarifying small volumes of aqueous solutions.

Directions for Use

WARNINGS:


- Do not use this product as an in-line filter for intravenous administration; it was not designed for long-term continuous use.
- Do not use with syringes smaller than 10 mL because pressures in excess of the maximum pressure rating may be reached, potentially causing damage to the filter unit and/or personal injury.

CAUTIONS:

- Do not use the Millex® filter unit at temperatures above 45 °C (113 °F).
- Do not use the same Millex® filter unit to filter solutions in both directions.
- Do not use the Millex® filter unit to filter emulsions or suspensions.
- Use caution when filtering solutions with very low concentrations of critical care drugs.
- Do not reuse the Millex® filter unit, as Merck Millipore Ltd. cannot assure performance and integrity beyond a single use.

Sterilization Guidance

Non-sterile Millex® filter units are compatible with ethylene oxide (EO) sterilization. Sterilization by any other method (e.g., gamma irradiation, electron beam, steam) is not recommended.

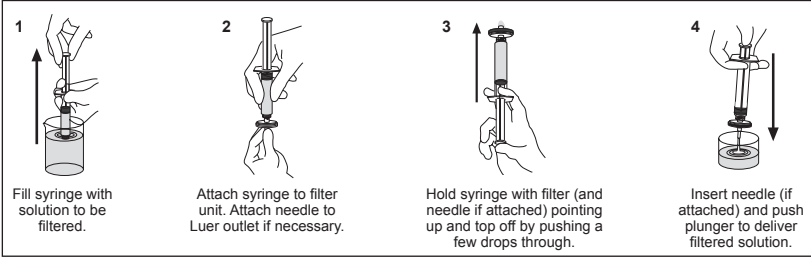Specifications

Materials	
Membrane	MF-Millipore™ membrane, mixed cellulose ester Millex®-OR filter unit: type GS, pore size 0.22 µm Millex®-GS filter unit: type GS, pore size 0.22 µm Millex®-HA filter unit: type HA, pore size 0.45 µm
Housing	
Dimensions	
Inlet to outlet	Millex®-OR filter unit: 23 mm (0.91 in.) Millex®-GS filter unit: 25 mm (0.98 in.) Millex®-HA filter unit: 25 mm (0.98 in.)
Diameter	29 mm (1.14 in.)
Filtration area	4 cm² (0.62 in²)
Temperature limit	45 °C (113 °F)
Pressure limit at 25 °C	5.2 bar (75 psi) inlet and differential
Filtration volume	1–100 mL
Hold-up volume	≤0.1 mL after air purge
Connections	Millex®-OR filter unit: Female Luer-Lok™ inlet, male Luer-Lok™ outlet Millex®-GS filter unit: Female Luer-Lok™ inlet, male Luer-slip outlet Millex®-HA filter unit: Female Luer-Lok™ inlet, male Luer-slip outlet
Flow rate at 2.1 bar (30 psi), 25 °C	Millex®-OR filter unit: ≥ 75 mL/min Millex®-GS filter unit: ≥ 75 mL/min Millex®-HA filter unit: ≥ 180 mL/min

Notice

The information in this document is subject to change without notice and should not be construed as a commitment by Merck Millipore Ltd. ("Millipore") or an affiliate. Neither Merck Millipore Ltd. nor any of its affiliates assumes responsibility for any errors that may appear in this document.

Technical Assistance

For more information, contact the office nearest you. In the U.S., call 1-800-MILLIPORE (1-800-645-5476). Outside the U.S., go to our web site at www.millipore.com/offices for up-to-date worldwide contact information. You can also visit the tech service page on our web site at www.millipore.com/techservice.

Standard Product Warranty

The applicable warranty for the products listed in this publication may be found at www.millipore.com/terms (within the "Terms and Conditions of Sale" applicable to your purchase transaction).

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