

43509 Legionella (GVPN) Selective Supplement

An antibiotic supplement recommended for selective isolation of Legionella species

Composition

(per vial, sufficient for 500 ml medium)

Glycine	1500 mg
Vancomycin hydrchloride	0.50 mg
Polymyxin B sulphate	40000 IU
Natamycin	20 mg

Directions

Rehydrate the contents of 1 vial aseptically with 10 ml of sterile distilled water. Mix well and aseptically add to 490 ml of sterile, molten Buffered Charcoal Yeast Extract Agar, Base (Cat. No. 86558) along with rehydrated contents of 1 vial of Legionella Supplement (Twin Pack; Cat. No. 89166).

Mix well and pour into sterile Petri plates.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

