



3050 Spruce Street  
Saint Louis, Missouri 63103 USA  
Telephone 800-325-5832 • (314) 771-5765  
Fax (314) 286-7828  
email: techserv@sia.com  
sigma-aldrich.com

## Product Information

### EGG YOLK TELLURITE EMULSION

Product Number **E 7774**

#### Product Description

Sterile stabilized tellurite emulsion of egg yolk recommended for identification of the *Staphylococcus* species.

#### Components

##### Item

Egg Yolk	30.0 mls
Sterile Saline	64.0 mls
Sterile 3.5% Potassium Tellurite Solution	6.0 mls

#### Preparation Instructions

Warm the refrigerated Egg Yolk Tellurite Emulsion to 45-50 °C and aseptically add 50 mls to 950 mls of sterile, molten B 1015 Baird Parker Agar Base. Mix well and pour into sterile petri plates.

#### Storage Temperature

Store at 2-8 °C.

**Precautions and Disclaimer** For laboratory use only.  
Not for drug, household or other uses.

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.