
MultiScreen₃₈₄-SEQ Filter Plate

Ready-to-use, automation-friendly, 384-well plate for dye terminator removal

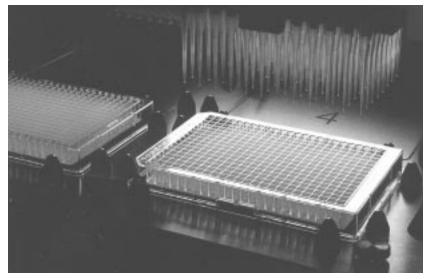

The MultiScreen₃₈₄-SEQ filter plate recovers highly purified sequencing products for long, accurate sequencing reads. The plates are fully automatable and provide a cost- and time-efficient alternative to ethanol precipitation for dye-terminator removal. They are ideal for use before electrokinetic injection or loading onto a slab gel-type DNA sequencer.

Designed for Automated Operations

The MultiScreen₃₈₄-SEQ filter plate is suited for use with most liquid handling systems. The filter plate operates by vacuum filtration enabling top-loading and recovery of sample. Our plate is designed with the needs of the high throughput laboratory in mind, meets SBS standards and is ideally suited for "walk away" automation.

Quick and Easy Protocol

Use of the MultiScreen₃₈₄-SEQ filter plate minimizes the procedural complexity involved in traditional precipitation or bind-and-elute methods. These plates eliminate the need for centifugation, time consuming incubations, ethanol-drying steps and manifold disassembly routines. Resuspension in electrokinetic injection buffer or water facilitates integration with DNA sequencers.

High Throughput Recovery of Even Miniaturized Sequencing Reactions

The MultiScreen₃₈₄-SEQ filter plate uses a size-selective membrane to separate sequencing products from contaminants. The plate retains sequencing products on the membrane surface, allowing salts and dye terminators to be filtered to waste. This timeand cost-effective method delivers highly purified sequencing products for long, accurate DNA sequencing reads.

MILLIPORE

Simply Wash, Reconstitute, and Recover



1. Introduce samples into wells



 Filter using MultiScreen₃₈₄ vacuum manifold for 6 minutes or until wells are empty. DNA is retained on the surface of the membrane while smaller contaminants are filtered to waste.



 Wash the sample once. Add resuspension buffer to each well (typically 10 – 25 μl) and recover purified DNA from the membrane surface by mixing or shaking. Retrieve purified samples by aspirition.

Features

- Automation friendly: Design and protocol saves time, money and improves reproducibility
- Efficient dye terminator & salt removal: Recovers highly purified sequencing products for long, accurate sequencing reads
- Eliminates dry-down step: Sequencing products may be resuspended in any aqueous electrokinetic injection buffer
- Saves reagents: plates suitable for use with 1/16 sequencing reactions

Performance

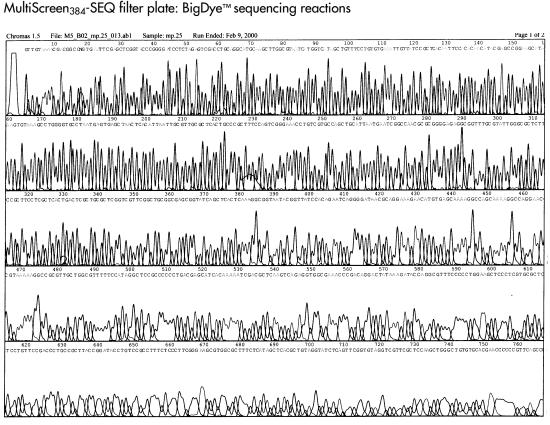


Figure 1. The MultiScreen₃₈₄-SEQ filter plate was used to purify BigDye sequencing reactions. Quarter reactions were cleaned up and injected onto the ABI Prism^M 3700. This example demonstrates the utility of the MultiScreen₃₈₄-SEQ filter plate for dye terminator removal, delivering long, high quality reads (Phred 20 \geq 627). Data courtesy of Duane Bartley, Johns Hopkins University.

MultiScreen₃₈₄-SEQ filter plate: ET Terminators

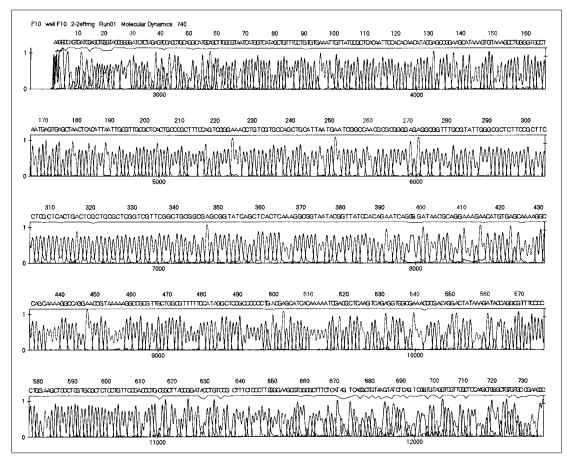


Figure 2. The MultiScreen₃₈₄-SEQ filter plate was used to purify ET terminator sequencing reactions. Quarter reactions were cleaned up and injected onto the APB MegaBACE[®] 1000 sequencer. Electrokinetic injection was performed out of water at 2 Kv for 10 seconds. Samples were run at 9 Kv for 120 minutes. This example further demonstrates the utility of the MultiScreen₃₈₄-SEQ filter plate for dye terminator removal, delivering long, high quality reads (Phred 20 \geq 607).

MultiScreen₃₈₄-SEQ filter plate

Phred 20 bases for 96 samples, MegaBACE 1000 sequencer, $_{\rm 1/8}$ reaction, plasmid template 54,445 Phred 20 bases for these 96 samples

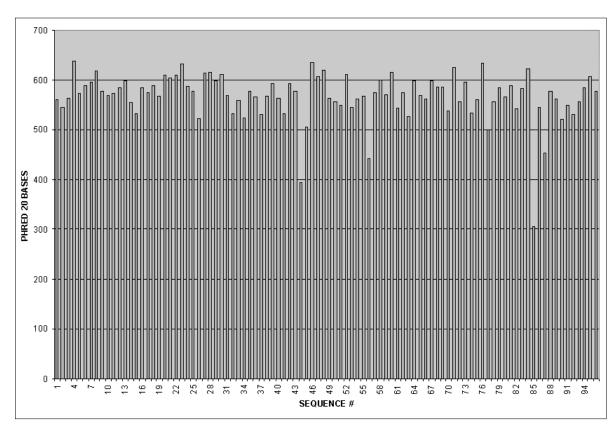


Figure 3. The MultiScreen₃₈₄-SEQ filter plate was evaluated for intra-plate reproducibility. Plasmid template was sequenced with ET Terminators (1/8 reactions) and MultiScreen₃₈₄-SEQ filter plates were used for dye terminator removal. The data show the reproducibility of read lengths from one quadrant of a 384 well plate. For these 96 samples, 54,445 Phred 20 bases were determined for an average read length of 567 bases. Similar experiments with BigDye sequencing chemistry yielded average Phred 20 read lengths of 630 bases (data not shown).

Specifications

Maximum well volume:130 μLWorking well volume:100 μLNon-recoverable volume:< 2 μL</td>Distance from bottom of plate to
membrane surface:3 mm

Materials of construction

Plastic housing: Clear Polystyrene Membrane: Proprietary

Automation Platform Recommendations

The MultiScreen₃₈₄-SEQ filter plate performs best with 96- or 384-tip liquid handlers capable of pipetting $\geq 25 \ \mu L$ /second or with plate shakers with 0.12 inch orbit at 1100 rpm. Other liquid handlers or shakers operating at slower speeds are equally effective, but resuspension times are longer.

Contact Millipore Technical Services or consult with one of the Millipore Application Scientists for help with your automation needs.

Chemical Compatibility

| Chemical | 1 | Chemical | |
|---------------|---|---------------------|---|
| EDTA | Е | Ethanol 100% | Q |
| Formamide 70% | Х | Ethanol 70% | E |
| Formamide 20% | E | Ethanol 20% | E |
| DMSO 100% | Х | Sodium Citrate | E |
| DMSO 50% | Х | 2-Pyrrolidinone 20% | Q |
| DMSO 20% | Q | | |

E= Excellcent Q= Questionable X= Not Recommended

Ordering Information

| Description | Pack Size | Catalogue No. |
|--|-----------|---------------|
| MultiScreen ₃₈₄ -SEQ filter plate | 10 | S384 SEQ 10 |
| | 50 | S384 SEQ 50 |
| MultiScreen ₃₈₄ Manifold | 1 | SAVM 384 01 |

Also available: Fully-automatable MultiScreen₃₈₄-PCR filter plate for one-step cleanup of PCR products. Request publication number PF1546EN00. Call the nearest Millipore office for information.

| MultiScreen ₃₈₄ -PCR filter plate | 10 | S384PCR10 |
|--|----|-----------|
| | 50 | S384PCR50 |

To Place an Order or Receive Technical Assistance

For additional information call your nearest Millipore office: In the U.S. and Canada, call toll-free 1-800-MILLIPORE (1-800-645-5476)In the U.S., Canada and Puerto

Rico, fax orders to 1-800-MILLIFX (1-800-645-5439)On the Internet http://www.millipore.com E-mail: tech service@millipore.com

pure e-commerce

Now you can buy Millipore products online @ www.millipore.com/purecommerce

MILLIPORE

Millipore and MultiScreen are trademarks of Millipore Corporation or an affiliated company MultiScreen₃₈₄-SEQ is patent pending. ABI Prism 3700 and BigDye are trademarks of PE Corporation or its subsidiaries. MegaBACE and DYEnamic are trademarks of Amersham Pharmacia BioTech, Inc. Millipore Lit. No. PF1545EN00 Printed in U.S.A. 10/00 00-185 Copyright 2000 Millipore Corporation or an affiliated company. All rights reserved.

Millipore Worldwide

AUSTRALIA C7ECH REPUBLIC Tel. 1 800 222 111 Tel. 02-2051 3841 02-2051 3842 or (02) 9888 8999 Fax 02-2051 4298 Fax (02) 9878 0788 AUSTRIA DENMARK Tel. (01) 877-8926 Tel. 70 10 00 23 Fax 70 10 13 14 Fax (01) 877-1654 IAPAN BALTIC COUNTRIES EASTERN EUROPE, C.I.S., Tel. +358 9 804 5110 AFRICA, MIDDLE EAST Fax +358 9 256 5660 AND GULF Analytical Division: **BELGIUM AND** Tel. (33) 3.88.38.9536 LUXEMBOURG Fax (33) 3.88.38.9539 Tel. (02) 726 88 40 BioProcess Division: Fax (02) 726 98 84 Tel. (43) 1 877-8926 Fax (43) 1 877-1654 BRA7II Laboratory Water Division: Tel. (33) 1.30.12.70.00 Tel. (011) 5548-7011 Fax (011) 5548-7923 Fax (33) 1.30.12.71.80 Fax 603-757 1711 CANADA Tel. 1-800-645-5476 FINLAND Tel. (09) 804 5110 Fax 1-800-645-5439 Fax (09) 256 5660 CHINA, PEOPLE'S **REPUBLIC OF** FRANCE Tel. (01) 30.12.70.00 Fax (01) 30.12.71.80 Beijing Tel. (8610) 8519-1250 (8610) 8519-1058 GFRMANY Fax (8610) 8519-1255 NORWAY Tel. (06196) 494-0 Guangzhou: Fax (06196) 43901 Tel. (8620) 8755-4021 Fax (8620) 8755-4350 HUNGARY POLAND Hong Kong: Tel. (852) 2803-9111 Tel. 1-205 9784 Fax 1-205 9792 Fax (852) 2513-0313 INDIA Shanghai Tel. (91) 80-839 46 57 Tel. (8621) 5306-9100 Fax (91) 80-839 63 45 Fax (8621) 5306-0838

ITALY Vimodrone (Milano): Tel. (02) 25.07.81 Fax (02) 26.50.324 Roma: Tel. (06) 52.03.600 Fax (06) 52.95.735

Tel. (03) 5442-9711 Fax (03) 5442-9736 Analytical 9737 BioProcess 9734 Lab Water

KOREA Tel. (822) 551-0990 Fax (822) 551-0228 MALAYSIA Tel. 603-757 1322

MEXICO Tel. (525) 576-9688 Fax (525) 576-8706

THE NETHERIANDS Tel. 076-5022000 Fax 076-5022436

Tel. 22 67 82 53 Fax 22 66 04 60

Tel. 22-669 12 25 22-663 70 31 Fax 22-663 70 33

PUERTO RICO Tel. (787) 273-8495 Fax (787) 747-6553

SINGAPORE Tel. (65) 842 1822

Fax (65) 842 4988

SPAIN AND PORTUGAL

Madrid: Tel. 917 283 960 Fax 917 292 909 Barcelona: Tel. 934 525 530 Fax 934 516 048

SWEDEN Tel. 08-628 6960 Fax 08-628 6457

SWITZERLAND

Volketswil: Tel. (01) 908-30-60 Fax (01) 908-30-80 Lausanne: Tel. (021) 641 2550 Fax (021) 641 2551

TAIWAN

Taipei: Tel. (886-2) 700-1742 Fax (886-2) 755-3267 Hsin Chu City: Tel. (886-3) 571-0178 Fax (886-3) 572-9520

U.K. AND IRELAND Tel. (01923) 816375 Fax (01923) 818297

U.S.A

Tel. (781) 533-6000 Fax (781) 533-3110

IN ALL OTHER

COUNTRIES Millipore Intertech (U.S.A.) Tel. +1 (781) 533-8622 Fax +1 (781) 533-8630