

Product Information

Phalloidin, Tetramethylrhodamine B isothiocyanate

Sequence from *Amanita phalloides* (synthetic peptide sequence)**P1951**

Product Description

Molecular Formula: C₆₀H₇₀N₁₂O₁₅S₂


Molecular Weight: 1231.40

Synonyms: Phalloidin-TRITC, TRITC-Phalloidin, Rhodamine-Phalloidin, Rhodamine Phalloidin-TRITC

Molar Extinction Coefficient:¹ 80,000 (545 nm)Excitation wavelength:¹ 540-545 nmEmission wavelength:¹ 570-573 nm

Storage temperature: -20 °C

Structure:

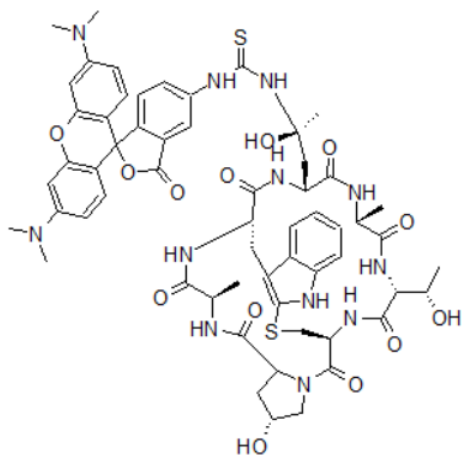

Phalloidin is a fungal toxin that occurs naturally in the poisonous mushroom *Amanita phalloides*.² Phalloidin toxicity is attributed to the ability to bind F actin in liver and muscle cells. As a result of binding phalloidin, actin filaments become strongly stabilized. Phalloidin has been found to bind only to polymeric and oligomeric forms of actin, and not to monomeric actin.³ The dissociation constant of the actin-phalloidin complex has been determined to be on the order of 3×10^{-8} M.⁴

Fluorescent conjugates of phalloidin, such as FITC (fluorescein isothiocyanate) or TRITC (tetramethylrhodamine B isothiocyanate) conjugates, have been used to probe actin filaments for histological applications.³⁻¹⁰ Some structural features of phalloidin are required for the binding to actin.³ However, the side chain of amino acid 7 (γ - δ -dihydroxyleucine) is accessible for chemical modifications without appreciable loss of affinity for actin. TRITC-phalloidin is considered to be less susceptible to photobleaching than FITC-phalloidin.⁶ The binding positions of phalloidin and of TRITC-phalloidin to F-actin have been studied by X-ray crystallography.¹¹

Precautions and Disclaimer

This product is for R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Storage/Stability

It is recommended to store this product at -20 °C.

Preparation Instructions

This product is tested for solubility in methanol at 2 mg/mL. In general, solutions of phalloidin compounds should be prepared fresh and protected from light whenever possible. One publication reports preparation of a ~0.5 mM stock solution of TRITC-phalloidin in DMSO.¹²

Several publications cite storage of stock solutions of TRITC-phalloidin in methanol in the dark at -20 °C,^{13,14} although we have ourselves not tested the solution stability of TRITC-phalloidin.

Procedure

The following procedure may serve as a general guideline for staining cells.¹⁵ Final staining solutions in aqueous physiological buffers have a phalloidin concentration range of 0.1-100 µM, with corresponding incubation times of 15 minutes to 72 hours.

1. Cells are washed with phosphate buffered saline (PBS).
2. Cells are fixed for 5 minutes in 3.7% formaldehyde solution in PBS, then washed extensively in PBS.
3. Cells may be dehydrated with acetone, permeabilized with 0.1% TRITON™ X-100 in PBS and washed again in PBS.
4. Cells are stained with a 50 µg/mL fluorescent phalloidin conjugate solution in PBS (containing 1% DMSO from the original stock solution) for 40 minutes at room temperature.
5. Wash several times with PBS to remove unbound phalloidin conjugate.

References

1. Faulstich, H. *et al.*, *J. Muscle Res. Cell Motil.*, **9(5)**, 370-383 (1988).
2. Abul-Haj, S. K. *et al.*, *New Eng. J. Med.*, **269**, 223-227 (1963).
3. Heidecker, M. *et al.*, *Biochemistry*, **34(35)**, 11017-11025 (1995).
4. Wulf, E. *et al.*, *Proc. Nat. Acad. of Sci. USA*, **76(9)**, 4498-4502 (1979).
5. Small, J. V. *et al.*, *J. Cell Sci.*, **89(Pt 1)**, 21-24 (1988).
6. Faulstich, H. *et al.*, *Exp. Cell Res.*, **144(1)**, 73-82 (1983).
7. Lawson, M. A., and Maxfield, F. R., *Nature*, **377(6544)**, 75-79 (1995).
8. Pinaev, G. *et al.*, *FEBS Lett.*, **369(2-3)**, 144-148 (1995).

9. Mahaffy, R. E., and Pollard, T. D., *Biochemistry*, **47(24)**, 6460-6467 (2008).
10. Wang, Y.-L., *J. Cell. Biol.*, **105(6 Pt 1)**, 2811-2816 (1987).
11. Oda, T. *et al.*, *Biophys. J.*, **88(4)**, 2727-2736 (2005).
12. Theesfeld, C. L. *et al.*, *J. Cell Biol.*, **146(5)**, 1019-1032 (1999).
13. Sund, S. E., and Axelrod, D., *Biophys J.*, **79(3)**, 1655-1659 (2000).
14. González-Gutiérrez, A. *et al.*, *Front. Plant Sci.*, **11**, 384 (2020).
15. Waggoner, A. *et al.*, *Methods Cell Biol.*, **30**, 449-478 (1989).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2009-2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.
P1951pis Rev 06/23 CS,RGB,RD,GCY,MAM

