

## Product Information

### *Instruction Guide for the*

# MINI-CENTRIFUGE

Congratulations on your purchase of a Sigma Personal Mini-Centrifuge!


**Catalog #C5730 – Mini-centrifuge – 220V**

**Catalog #C5065 – Mini-centrifuge – 110V**


Following are some recommended operating procedures:

### Operation


- After unpacking, place the Mini-Centrifuge on a flat, clear surface.
- To open, press firmly down on release tab while lifting dome.
- Insert tubes into rotor completely. Be sure that each tube is securely closed and of equal volume. If necessary, add dummy-tubes to balance.
- Load rotor as indicated below. Properly balanced loads are essential for safe and quiet operation.

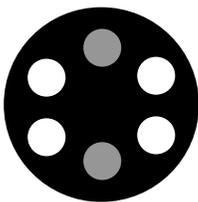


**2 Tubes**

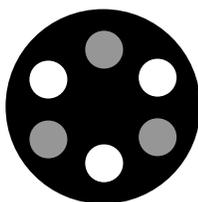

**3 Tubes**

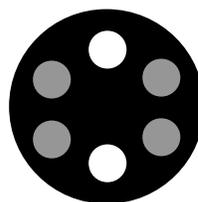

**4 Tubes**

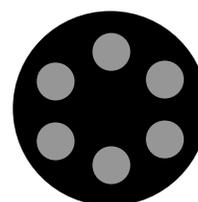

**6 Tubes**

- Close dome securely and activate side switch. Device will not operate with dome open.
- To stop device, deactivate side switch. Wait until device has completely stopped to release dome. Remove samples.

## Safeguard

- Do not immerse device in water. Injury due to shock or fire may result.
- Do not operate device with dome open, or open dome before rotor has ceased spinning.
- Do not insert non-standard tubes or other foreign articles into rotor.
- Do not operate under uneven loads. Destabilized operation may result in damage to the device or injury to the user.
- Store item at room temperature in a dry area. Do not expose to sun, moisture or extreme temperatures for prolonged periods of time.



**Our one-year warranty assures you reliable quality at an unbeatable value!**

Sigma-Aldrich • P.O. Box 14508 • St. Louis, Missouri 63178  
Phone: (800) 325-3010 • Fax: (800) 325-5052 • [www.sigma-aldrich.com](http://www.sigma-aldrich.com)