Vented Millex®-GS Sterilizing Filter Unit

SLGSV255F

- Single use only
 - Sterile
 - Non-pyrogenic
- Introduction**
- The Vented Millex®-GS filter unit is a sterilizing filter for aqueous solutions dispensed with a syringe. It will remove microorganisms, particles, precipitates, and undissolved powders larger than 0.22 micron (µm). The sterile Vented Millex®-GS filter unit is non-pyrogenic and non-toxic. This single-use product consists of a 0.22 µm membrane filter sealed in a plastic housing. A 0.03 µm hydrophobic filter in the dome prevents air locking by automatically venting air introduced upstream. Typical applications include the sterile filtration of antibodies and critical care solutions (TPN without lipids or additives).
- Chemical Compatibility**
- The Vented Millex®-GS filter unit is compatible with most aqueous solutions. It may be used to filter the agents listed below. This guide has been developed from technical publications, materials suppliers, and laboratory tests, and is believed to be reliable. However, because of variability in temperature, concentrations, duration of exposure, and other factors outside of our control which may affect the use of the unit, no warranty is given or is to be implied with respect to such information. Agents not listed below should be tested with the Vented Millex®-GS filter prior to use.

Acetic acid (5%)	Butyl alcohol	Hypo (photo)	Pentane
Aliphatic ethers	Freon® solvent, TF or PCA	Isobutyl alcohol	Petroleum based oils
Ammonium hydroxide (6 N)	Glycerine (glycerol)	Isopropyl alcohol	Phenol (0.5%)
Benzyl alcohol (1%)	Helium	Kerosene	Silicone oils
Boric acid	Hexane	Mineral spirits	
Brine (sea water)	Hydrogen (gas)	Nitrogen	

- Directions for Use**
- WARNINGS:**
- To ensure sterility, do not use this product if the package is damaged.
 - Do not use with syringes smaller than 10 mL because pressures in excess of the maximum pressure rating may be reached, potentially causing damage to the filter unit and/or personal injury.
 - Ne pas utiliser avec des seringues de moins de 10 ml car des pressions supérieures à la pression nominale maximale pourraient être atteintes et celles-ci pourraient endommager le matériel et causer des blessures graves.
- CAUTIONS:**
- Do not resterilize or reuse this filter unit, as Merck Millipore Ltd. cannot assure the sterility, integrity, and performance beyond a single use.
 - Do not use the Vented Millex®-GS unit at temperatures above 45 °C (113 °F).
 - Do not use the Vented Millex®-GS unit to filter emulsions or suspensions.
 - Do not use the Vented Millex®-GS unit to filter 5 mg or less of active drug materials unless binding studies have been performed.

1	2	3	4	5
Fill syringe with solution to be filtered.	Aseptically remove cover from package.	Attach syringe to filter unit and remove assembly from package. Attach needle to Luer-slip outlet if necessary.	Hold syringe with filter (and needle if attached) pointing up and top off plunger to deliver filtered solution.	Insert needle (if attached) and push down plunger to deliver filtered solution.

Specifications		
Materials		
Filter membrane	Mixed cellulose ester, type GS, pore size 0.22 µm	
Vent membrane	Hydrophobic PTFE (polytetrafluoroethylene), 0.03 µm	
Housing	PVC (polyvinyl chloride)	
Dimensions		
Inlet to outlet	27 mm (1.06 in.)	
Diameter	29 mm (1.14 in.)	
Filtration area	4 cm² (0.62 in²)	
Temperature limit	45 °C (113 °F)	
Pressure limit at 25 °C	5.2 bar (75 psi) inlet and differential	
Filtration volume	1–100 mL	
Hold-up volume	≤ 0.1 mL after air purge	
Sterilization method	Ethylene oxide gas	
Connections	Female Luer-Lok™ inlet, male Luer-slip outlet	
Flow rate at 0.7 bar (10 psi), 25 °C	≥ 30 mL/min	
Notice		
The information in this document is subject to change without notice and should not be construed as a commitment by Merck Millipore Ltd. ("Millipore") or an affiliate. Neither Merck Millipore Ltd. nor any of its affiliates assumes responsibility for any errors that may appear in this document.		
Technical Assistance		
For more information, contact the office nearest you. In the U.S., call 1-800-MILLIPORE (1-800-645-5476). Outside the U.S., go to our web site at www.millipore.com/offices for up-to-date world-wide contact information. You can also visit the tech service page on our web site at www.millipore.com/techservice .		
Standard Product Warranty		
The applicable warranty for the products listed in this publication may be found at www.millipore.com/terms (within the "Terms and Conditions of Sale" applicable to your purchase transaction).		

© 2012 EMD Millipore Corporation. All rights reserved.

The M logo is a trademark of Merck KGaA, Darmstadt, Germany.

Millipore and Millex are registered trademarks of Merck KGaA.

Freon is a registered trademark of E. I. du Pont de Nemours and Co.

Luer-Lok is a trademark of Becton, Dickinson and Co.

PR04005, Rev. A, 02/12

English	Ελληνικά	Magyar
Français	Nederlands	Česky
Italiano	Dansk	Polski
Deutsch	Svenska	日本語
Español	Latviski	
Português	Lietuviškai	

English	Ελληνικά	Magyar
Français	Nederlands	Česky
Italiano	Dansk	Polski
Deutsch	Svenska	日本語
Español	Latviski	
Português	Lietuviškai	

1	2	3	4	5
Refill the syringe with the solution to be filtered.	Remove aseptically the cap from the package.	Attach the syringe to the filter unit and remove the assembly from the package. Attach the needle to the Luer-lock outlet if necessary.	Hold the syringe with the filter (and needle if attached) pointing up and push the plunger down to deliver the filtered solution.	Insert the needle (if attached) and push down the plunger to deliver the filtered solution.

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PR04005, Rev.A.ind 1

Vented Millex®-GS Sterilizing Filter Unit

Cat. No. SLGSV255F (50/box)

- Single use only
 - Sterile
 - Non-pyrogenic
- Introduction**
- The Vented Millex®-GS filter unit is a sterilizing filter for aqueous solutions dispensed with a syringe. It will remove microorganisms, particles, precipitates, and undissolved powders larger than 0.22 micron (µm). The sterile Vented Millex®-GS filter unit is non-pyrogenic and non-toxic. This single-use product consists of a 0.22 µm membrane filter sealed in a plastic housing. A 0.03 µm hydrophobic filter in the dome prevents air locking by automatically venting air introduced upstream. Typical applications include the sterile filtration of antibodies and critical care solutions (TPN without lipids or additives).

Acetic acid (5%)	Butyl alcohol	Hypo (photo)	Pentane
Aliphatic ethers	Freon® solvent, TF or PCA	Isobutyl alcohol	Petroleum based oils
Ammonium hydroxide (6 N)	Glycerine (glycerol)	Isopropyl alcohol	Phenol (0.5%)
Benzyl alcohol (1%)	Helium	Kerosene	Silicone oils
Boric acid	Hexane	Mineral spirits	
Brine (sea water)	Hydrogen (gas)	Nitrogen	

- Directions for Use**
- WARNINGS:**
- To ensure sterility, do not use this product if the package is damaged.
 - Do not use with syringes smaller than 10 mL because pressures in excess of the maximum pressure rating may be reached, potentially causing damage to the filter unit and/or personal injury.
 - Ne pas utiliser avec des seringues de moins de 10 ml car des pressions supérieures à la pression nominale maximale pourraient être atteintes et celles-ci pourraient endommager le matériel et causer des blessures graves.
- CAUTIONS:**
- Do not resterilize or reuse this filter unit, as Merck Millipore Ltd. cannot assure the sterility, integrity, and performance beyond a single use.
 - Do not use the Vented Millex®-GS unit at temperatures above 45 °C (113 °F).
 - Do not use the Vented Millex®-GS unit to filter emulsions or suspensions.
 - Do not use the Vented Millex®-GS unit to filter 5 mg or less of active drug materials unless binding studies have been performed.

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